

J-1000 series

CD Spectrometer

Software Manual

Spectra Manager™ Ver2.5 CFR



Preface

This instruction manual serves as a guide for using this software. It is intended to instruct first-time users on how to properly use the software, and to serve as a reference for experienced users.

Before using the software, read this instruction manual carefully, and make sure you fully understand its contents. This manual should be easily accessible to the operator at all times during software operation. When not using the software, keep this manual stored in a safe place. Should this instruction manual be lost, order a replacement from your local JASCO distributor.

Servicing

Contact your local JASCO distributor for instrument servicing. In addition, contact your JASCO distributor before moving the instrument to another location. Consumable parts should be ordered according to part number from your local JASCO distributor. If a part number is unknown, give your JASCO distributor the model name and serial number of your instrument.

Notices

- (1) JASCO shall not be held liable, either directly or indirectly, for any consequential damage incurred as a result of product use.
- (2) Prohibitions on the use of JASCO software
 - Copying software for purposes other than backup
 - Transfer or licensing of the right to use software to a third party
 - Disclosure of confidential information regarding software
 - Modification of software
 - Use of software on multiple workstations, network terminals, or by other methods (not applicable under a network licensing agreement concluded with JASCO)
- (3) The contents of this manual are subject to change without notice for product improvement.
- (4) This manual is considered complete and accurate at publication.
- (5) This manual does not guarantee the validity of any patent rights or other rights.
- (6) If a JASCO software program has failed causing an error or improper operation, this may be caused by a conflict from another program operating on the PC. In this case, take corrective action by uninstalling the conflicting product(s).
- (7) *Windows* is a registered trademark of *Microsoft Corporation* in the United States and other countries. In general, company names and product names are trademarks or registered trademarks of the respective companies. However, the TM and [®] marks are not used in all cases in this manual.
- (8) JASCO and the JASCO logo are registered trademarks of JASCO Corporation in Japan and other countries.

Limited Warranty

Products sold by JASCO, unless otherwise specified, are warranted for a period of one year from the date of shipment to be free of defects in materials and workmanship. If any defects in the product are found during this warranty period, JASCO will repair or replace the defective part(s) or product free of charge.

THIS WARRANTY DOES NOT APPLY TO DEFECTS RESULTING FROM THE FOLLOWING:

- (1) IMPROPER OR INADEQUATE INSTALLATION
- (2) IMPROPER OR INADEQUATE OPERATION, MAINTENANCE, ADJUSTMENT OR CALIBRATION
- (3) UNAUTHORIZED MODIFICATION OR MISUSE
- (4) USE OF CONSUMABLE PARTS NOT SUPPLIED BY AN AUTHORIZED JASCO DISTRIBUTOR
- (5) CORROSION DUE TO THE USE OF IMPROPER SOLVENTS, SAMPLES, OR DUE TO SURROUNDING GASES
- (6) ACCIDENTS BEYOND JASCO'S CONTROL, INCLUDING NATURAL DISASTERS

This warranty does not cover the consumable parts listed below:

- (1) Tungsten lamp, and other light sources
- (2) Mirrors in the light source section, and cell windows
- (3) Fuses, batteries, glassware, chart paper and ink

THE WARRANTY FOR ALL PARTS SUPPLIED AND REPAIRS PROVIDED UNDER THIS WARRANTY EXPIRES ON THE WARRANTY EXPIRATION DATE OF THE ORIGINAL PRODUCT. FOR INQUIRIES CONCERNING REPAIR SERVICE, CONTACT YOUR JASCO DISTRIBUTOR AFTER CONFIRMING THE MODEL NAME AND SERIAL NUMBER OF YOUR INSTRUMENT.

JASCO Corporation
2967-5, Ishikawa-machi, Hachioji
Tokyo 192-8537
Japan

Notation Used

The following notational conventions are used throughout this manual:

General Notation

Notation	Meaning
[Measurement] menu [Parameters...] command	Names of menus, commands, and text boxes are enclosed in square brackets '[]', followed by a description indicating whether the function is a menu, command, text box, etc. Shortcut keys used to select menus or commands are underlined.
<OK>, <Cancel>	Names of buttons are enclosed in angular brackets '< >'.

Keyboard Operations

Notation	Meaning
Shift Ctrl	The key is enclosed in a square and shown in boldface.
Alt , F	Keys that are to be pressed in succession are separated by commas. In the example shown on the left, the Alt key is to be pressed and released, followed by the F key.
Shift + →	Keys that are pressed simultaneously are separated by a "plus" sign. In the example shown on the left, press the → key while holding down the Shift key.

Mouse Operations

Notation	Meaning
Point	Move the mouse pointer to the specified item.
Click	Quickly press and release the mouse button.
Double-click	Click the mouse button twice in rapid succession.
Drag	Point to an item, click and hold down the mouse button. Move the mouse with the button held down, and release the button when the pointer is in the desired position.

Contents

Preface.....	i
Servicing	i
Notices.....	ii
Limited Warranty	iii
Notation Used	iv
Contents	v
1. Introduction	1
1.1 How to Use This Manual.....	1
1.2 FDA 21CFR Part 11.....	2
1.3 Features of [Spectra Manager] Compatible with FDA 21CFR Part 11	2
1.4 Overview of [Spectra Manager].....	6
2. Starting and Exiting Programs and the [Spectra Manager] Program Reference	7
2.1 Starting the program	7
2.1.1 Turning on the CD Spectrometer	7
2.1.2 Starting the PC and Windows and Registering a J-1000 Series CD spectrometer	8
2.1.3 Starting [Spectra Manager CFR].....	23
2.1.4 Components of the Program Window	25
2.2 Exiting the Program	27
2.2.1 Exiting [Spectra Manager CFR]	27
2.2.2 Turning off the PC and CD Spectrometer	28
2.3 [Spectra Manager CFR] Program Reference	28
3. Accessory Registration Reference.....	30
3.1 Installing/Attaching an Accessory	30
3.2 Precautions for Peltier Thermostatic Cell Holder.....	31
3.2.1 Removing the Peltier Thermostatic Cell Holder.....	32
3.3 Registering Accessories	33
3.3.1 Registering an Automatically-Recognized Accessory	33
3.3.2 Registering an Accessory Requiring Manual Registration.....	33
3.4 Actions of [Spectra Manager] When Recognizing an Accessory	36
3.4.1 When an Automatically-Recognized Accessory is Attached.....	36
3.4.2 When the Automatically-Recognized Accessory is Removed	37
3.5 Status Display for the CD Spectrometer	38
4. Introduction to the [Spectra Measurement] Program	39
4.1 Procedures for the [Spectra Measurement] Program	39
4.2 Starting the [Spectra Measurement] Program	40
4.3 Specifying and Saving Measurement Parameters.....	41
4.4 Reviewing and Approving the Parameter File	48
4.5 Measuring the Sample	50
4.6 Saving Spectra	52
4.7 Printing Results	53

4.8	Exiting the Program	54
5.	Introduction to the [Time Course Measurement] Program.....	55
5.1	Procedures for the [Time Course Measurement] Program	55
5.2	Starting the [Time Course Measurement] Program	56
5.3	Specifying and Saving Measurement Parameters.....	57
5.4	Reviewing and Approving the Parameter File	64
5.5	Measuring the Sample	66
5.6	Saving Time-Course Data.....	68
5.7	Printing Results	69
5.8	Exiting the Program	70
6.	Measurement Program Reference	71
6.1	Overview of Measurement Programs.....	71
6.2	[File] Menu.....	76
6.2.1	[Save Data...].....	76
6.2.2	[Send to Analysis].....	77
6.2.3	[Open Parameters...]	78
6.2.4	[Save Parameters...]	79
6.2.5	[Open Baseline...]	80
6.2.6	[Save Baseline...].....	82
6.2.7	[Print Setup...].....	83
6.2.8	[Exit]	83
6.3	[Measure] Menu.....	84
6.3.1	[Cancel]	84
6.3.2	[Sequence Measurement].....	84
6.3.3	[Sample Measurement].....	84
6.3.4	[Baseline Measurement]	85
6.3.5	[Blank Measurement].....	85
6.3.6	[Dark Measurement]	86
6.3.7	[Parameters...].....	86
6.4	[Control] menu	87
6.4.1	[Zero Clear]	87
6.4.2	[Shutter].....	87
6.4.3	[Move Wavelength...].....	87
6.4.4	[Detector ...].....	88
6.4.5	[Light Source...]	89
6.4.6	[Nitrogen Flowrate...]	89
6.4.7	[Select Detector...]	89
6.4.8	[Select Accessory...]	90
6.4.9	[Accessory Model Number].....	91
6.5	[View] Menu.....	92
6.5.1	[Scales...].....	92
6.5.2	[Patterns...]	93
6.5.3	[Fonts...]	94
6.5.4	[Gridlines...]	95
6.5.5	[Styles...].....	96
6.5.6	[Overlay].....	97
6.5.7	[Select Graph...].....	97
6.5.8	[Information Bar]	97

6.5.9	[Accessory Bar]	97
6.5.10	[Toolbars]	97
6.5.11	[Status Bar].....	97
6.5.12	[View Mode].....	98
6.5.13	[Graph View].....	99
6.6	[Settings] Menu.....	99
6.6.1	[Default Parameters...].....	99
6.6.2	[Customize Toolbar...].....	100
6.6.3	[Monitor Bar...].....	102
6.7	[Help] Menu	102
6.7.1	[Contents].....	102
6.7.2	[Search Topic]	102
6.7.3	[About...].....	102
7.	[Spectra Measurement] Program Reference.....	103
7.1	[Parameters...].....	103
7.1.1	[General] tab.....	104
7.1.2	[Control] tab.....	108
7.1.3	[Information] tab.....	109
7.1.4	[Data] tab.....	110
8.	[Time Course Measurement] Program Reference	112
8.1	[Parameters...].....	112
8.1.1	[General] tab.....	113
8.1.2	[Control] tab.....	116
8.1.3	[Information] tab.....	117
8.1.4	[Data] tab.....	118
9.	[Data Monitor] Program Reference.....	120
9.1	[Parameters...]	121
9.1.1	[Parameters].....	121
10.	Appendix.....	124
10.1	Setting the Measurement Parameters.....	124
10.1.1	Measurement Wavelength Range.....	124
10.1.2	Data Acquisition Interval (Resolution)	124
10.1.2.1	Spectra measurement.....	124
10.1.2.2	Time-course measurement	124
10.1.3	Scanning Speed	124
10.1.4	Number of Accumulations.....	125
10.1.5	Response (D.I.T.)	125
10.1.5.1	Spectra measurement.....	125
10.1.5.2	Time-course measurement	126
10.1.6	Bandwidth and slit width	126

1. Introduction

1.1 How to Use This Manual

This manual is for the JASCO J-1000 series CD spectrometer measurement program that runs on *Microsoft Windows*. The J-1000 series CD spectrometer measurement program consists of [Spectra Manager CFR-Compliant Program], which conforms to FDA regulation 21 CFR Part 11 concerning electronic records and electronic signatures, and [Spectra Manager v.2-Compliant Program], which does not conform to these standards. This instruction manual refers to [Spectra Manager CFR-Compliant Program] and [Spectra Manager v.2-Compliant Program] as [CFR Version Program] and [Non-CFR Version Program], respectively.

This chapter describes the structure and use of this manual. The manual consists of 7 chapters including this one. An explanation of each chapter is given below. For this manual, *Microsoft Windows* is referred to as *Windows* and a personal computer as a PC.

1. Introduction

This chapter describes how to use this manual and gives an overview of each program and menu. Read this chapter first.

2. Starting and Exiting Programs and the [Spectra Manager] Program Reference

This chapter describes detailed procedures ranging from turning on/off the CD spectrometer and PC to starting/exiting the programs. Refer to Chapter 3 for details about program operations. In addition, this chapter provides the [Spectra Manager] program reference.

3. Accessory Registration Reference

This chapter describes how to register accessories used with the CD spectrometer.

4. Introduction to the [Spectra Measurement] Program

This chapter describes the [Spectra Measurement/Analysis] programs with examples of basic procedures. Users unfamiliar with *Windows* or first-time users of the CD spectrometer should follow the procedures described in this chapter for a general overview of how to operate the program.

5. Introduction to the [Time Course Measurement] Programs

This chapter describes the [Time Course Measurement] programs with examples of basic procedures.

6 to 9. Measurement Program References

These chapters describe common toolbar and menu commands for the [Spectra Measurement], [Time Course Measurement] and [Data Monitor] programs, in addition to the parameters used in these programs. Read the relevant chapters as required.

6. Toolbar and menu reference.
7. [Spectra Measurement] program reference
8. [Time Course Measurement] program reference
9. [Data Monitor] program reference

10. Appendix

This chapter provides useful suggestions for specifying parameters for CD measurements.

1.2 FDA 21CFR Part 11

FDA 21 CFR Part 11 is a set of regulations regarding digital archiving of data and data records for GLP, GCP, and GMP procedures. These regulations cover a number of areas, including: 1) access control and electronic signature requirements for data recorded by any computer-controlled analytical system in which the results are digitally archived; 2) the provision of security functions that can only be accessed by authorized personnel to ensure the security and integrity of data; and 3) a data auditing mechanism with the automatic creation of an Audit Trail to maintain a record of any creation, modification or deletion of instrument data.

1.3 Features of [Spectra Manager] Compatible with FDA 21CFR Part 11

JASCO Spectra Manager™ is an integrated software package that acts as a common platform for the range of JASCO analytical instruments. The Spectra Manager™ software provides functionality ranging from analytical instrument control and spectral data processing to sophisticated and specialized data analysis programs. JASCO Spectra Manager™ Ver.2.5 CFR provides security and auditing functionality that ensures the security, integrity and confidentiality of electronic records by enforcing the use of electronic signatures and other security measures as described in the regulations outlined in 21 CFR Part 11.

In addition, in order to ensure strict data integrity, functions are provided for electronically recording intermediate results even in the case where a measurement is stopped, either for a single spectrum or time-dependent spectral data. Also, saving of data before and after changing metadata such as the sample name or comments is strictly enforced.

JASCO Spectra Manager™ CFR consists of three core modules: Administrative Tools and Security Manager modules; intrinsic components of the Spectra Manager™ CFR software; and a database, which is an external component to the Spectra Manager™ CFR software. In addition, there are two broad classes of add-on modules: instrument drivers, which are spectrometer, or accessory-specific; and spectral analysis modules, which provide a range of sophisticated analytical and data processing tools. These add-on modules integrate seamlessly with the three core components of the Spectra Manager system to provide compliance with 21 CFR Part 11 regulations over the entire range of JASCO products. Figure

1.1 illustrates the system organization.

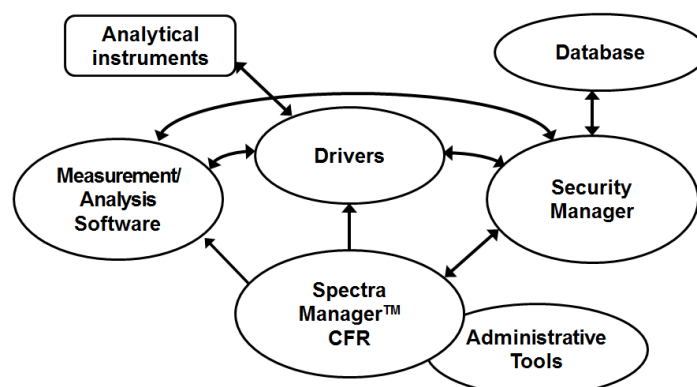


Figure 1.1 Structure of Spectra Manager™ CFR

All of the security management functions of JASCO Spectra Manager™ CFR are accessed through the Administrative Tools interface. The Security Manager components of the system are not directly accessible by a user, and thus the operation of these modules is not described. For secure management of data and data records, JASCO Spectra Manager™ CFR has two levels of security: system-level security and workgroup-level security.

When a user is registered in JASCO Spectra Manager™ CFR, the user must be assigned to one of four possible System Access Levels: “Administrators”, “Power Users”, “Users”, or “Limited Users”. The level of control that a user has over the management functions of the software system is determined by the System Access Level assigned to that user. Table 1.1 outlines the different rights assigned to the different types of users depending on their System Access Levels. Users with an “Administrators” system access level are referred to as “[Spectra Manager CFR] administrators”. In contrast, users belonging to the “Administrators” group in *Windows* are referred to as “*Windows* administrators”.

Note: An “Administrators” access level has no relation to “Windows administrator” and means administrator privileges for the [Spectra Manager CFR] program.

Table 1.1 Rights of each System Access Level

	Administrators	Power Users	Users	Limited Users
Display/modify system policy	OK			
Add/delete users	OK			
Enable/disable user account	OK			
Modify user properties (incl. System Access Level)	OK			
Add a workgroup	OK			
Enable/disable a workgroup	OK			
Modify workgroup users	Workgroup Managers only			
Add/delete instruments	OK			
Modify instrument configuration	OK			
Add/delete programs	OK			
Display log management	OK	OK		

Workgroups correspond to actual groupings of users within the organization and may represent divisions such as projects, departments, or research groups. Each workgroup has a set of users, instruments, accessories, and analysis programs associated with it; only users registered with a specific workgroup may access the instruments and programs of that workgroup. Within a workgroup, users are granted access as either a “Manager”, “Analyst”, or “Operator”, with authorities to modify instrument and analysis settings and use resources within the workgroup determined by the Workgroup User Rights granted. Selectable Workgroup User Rights vary depending on the system access level of the user. Table 1.2 outlines the relationship between the System Access Levels and Workgroup User Rights.

Table 1.2 Qualified access levels by Workgroup User Rights

	Administrators	Power Users	Users	Limited Users
Managers	OK	OK		
Analysts	OK	OK	OK	
Operators	OK	OK	OK	OK

By employing a two-level security model, global security authorization is independent of instrument, accessory, and application authorization, resulting in a highly flexible security system. Table 1.3 outlines the different rights assigned to the different types of users depending on their Workgroup User Rights.

Table 1.3 Rights of each Workgroup User Rights

	Managers	Analysts	Operators
Register/unregister workgroup users	OK		
Change Workgroup user authorities	OK		
Register/unregister instruments for the workgroup	OK		
Register/unregister programs for the workgroup	OK		
Use a registered instrument	OK	OK	OK
Use a registered instrument application	OK	OK	OK
Modify measurement parameters	OK	OK	
Use a registered analysis program	OK	OK	OK
Modify analysis parameters	OK	OK	
Electronically sign data	OK	OK	OK

Figure 1.2 shows the typical operational flow of the JASCO Spectra Manager™ CFR. First, a user is created (registered) and their System Access Level is assigned using Administrative Tools. A Workgroup is created (registered), and individual users and analytical instruments are assigned to the Workgroup. Detailed access to instrument control and analysis programs can be set at the User and Workgroup levels. Access control levels as required by FDA 21 CFR Part 11 regulations are determined by the Administrative Tools settings.

Routine analyses can then be conducted by the operator using the measurement/analysis programs in JASCO Spectra Manager™ CFR. Since all levels of access are managed by

Security Manager, individual users are only able to execute procedures that are allowed by their assigned access level. Access control ranges from declaring measurement parameters to conducting measurements and performing data analysis.

Analytical data obtained using any of the procedures described above can be saved as an electronic record with an electronic signature.

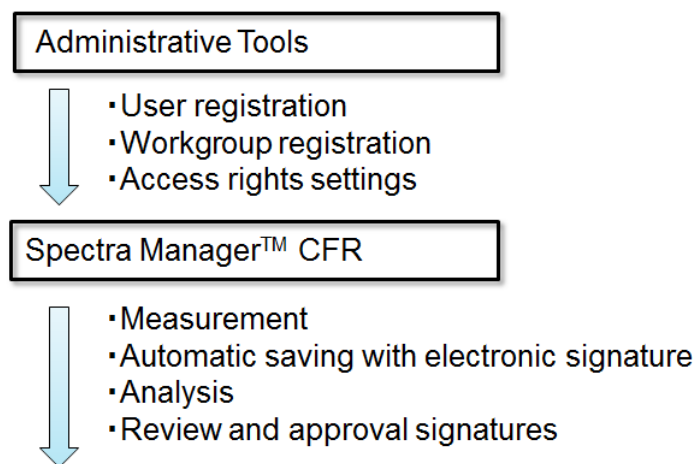


Figure 1.2 Operational flow

1.4 Overview of [Spectra Manager]

[Spectra Manager] refers to the entire suite of measurement/analysis and administrative programs for JASCO CD spectrometers.

Specifically, [Spectra Manager] is a software program that communicates with the CD Spectrometer to start various measurement and analysis programs. Common programs that are used include programs for performing measurements using the CD Spectrometer, the [Validation] program for checking the performance of the CD spectrometer, and the [Spectra Analysis], [JASCO Canvas] and [Administrative Tools] programs.

Spectra Measurement	Controls the CD spectrometer and measures samples.
Validation	Checks the performance of the CD spectrometer.
Spectra Analysis	Displays, processes, and prints the data obtained by [Spectra Measurement].
JASCO Canvas	Customizes the format of spectral data for printing.
Administrative Tools	Controls the administration of the CD spectrometer and measurement/analysis programs including the assignment of user authority.

This manual mainly explains how to use several of the measurement programs. Refer to [Hardware Manual] for details about how to use the CD spectrometer. Refer to the individual instruction manuals for details about the [Validation], [Interval Scan Measurement], [Spectra Analysis], [Interval Data Analysis], and [JASCO Canvas] programs.

2. Starting and Exiting Programs and the [Spectra Manager] Program Reference

2.1 Starting the program

2.1.1 Turning on the CD Spectrometer

- (1) Confirm that the standard detector is installed in the J-1000 series CD spectrometer. If another detector has been installed, replace it with the standard detector. Refer to the hardware manual "Model J-1100/1500 CD Spectrometer" or "Model J-1700 CD Spectrometer" for instructions on removing or installing the detector.
- (2) Supply nitrogen gas at a flow rate of at least 2 L/min. Absorption due to oxygen increases with decreasing wavelength, which reduces the light intensity reaching the detector. Please increase the flow rate to optimize the intensity within the desired wavelength range.

Nitrogen flow rate for different measurement wavelength ranges:

Up to 185 nm	2 L/min
Up to 180 nm	5 L/min
Up to 175 nm	15 L/min
Less than 175 nm	More than 15 L/min

Note: Depending on the residual amount of nitrogen gas present when the spectrometer is first turned on, it is possible for absorption due to oxygen to be high even if the recommended nitrogen flow rate is used. In this case, pre-purge at a nitrogen flow rate of 15 L/min before performing measurements. The recommended pre-purge times are: 10 min for wavelengths up to 185 nm, 30 min for wavelengths up to 180 nm and 60 min for wavelengths up to 175 nm.

- (3) When using the 450 W Xe lamp, feed cooling water to the light source at a flow rate of about 2 L/min.
- (4) Turn on the "Power" switch at the left side of the CD spectrometer.

WARNING!: Running the xenon (Xe) lamp in air will generate ozone gas, which is harmful to humans. Always supply nitrogen gas for a few minutes before turning on the light source.

Note 1: When performing measurements at wavelengths shorter than 180 nm, to maintain the effect of the nitrogen purge, after inserting or removing samples close the sample compartment lid as soon as possible.

Note 2: A high nitrogen flow rate is necessary when performing measurements at wavelengths shorter than 175 nm because of the sharp increase in absorption due to oxygen.

Note 3: Ozone gas absorbs light at wavelengths of 200 to 300 nm and 600 to 700 nm. This makes it difficult to perform measurements in these wavelength regions and can also damage optical parts due to corrosion. Always supply nitrogen gas before turning on the light source.

Note 4: Never use helium gas. It will permeate into the photomultiplier tube, reducing the vacuum level and the functionality of the tube.

Note 5: If a PC-controlled nitrogen gas flow meter (an optional accessory) is installed, all that is necessary to do is to simply open the main valve of the gas cylinder. Nitrogen will start flowing when the spectrometer is started. The flow rate can be controlled via the measurement program.

Note 6: Step (3) is not necessary with the 150 W Xe lamp, which is cooled with air.

Note 7: Just turning ON the "Power" switch does not result in the light source being turned on. The light source will be turned on automatically when the measurement program is started.

2.1.2 Starting the PC and Windows and Registering a J-1000 Series CD spectrometer

Turn on the PC and monitor to start *Windows*. When a J-1000 series CD spectrometer is connected for the first time, registration of the CD spectrometer is required in [Spectra Manager].

Note: The operations described in this section can be performed only by Administrators.

- (1) Start [Administrative Tools], right-click [Instrument] and select [Register Instrument...] from the shortcut menu to display the window in Figure 2.1.

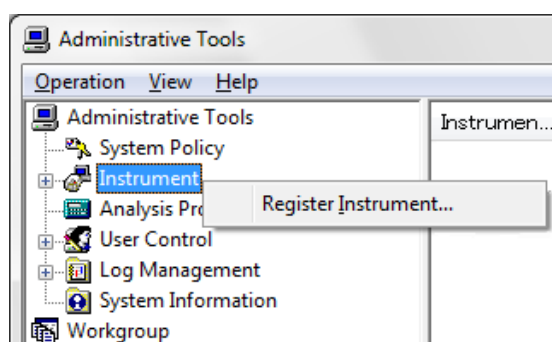


Figure 2.1 [Administrative Tools] window

- (2) Select the check box for the JASCO J-1000 series control driver and click the <Next> button to display the dialog box in Figure 2.2.


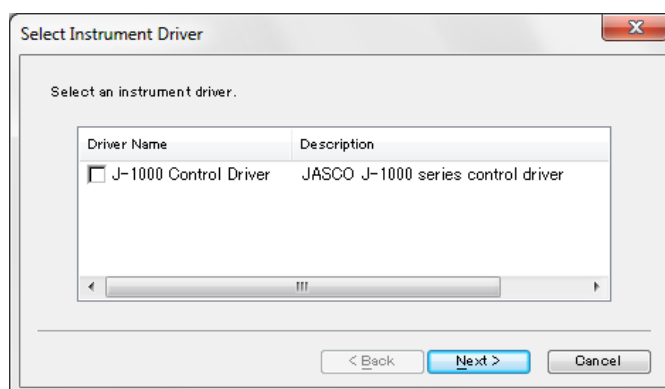


Figure 2.2 [Select Instrument Driver] dialog box

- (3) Enter the CD spectrometer name, model name and serial number and then click the <Finish> button to register the CD spectrometer. The model name and serial number are inscribed on the name plate on the rear panel of the CD spectrometer.

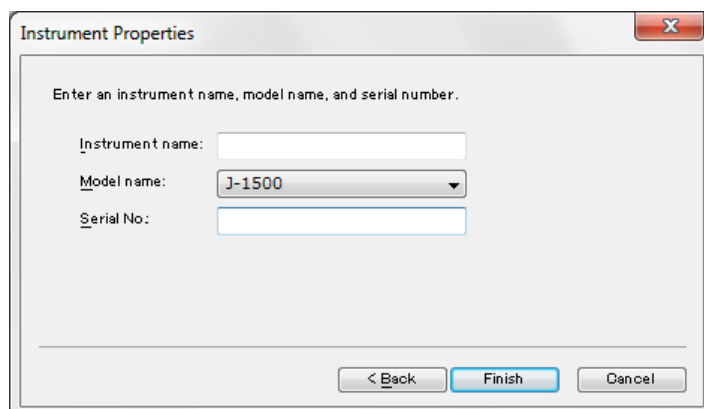

The 'Instrument Properties' dialog box contains the following fields and controls:

- Instruction: "Enter an instrument name, model name, and serial number."
- Instrument name:
- Model name:
- Serial No.:
- Buttons: "< Back", "Finish", "Cancel"

Figure 2.3 [Instrument Properties] dialog box

The CD spectrometer is registered and the name is displayed in the right pane of the [Administrative Tools] window.

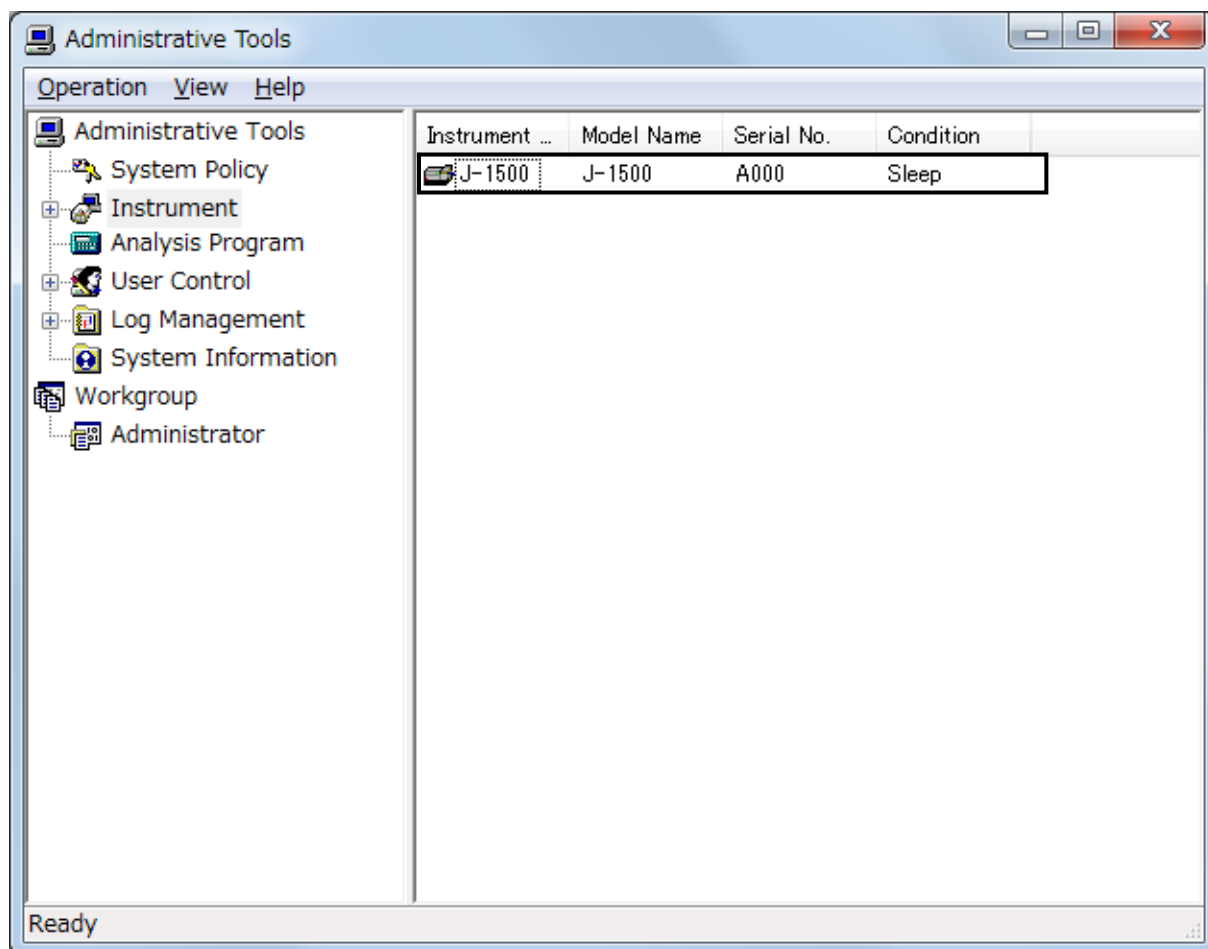


Figure 2.4 [Administrative Tools] window

Note: After registering the CD spectrometer, exit the program and perform the following operations after logging on again.

- (4) Restart [Spectra Manager] and start [Administrative Tools]. Right-click the registered CD spectrometer name and select [Properties...] to display the dialog box in Figure 2.5 . First, check that [Model name], [Display name], [Serial number], and [Driver version] are displayed in the [General] tab.

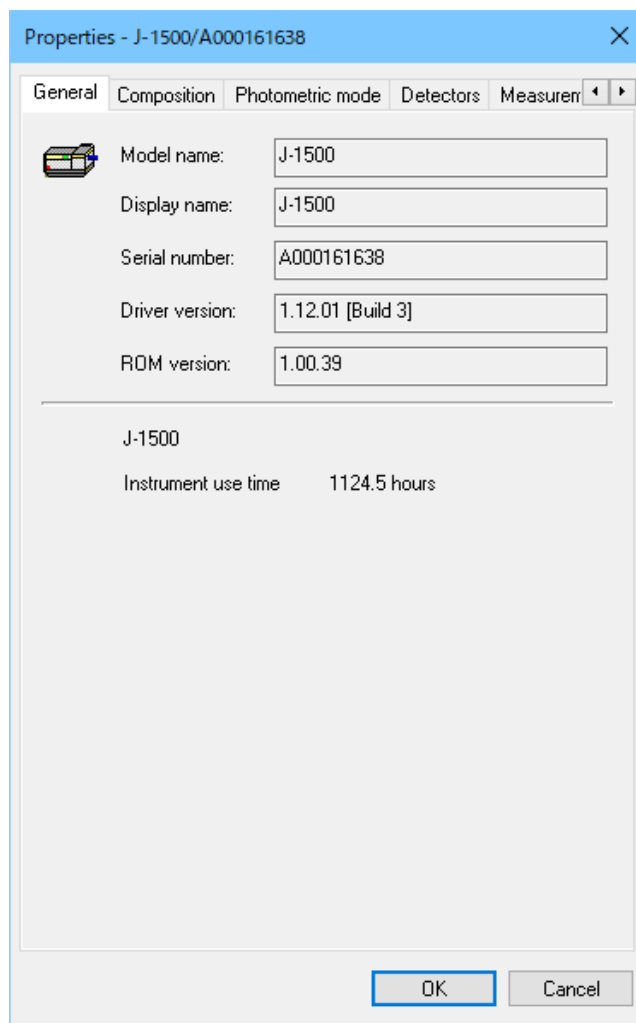


Figure 2.5 [General] tab in [Properties] dialog box

- (5) Click the [Composition] tab to set parameters for the nitrogen gas flow meter and the light source.

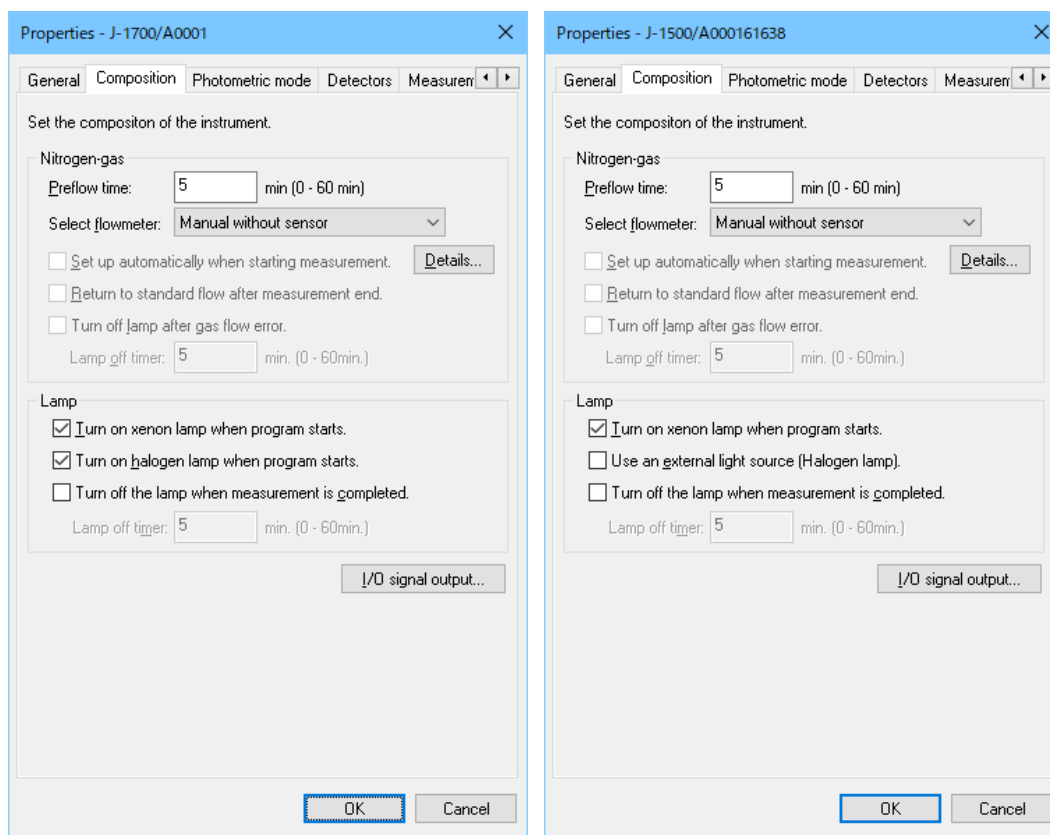


Figure 2.6 [Composition] tab in [Properties] dialog box J-1700 (left), J-1500 (right)

Nitrogen-gas	
[Prewflow time]	Specifies the nitrogen gas preflow time before the xenon light source is turned on.
[Select flowmeter]	<p>Selects the nitrogen gas flow meter.</p> <p>[Manual without sensor]: Select this when using a manual flow meter with no sensor.</p> <p>[Manual with sensor]: Select this when using a manual flow meter with a sensor.</p> <p>[Automatic (PC control)]: Select this when using a PC-controlled automatic flow meter.</p>
[Details]	<p>Specifies the flow rate of nitrogen gas under each condition.</p> <div style="border: 1px solid black; padding: 5px; margin-top: 10px;"> <p><i>Note: Refer to the "J-1000 Nitrogen Flow Meter Instruction Manual" for details.</i></p> </div>
[Set up automatically at measurement]	Specifies the nitrogen gas flow rate for the idling or measurement state. The flow rate is automatically controlled. This function is only available when the [Automatic (PC control)] flow meter is selected.

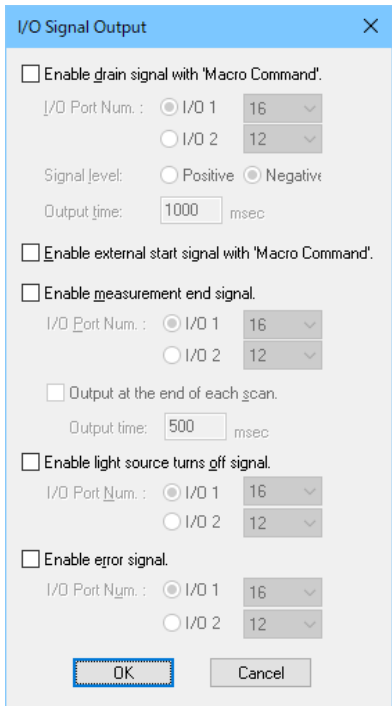
[Return to standard flow after a measurement end]	Specifies that the flow rate is automatically returned to the standard flow rate after measurement. This function is only available when the [Automatic (PC control)] flow meter is selected.
[Turn off lamp after gas flow error]	Specifies that the lamp is automatically turned off after a gas flow error has occurred. This function is only available when the [Automatic (PC control)] or [Manual with sensor] flow meter is selected.
[Lamp off timer]	Specifies the time before the light source is automatically turned off after a nitrogen gas flow error has occurred. This function is only available when the [Automatic (PC control)] or [Manual with sensor] flow meter is selected.
Lamp	
[Turn on xenon lamp when program starts]	Specifies that the Xe lamp is automatically turned on when the measurement program is started.
[Turn on halogen lamp when program starts]	Specifies that the Wl lamp is automatically turned on when the measurement program is started. (only for J-1700)
[Use an external light source (Halogen lamp)]	Specifies that the external halogen lamp is used. (only for J-1500)
<I/O signal output...>	<p>Specifies that each signal (drain, external start, measurement end, light source turns off and error) from the CD spectrometer is sent to the accessory attached to the I/O ports. Select the desired check boxes. When the accessory is made by JASCO, these settings are unnecessary.</p> 

Figure 2.7 [I/O Signal Output] dialog box

- (6) Click the [Photometric mode] tab to register the photometric mode.

The electric signals except other than CD signals can be detected using the J-1000 series CD spectrometer. The photometric modes selected in this dialog box are displayed in the measurement program. For J-1100, the available photometric modes are limited.

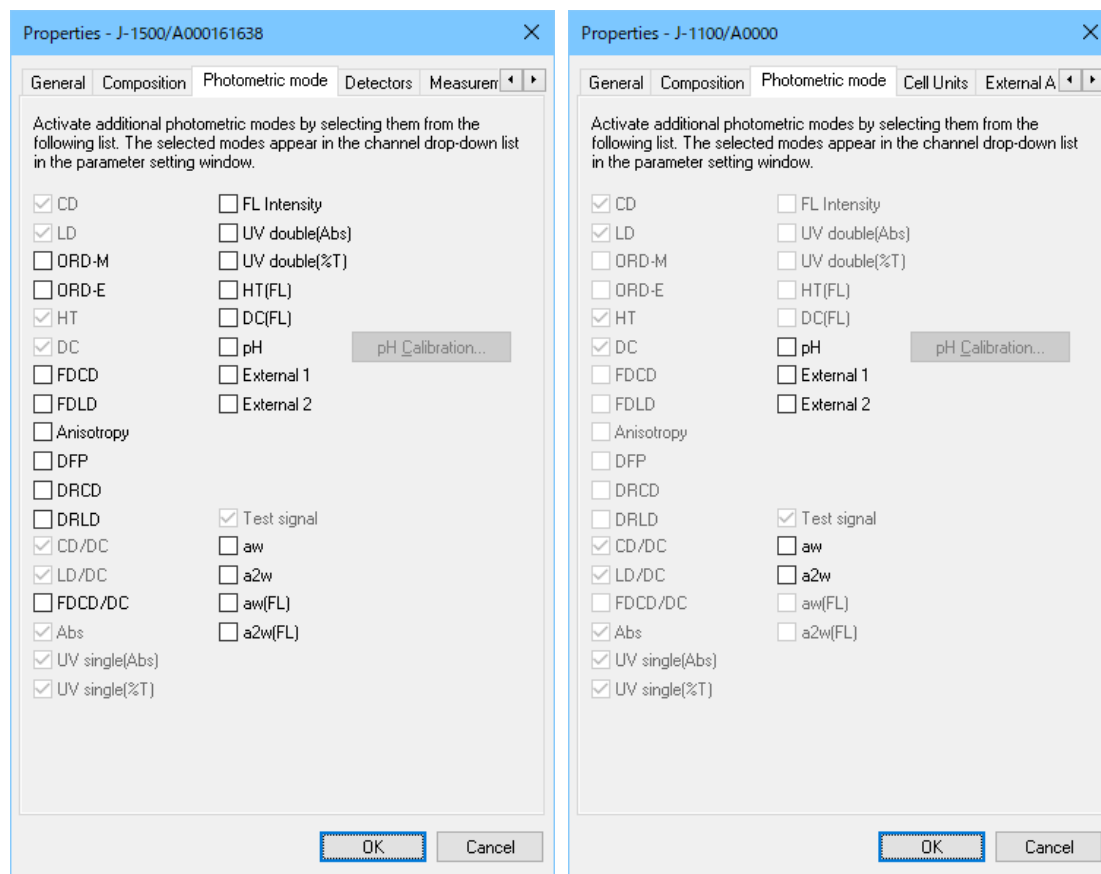


Figure 2.8 [Photometric mode] tab in [Properties] dialog box J-1500/1700 (left), J-1100 (right)

CD	Measures the circular dichroism signal
LD	Measures the linear dichroism signal
ORD-M	Measures the optical rotational dispersion signal. The ORD-M accessory is required. (Not available for J-1100)
ORD-E	Measures the optical rotational dispersion signal. The ORD-E accessory is required. (Not available for J-1100)
HT	Measures the high-tension (HT) voltage applied to the CD photomultiplier tube in the transmission configuration.
DC	Measures the direct current (DC) voltage applied to the CD photomultiplier tube in the transmission configuration.

FDCD	<p>Measures the fluorescence detected circular dichroism signal. The FDCD accessory is required. (Not available for J-1100)</p> $\Delta F = F_l - F_r$ <p>F_l: Fluorescence intensity excited by left circularly polarized light</p> <p>F_r: Fluorescence intensity excited by right circularly polarized light</p>
FDLD	<p>Measures the fluorescence detected linear dichroism signal. The FDCD accessory is required. (Not available for J-1100)</p> $\Delta F = F_{//} - F_{\perp}$ <p>$F_{//}$: Fluorescence intensity excited by horizontally linear polarized light</p> <p>F_{\perp}: Fluorescence intensity excited by vertically linear polarized light</p>
Anisotropy	Measures the fluorescence anisotropy signal. (Not available for J-1100).
D.F.P.	Measures the degree of fluorescence polarization signal. (Not available for J-1100).
DRCD	Measures the circular dichroism signal using detector for fluorescence configuration
DRLD	Measures the linear dichroism signal using detector for fluorescence configuration
CD/DC	Measures the CD and DC signals and displays the CD signal divided by the DC signal.
FDCD/DC	Measures the FDCD and DC signals and displays the FDCD signal divided by the DC signal. (Not available for J-1100).
Abs.	Measures the absorption spectrum. The absorption signal is obtained by conversion of the HT signal. Select this when the HT voltage for the photomultiplier tube in the transmission configuration is set to Auto.
UV single (Abs.)	Measures the absorption spectrum. The absorption signal is obtained by conversion of the DC signal. Select this when the HT voltage for the photomultiplier tube in the transmission configuration is set to Manual.
UV single (%T)	Measures the transmittance (%T) spectrum. The %T signal is obtained by conversion of the DC signal. Select this when the HT voltage for the photomultiplier tube in the transmission configuration is set to Manual.

Intensity	Measures the fluorescence intensity. Select this when the HT voltage for the photomultiplier tube in the fluorescence configuration (90 degrees with respect to the measurement beam direction) is set to Auto. (Not available for J-1100)
UV double (Abs.)	Measures the absorption spectrum using a UV double beam attachment. (Not available for J-1100)
UV double (%T)	Measures the transmittance spectrum using a UV double beam attachment. (Not available for J-1100)
HT(FL)	Measures the applied voltage (HTV) to the photomultiplier tube in the fluorescence configuration (90 degrees with respect to the measurement beam direction). (Not available for J-1100)
DC(FL)	Measures the direct current (DC) voltage to the photomultiplier tube in the fluorescence configuration (90 degrees with respect to the measurement beam direction). (Not available for J-1100)
pH	Measures the direct current signal from the pH meter. The pH meter is required.
External 1	Measures the signal from external signal input 1 ("EXT 1" terminal on the rear panel of the amplifier).
External 2	Measures the signal from external signal input 2 ("EXT 2" terminal on the rear panel of the amplifier).
Test signal	This is mainly used for checking the instrument.
aw	Measures the 50 kHz signal from the detector in the transmission configuration for the measurement beam. (Used for JASCO service engineers.)
a2w	Measures the 100 kHz signal from the detector in the transmission configuration for the measurement beam. (Used for JASCO service engineers.)
aw(FL)	Measures the 50 kHz signal from the detector in the fluorescence configuration (90 degrees with respect to the measurement beam direction) for the measurement beam. (Used for JASCO service engineers. Not available for J-1100)
a2w(FL)	Measures the 100 kHz signal from the detector in the fluorescence configuration (90 degrees with respect to the measurement beam direction) for the measurement beam. Used for JASCO service engineers. (Not available for J-1100)

The example of combination in accessories and photometric mode

Accessory	Potometric mode
FDCD unit	FDCD, FDL D, DC(FL) Set the FL detector HT voltage mode to Manual
Integrating sphere	DRCD, DRLD, HT(FL)
Emission monochromator (FDT-538)	Intensity Set the FL detector HT voltage mode to Manual
Stopped flow	CD/DC, DC Set the CD detector HT voltage mode to Manual
ORDM-520	ORD-M, HT
ORDE-521	ORD-E, HT

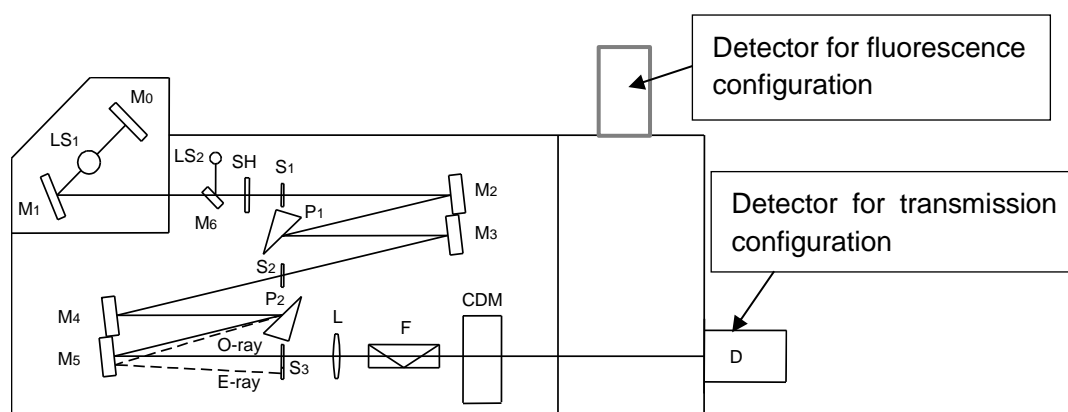


Figure 2.9 Configuration of detectors

- (7) Click the [Detectors], [Measurement Units], [Cell Units] and [External Accessories] tabs to register detectors, measurement units, cell units, and external accessories, respectively. Refer to Chapter 3 for details about registering accessories.

(8) Click the [Adjustments] tab to perform CD spectrometer adjustments.

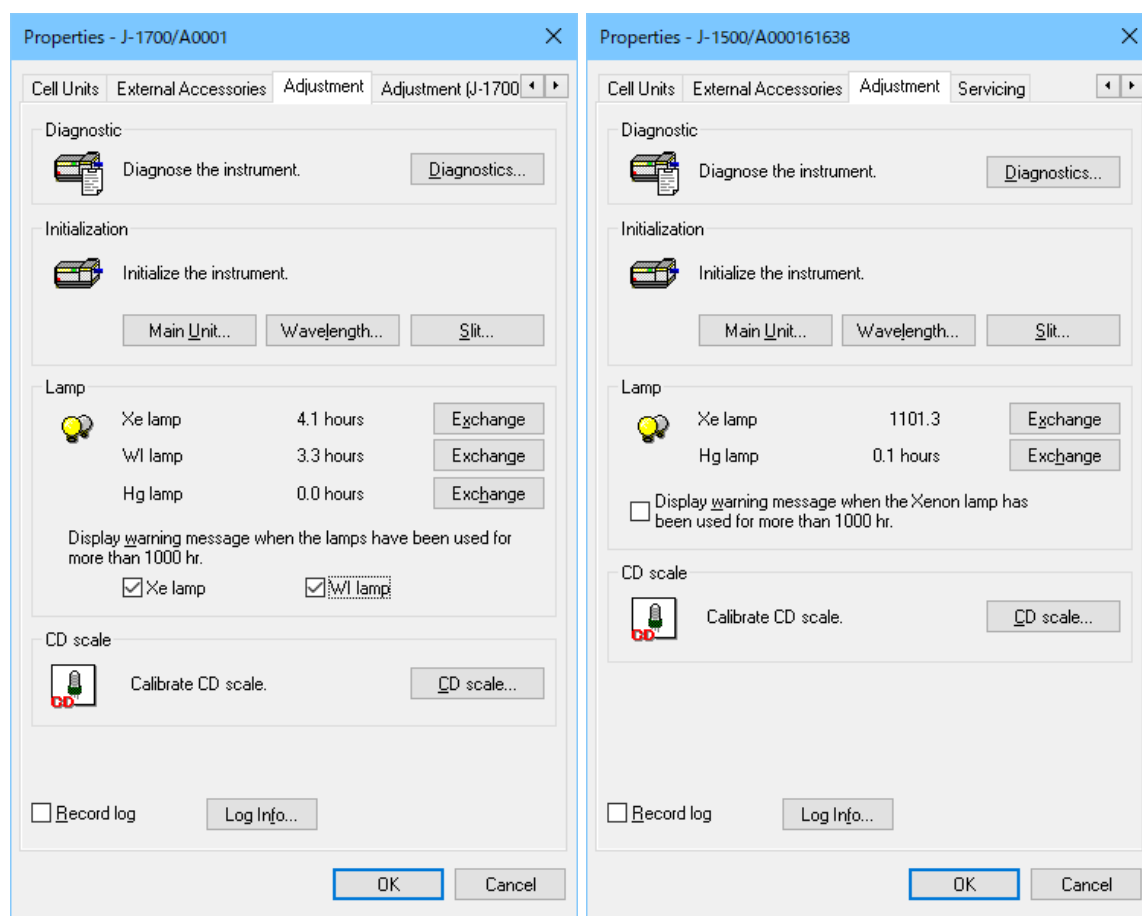
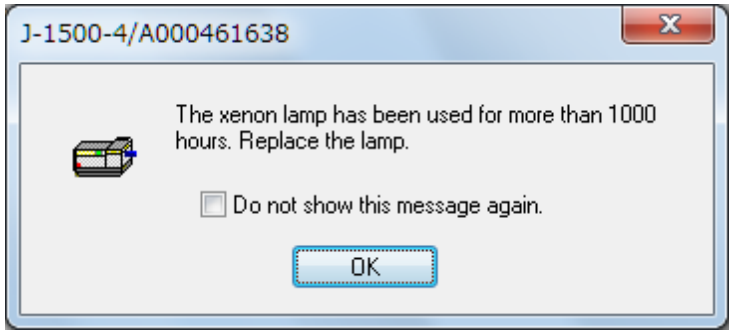
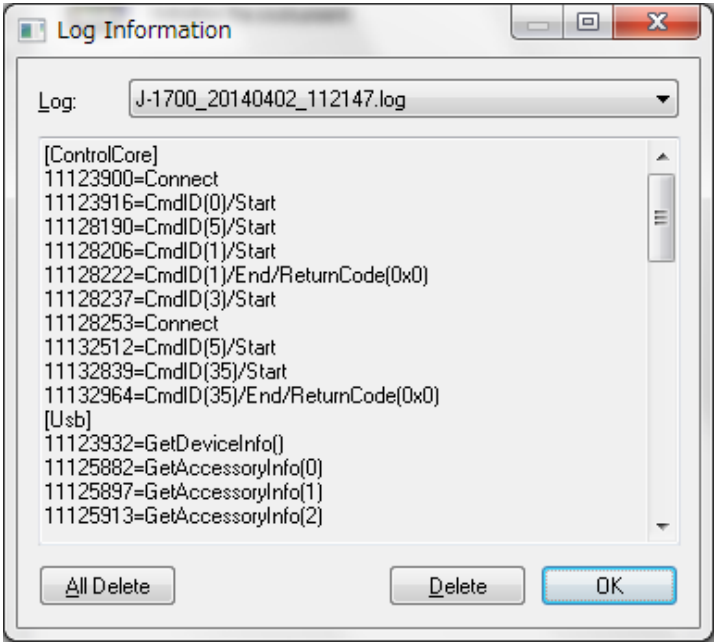


Figure 2.10 [Adjustments] tab in [Properties] dialog box 1700 (left), J-1100/1500 (right)

Diagnostic	
<Diagnostics...>	Confirms whether the CD spectrometer is functioning correctly.
Initialization	
<Main Unit>	Initializes the instrument settings.
<Wavelength>	Initializes the wavelength settings.
<Slit>	Initializes the slit width settings.
Lamp	
[Xe lamp]	Displays the number of hours the xenon lamp has been used. When the xenon lamp is exchanged, click the <Exchange> button to reset the number of hours used to 0.
[Wl lamp]	Displays the number of hours the halogen lamp has been used. When the halogen lamp is exchanged, click the <Exchange> button to reset the number of hours used to 0. (only for J-1700)

[Hg lamp]	Displays the number of hours the mercury lamp has been used. When the mercury lamp is exchanged, click the <Exchange> button to reset the number of hours used to 0.
[Display warning message when the lamps have been used for more than 1000 hr]	<p>Displays a warning message when the xenon and/or halogen lamp have been used for more than 1000 hours (see Figure 2.11). Halogen lamp can be selected only for J-1700.</p>  <p>Figure 2.11 Warning message for xenon lamp usage time</p> <div data-bbox="651 943 1417 1128" style="border: 1px solid black; padding: 5px;"> <p><i>Note: The lifetime of the xenon lamp is approximately 1000 hours. When it has been used for more than 1000 hours, the decrease in intensity of the light source will effect measurements.</i></p> </div>
CD scale	Calibrates the scale for CD signal. Used for users.
Record log	<p>Records a log of J-1000. When error occurs, send a log to JASCO service engineers.</p>  <p>Figure 2.12 [Log Information] dialog box</p>

Click the <Diagnostics...> button to display the [Diagnostics] dialog box. This function confirms whether the CD spectrometer is functioning correctly.

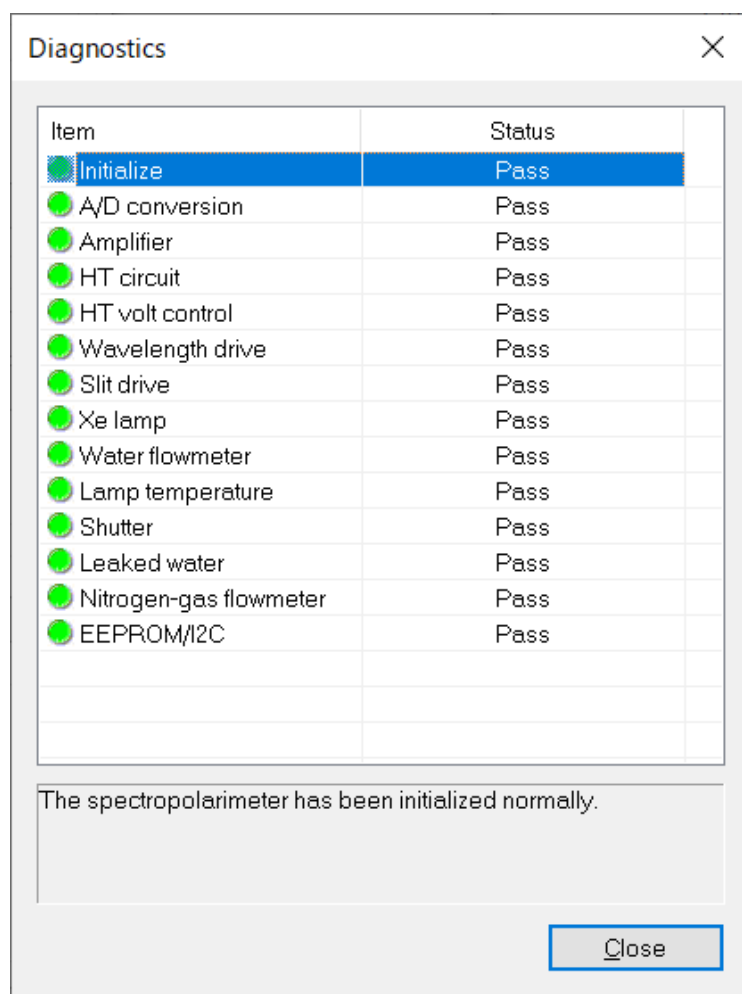


Figure 2.13 [Diagnostics] dialog box

If an error message is displayed in the message area, contact your local JASCO distributor.

- (9) Click the [Adjustment (J-1700)] tab to perform CD spectrometer adjustments. (Only J-1700)

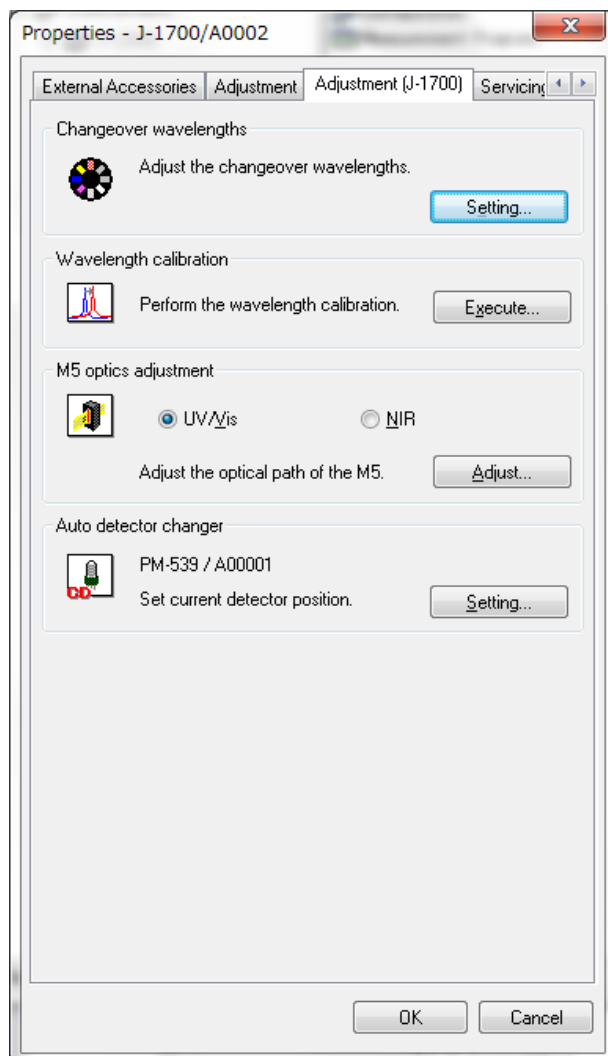


Figure 2.14 [Diagnostics] dialog box

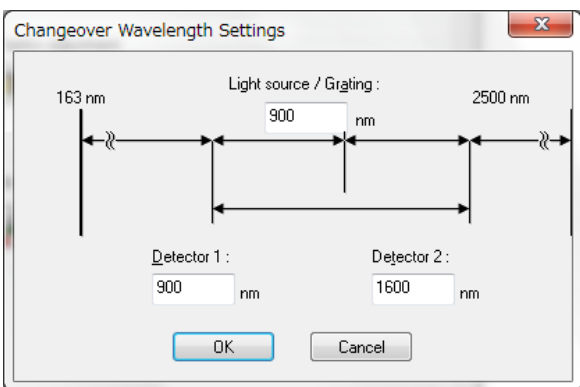
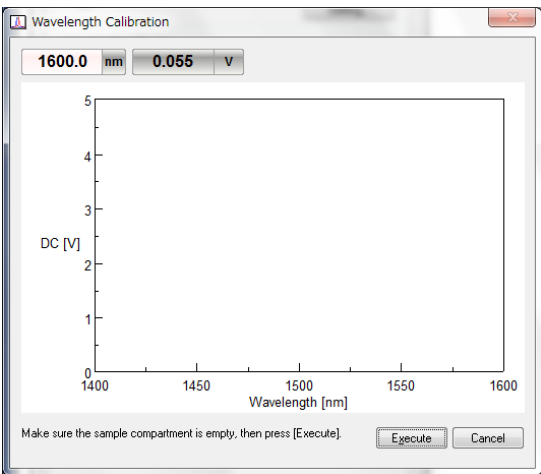
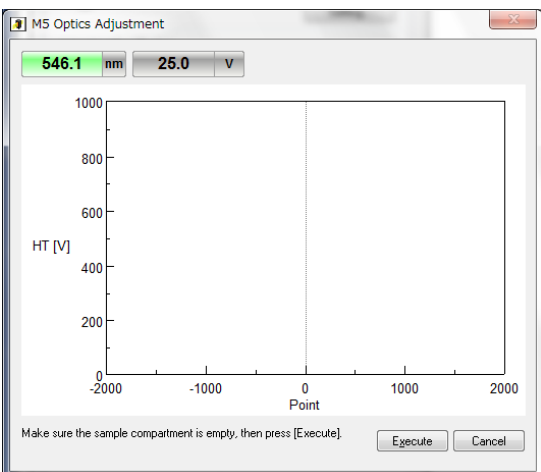
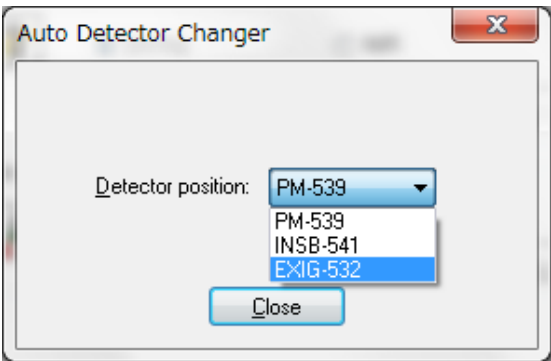
Changeover wavelengths	
<Setting...>	<p>Sets the changeover wavelengths for light source, grating and detectors.</p> 

Figure 2.15 [Changeover Wavelength Settings] dialog box

Wavelength calibration	
<Execute...>	<p>Calibrates the wavelength for grating in NIR region.</p>  <p>The dialog box titled 'Wavelength Calibration' features two input fields at the top: '1600.0 nm' and '0.055 V'. Below these is a graph with 'DC [V]' on the y-axis (0 to 5) and 'Wavelength [nm]' on the x-axis (1400 to 1600). At the bottom, there is a text instruction: 'Make sure the sample compartment is empty, then press [Execute].', followed by 'Execute' and 'Cancel' buttons.</p> <p>Figure 2.16 [Wavelength Calibration] dialog box</p>
M5 optics adjustment	
<Adjust>	<p>Adjusts the M5 mirror in UV/Vis or NIR region.</p>  <p>The dialog box titled 'M5 Optics Adjustment' has two input fields at the top: '546.1 nm' and '25.0 V'. Below is a graph with 'HT [V]' on the y-axis (0 to 1000) and 'Point' on the x-axis (-2000 to 2000). At the bottom, there is a text instruction: 'Make sure the sample compartment is empty, then press [Execute].', followed by 'Execute' and 'Cancel' buttons.</p> <p>Figure 2.17 [M5 Optics Adjustment] dialog box</p>
Auto detector changer	
<Setting>	<p>Sets current detector positions.</p>  <p>The dialog box titled 'Auto Detector Changer' contains a label 'Detector position:' next to a dropdown menu. The menu is open, showing three options: 'PM-539', 'INSB-541', and 'EXIG-532'. Below the menu is a 'Close' button.</p> <p>Figure 2.18 [Auto Detector Changer] dialog box</p>

- (10) Click the [Servicing] tab to calibrate CD spectrometer conditions such as the detector sensitivity. If calibration is required, contact your local JASCO distributor.

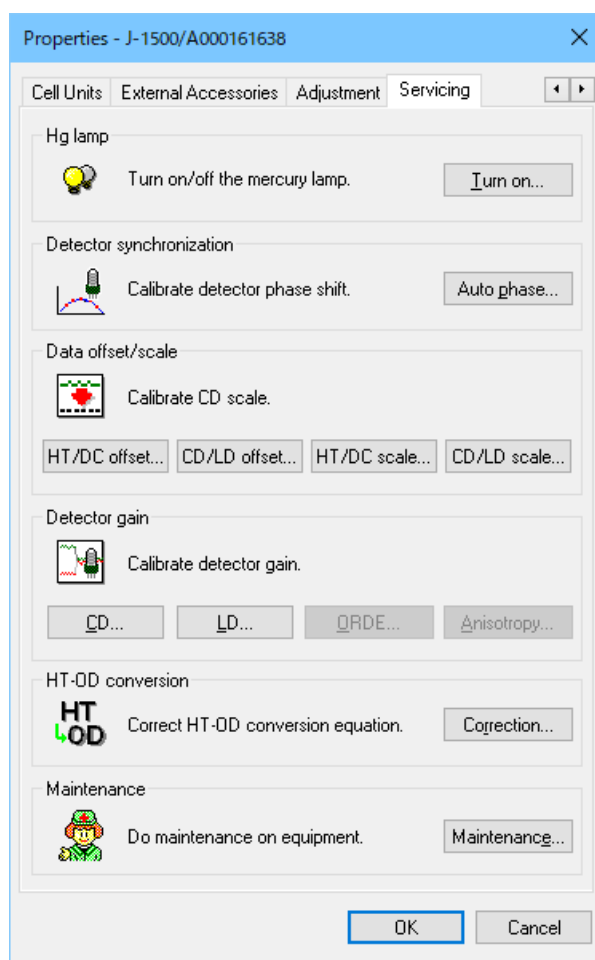


Figure 2.19 [Calibrations] tab in [Properties] dialog box

Hg lamp	Turns the mercury lamp on or off.
Detector synchronization	Calibrates the phase shift of the CD and FL detectors. Used for JASCO service engineers.
Data offset/scale	Calibrates the offset and scale of Main unit (J-1000). Used for JASCO service engineers.
Detector gain	Calibrates the scale for CD, LD, ORDE and anisotropy signals. Used for JASCO service engineers.
HT-OD conversion	Calibrates the HT-OD conversion equation. Used for JASCO service engineers.
Maintenance	Used for JASCO service engineers.

Note: In the CFR version, J-1000 must be registered to Workgroup before use. Refer to the [Spectra Manager CFR Administrative Tools] manual for details about registering the CD spectrometer to Workgroup.

2.1.3 Starting [Spectra Manager CFR]

- (1) After starting *Windows*, double-click the [Spectra Manager] icon located on the *Windows* desktop. Type the username and password in the logon dialog box to display the window in Figure 2.20.

Note: [Spectra Manager] can also be started by selecting [Start] - [All Programs] - [JASCO] - [Spectra Manager CFR] from the Windows [Start] menu.

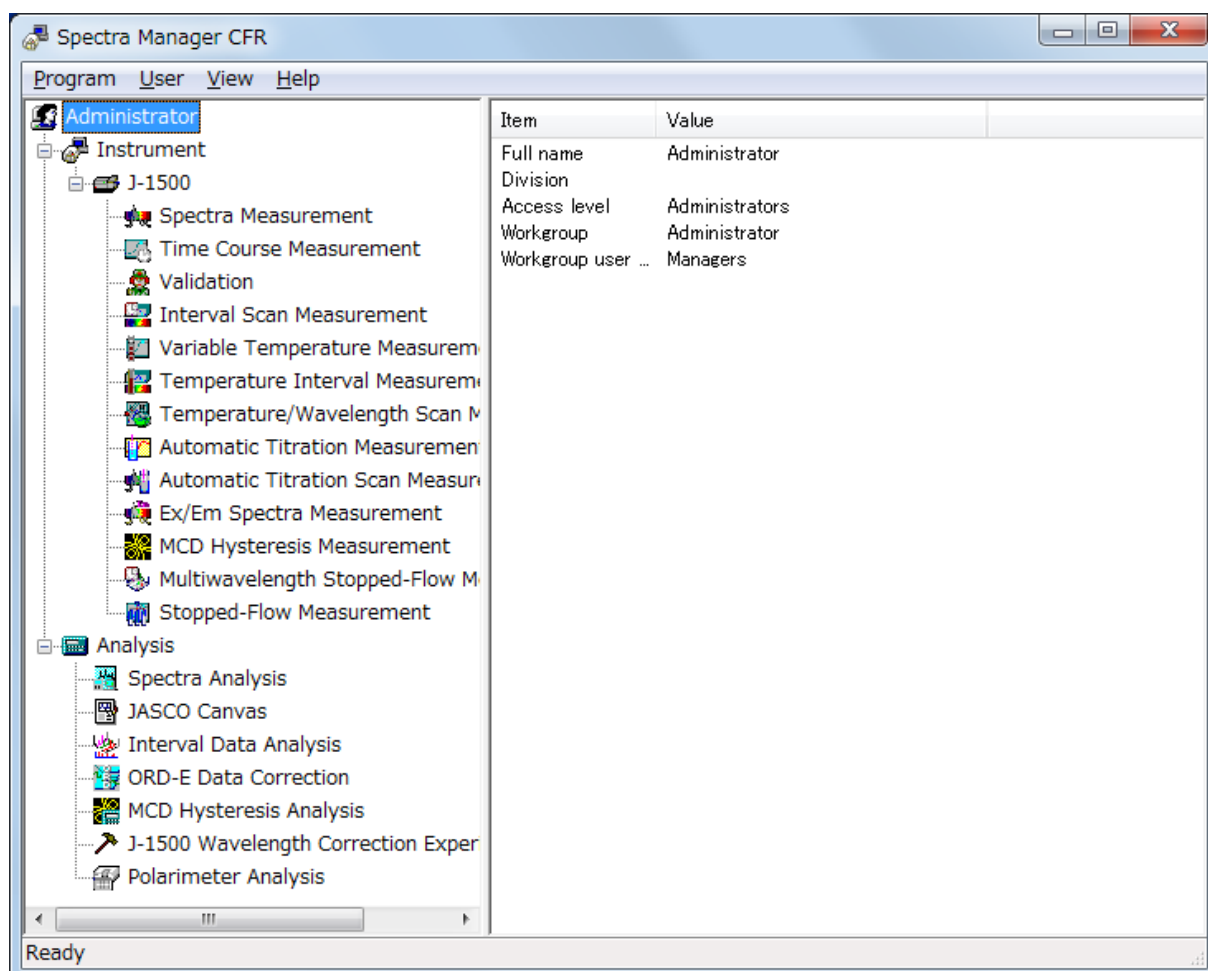


Figure 2.20 [Spectra Manager CFR] window

- (2) Start the [Spectra Measurement] program to measure a sample.

Double-click [Spectra Measurement] from the [Instrument] list of [Spectra Manager] to start the program. The window in Figure 2.21 is displayed.

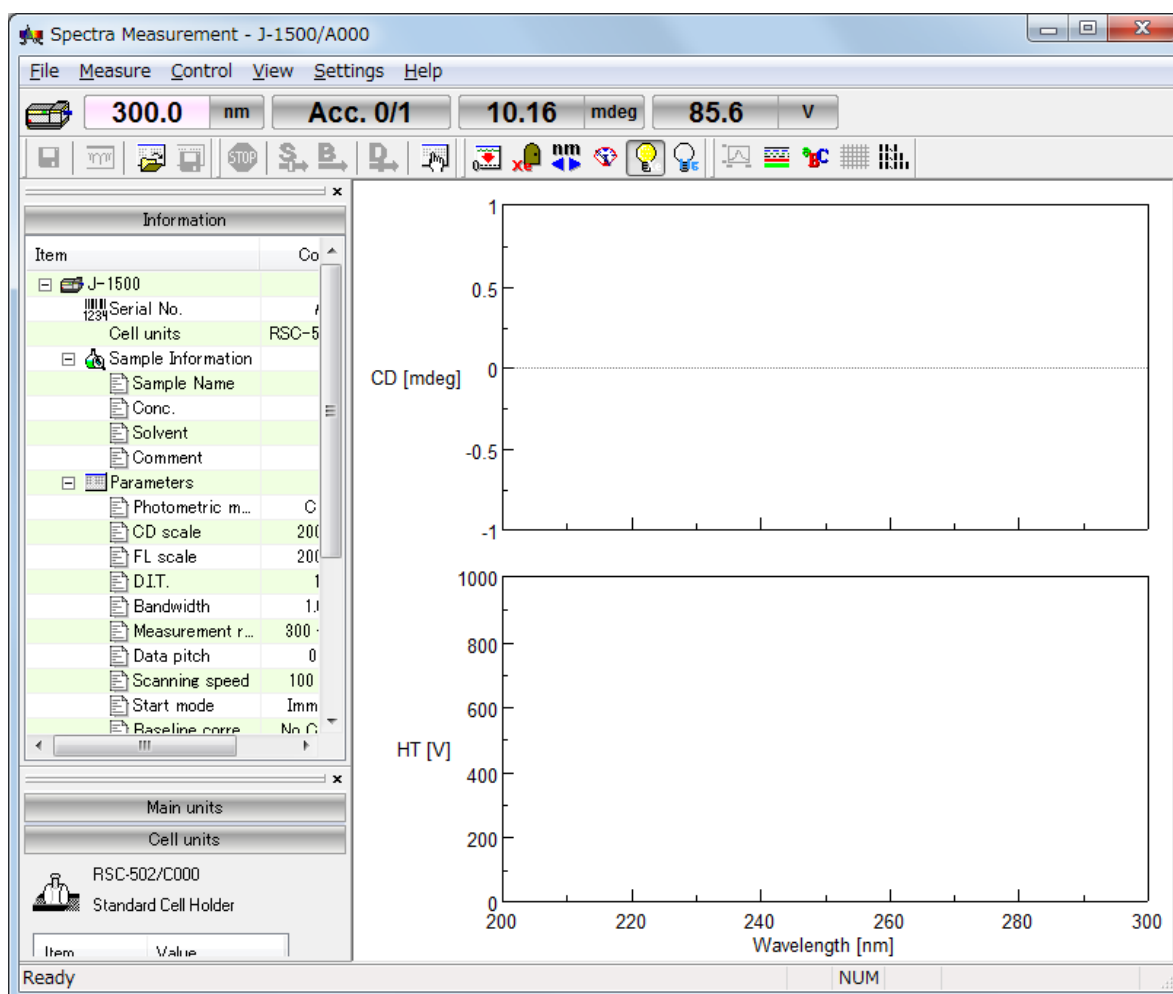


Figure 2.21 [Spectra Measurement] window

2.1.4 Components of the Program Window

In this section, the [Spectra Analysis] program window is used as an example to explain program operation components. The names of the various components (windows and dialog boxes) are also described.

View window

[Spectra Analysis] shows multiple spectrum display windows (called “Views”) in the main program window. Views have no dedicated menu bar, toolbar or status bar. Instead, the toolbar buttons and other components in the [Spectra Analysis] program window can be used for the active View window.

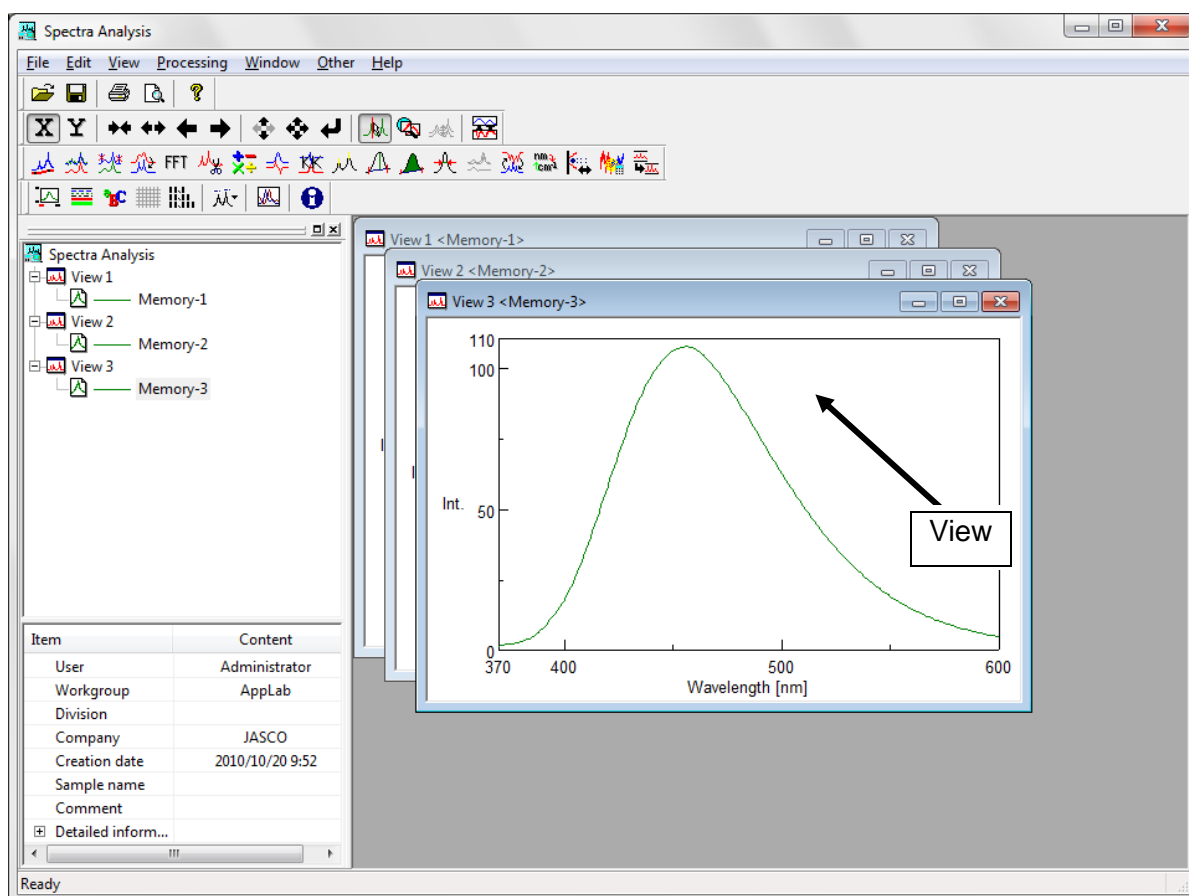


Figure 2.22 “Views” in [Spectra Measurement] program window

Dialog boxes

Commands and menu items ending with an ellipsis (...) open a dialog box when clicked. The opened dialog box usually contains parameters that must be specified. The [Styles] dialog box from the [View] menu is used as an example to explain the names and functions, as well as the operational rules for various sections of the dialog box.

To display the [Styles] dialog box in Figure 2.23, select [Styles...] from the [View] menu in the [Spectra Analysis] program. The names of various sections of the dialog box are indicated below.

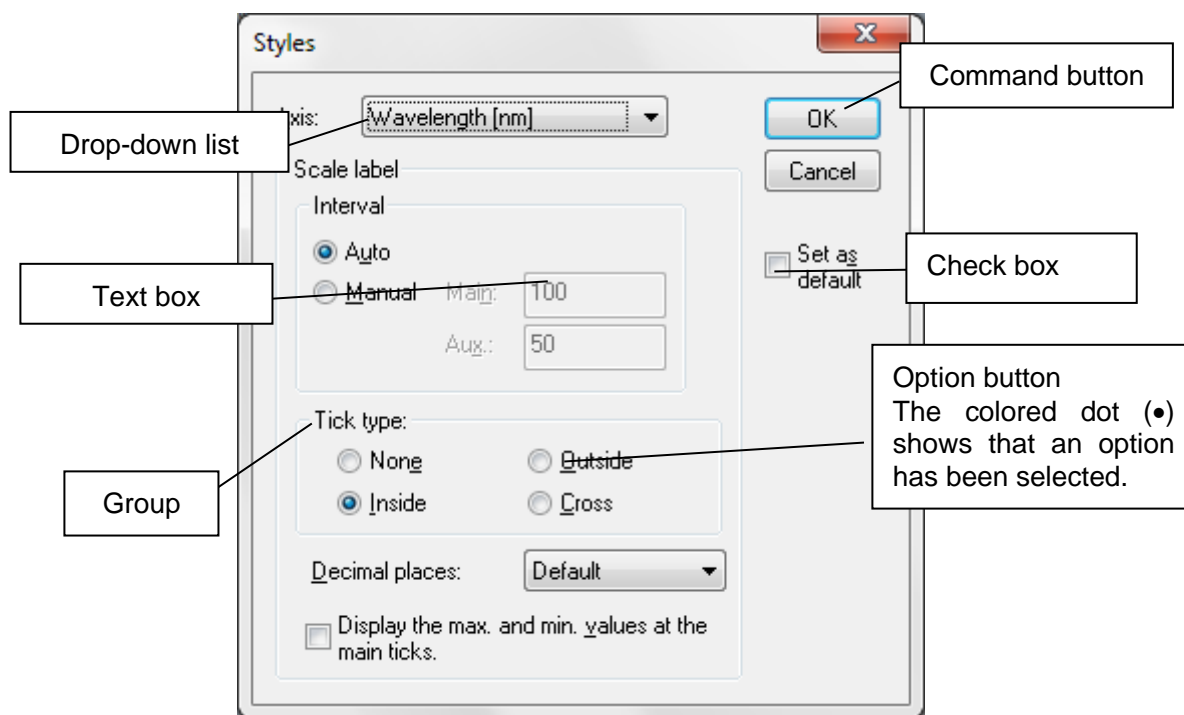


Figure 2.23 [Styles] dialog box

2.2 Exiting the Program

2.2.1 Exiting [Spectra Manager CFR]

- (1) Select [Exit] from the [Program] menu to close [Spectra Manager CFR].

Note: [Spectra Manager CFR] can also be exited by clicking the <Close> button.

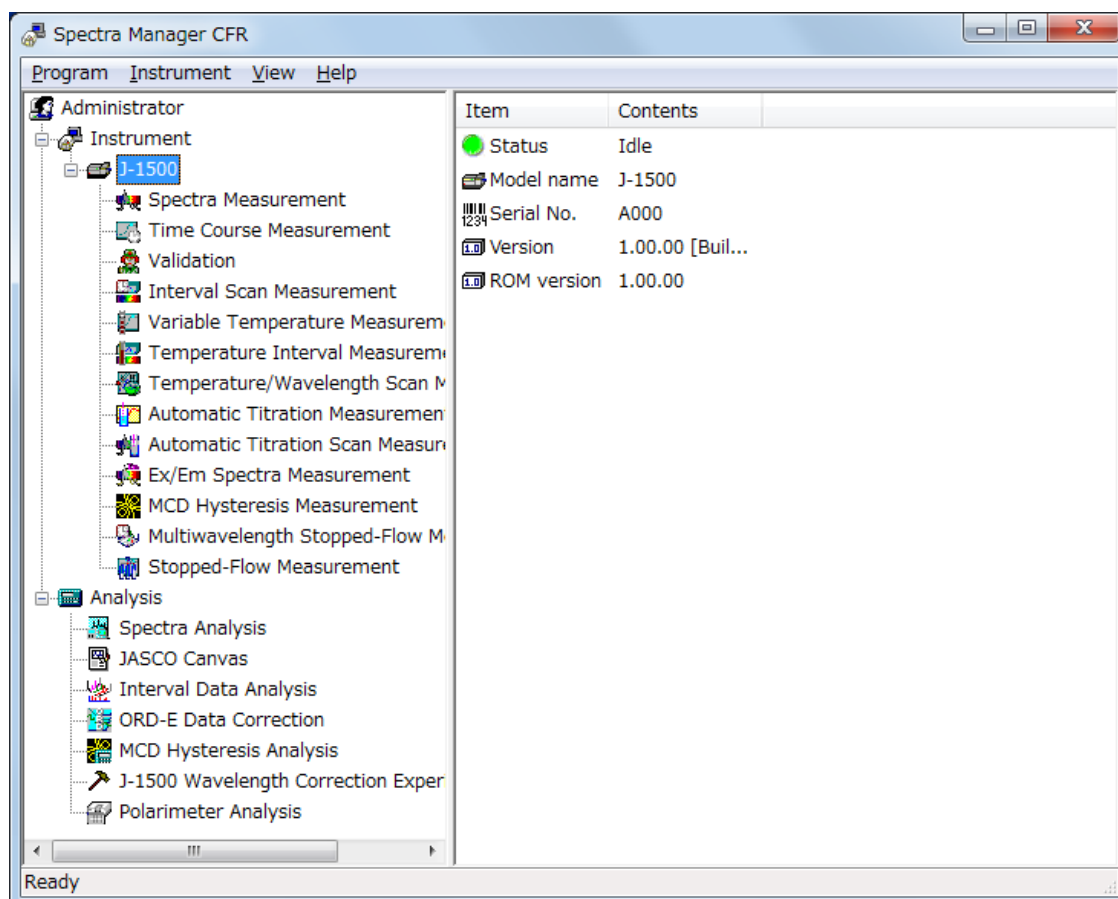


Figure 2.24 [Spectra Manager CFR] window

[Spectra Manager CFR] cannot be exited before closing the running measurement/analysis programs. If programs are running, the dialog box in Figure 2.25 is displayed. After clicking the <OK> button, close the running programs.

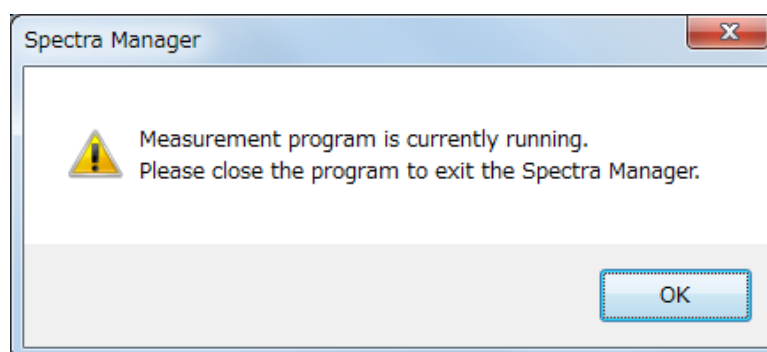


Figure 2.25 Message when a program is still running

(2) Exiting *Windows*

Exit *Windows* in accordance with standard *Windows* operation procedures.

2.2.2 Turning off the PC and CD Spectrometer

- (1) First, turn off the PC and monitor. Do not forget to turn off the monitor.
- (2) Ensure that the sample compartment is empty, and then turn off the CD spectrometer.

2.3 [Spectra Manager CFR] Program Reference

[Spectra Manager CFR] is an integrated software package that acts as a common platform for the entire range of JASCO analytical instruments.

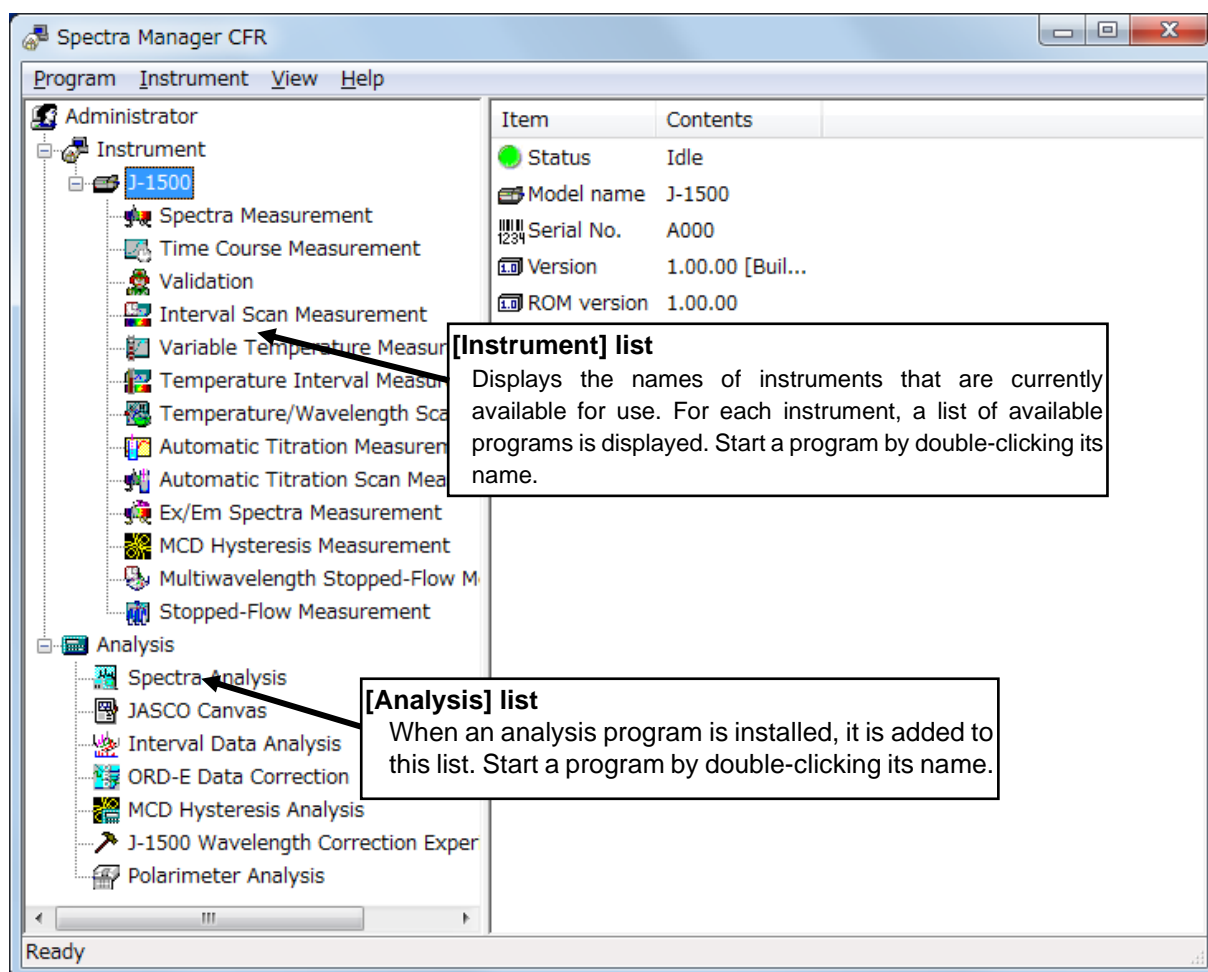


Figure 2.26 [Spectra Manager CFR] window

The standard measurement and analysis programs are displayed in Figure 2.26. When an optional program is installed, it is added to one of these lists.

Note 1: The CD Spectrometer is also referred to as the “instrument” in this manual.

Note 2: “Communication” refers to sending a control signal from the PC to the CD spectrometer or transferring measured data from the CD spectrometer to the PC.

Menu

[Program] menu	
[Administrative Tools]	Starts the [Administrative Tools] program.
[Change Password...]	Changes the current password.
[Logoff...]	Logs off the current user.
[Exit]	Exits [Spectra Manager].
[User] menu	
[Change Workgroup...]	Changes the Workgroup.
[View] menu	
[Status Bar]	Shows/hides the status bar.
[Help] menu	
[Contents]	Displays the Help contents window.
[Search Topic]	Displays the Help search function window.
[About...]	Displays version information about the CD spectrometer control driver.

3. Accessory Registration Reference

The J-1000 series CD spectrometers and software automatically recognize an attached accessory, display information about the accessory, and start registered programs based on recognized accessories. This chapter provides information of installing and registering accessories, the program that is started when an accessory is recognized, and the actions of [Spectra Manager CFR] when recognizing an accessory.

3.1 Installing/Attaching an Accessory

Accessories for J-1000 series CD spectrometers are automatically recognized. Information about the respective accessory is contained within the accessory itself. The accessory is recognized when its cable is connected to the sample compartment connector. When installing a J-1000 accessory, confirm that the accessory cable is properly connected to the CD spectrometer connector.

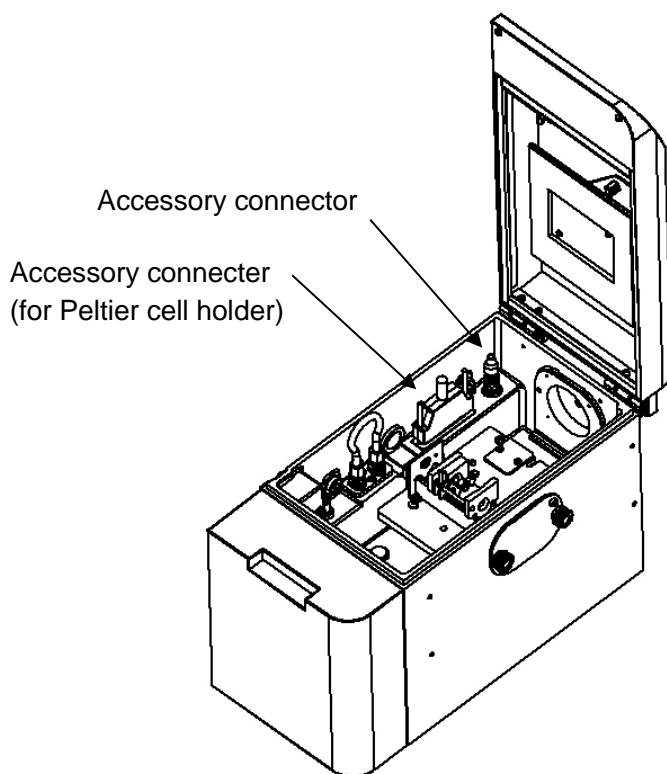


Figure 3.1 Location of J-1000 accessory connectors

3.2 Precautions for Peltier Thermostatic Cell Holder

The electrical power for the Peltier thermostatic cell holder is supplied from the J-1000 series CD spectrometer when this accessory is connected and detected by the instrument. When power is being supplied, an icon is displayed on the status bar.

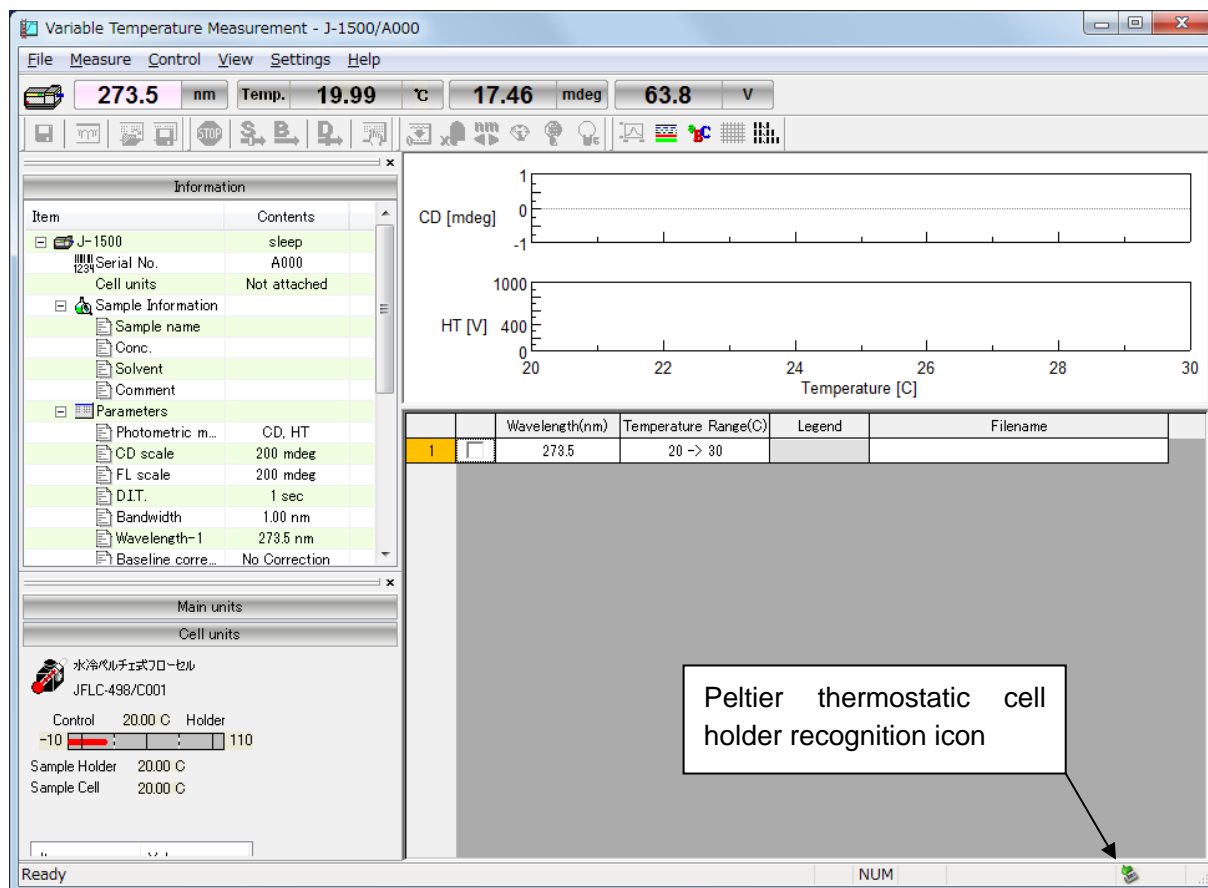


Figure 3.2 Icon displayed on the status bar after the Peltier thermostatic cell holder has been recognized

3.2.1 Removing the Peltier Thermostatic Cell Holder

This section describes how to remove the Peltier thermostatic cell holder while the J-1000 series instrument is running.

- (1) Click the cursor on the accessory icon to display the [Safely Remove Temp unit] pop-up menu.

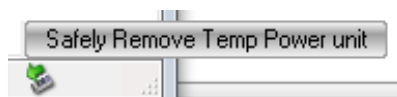


Figure 3.3 Pop-up prompt for safely removing the thermostatic unit

- (2) Left-click the command to display the [Safely Remove Thermostatic Unit] dialog box.

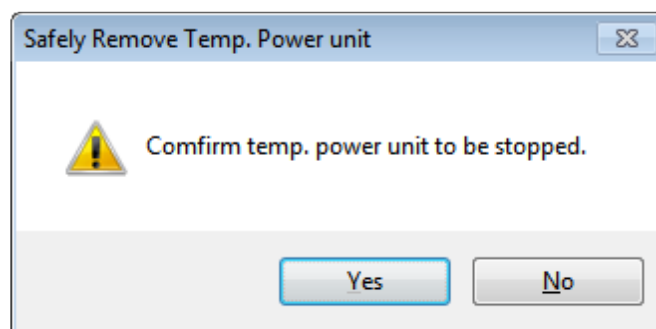


Figure 3.4 [Safely Remove Thermostatic Unit] dialog box

- (3) Click the <Yes> button. The electrical power supply for the Peltier thermostatic cell holder will be stopped and the cell holder can now be safely removed.

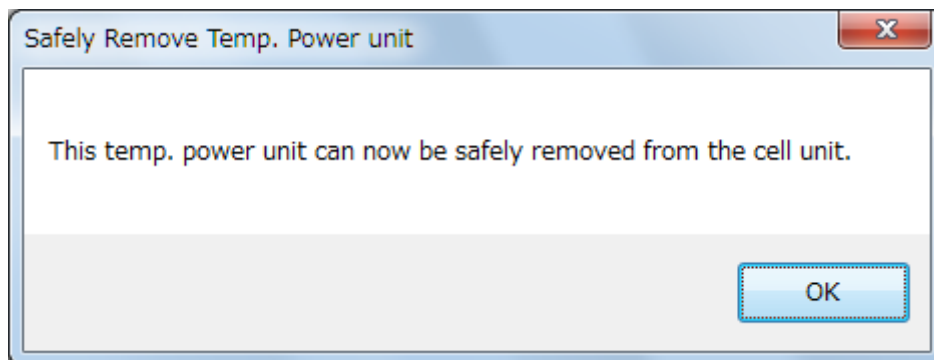


Figure 3.5 Confirmation message that thermostatic unit can be safely removed

To use the Peltier thermostatic cell holder again, reconnect its accessory cable to the J-1000 series instrument.

3.3 Registering Accessories

This section describes the procedure for registration of accessories used with the CD spectrometer.

3.3.1 Registering an Automatically-Recognized Accessory

- (1) Confirm that the instrument is turned on and that [Spectra Manager] is running.
- (2) Connect the automatically-recognized accessory to the sample compartment. The accessory name, accessory ID and serial number are automatically registered.

3.3.2 Registering an Accessory Requiring Manual Registration

Note: This operation can be performed only by Administrators.

- (1) Confirm that the CD spectrometer is turned on and that [Spectra Manager] is running.
- (2) After starting [Administrative Tools], right-click the instrument name and select [Properties...] (see Figure 3.6).

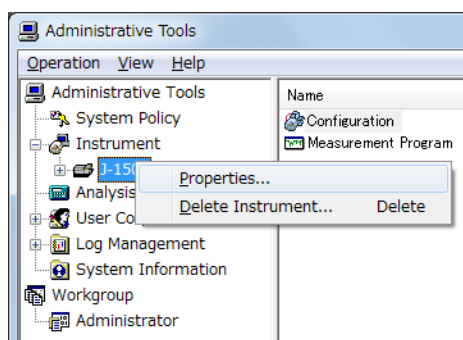


Figure 3.6 J-1000 instrument shortcut menu

- (3) Click the [Detectors] tab, [Measurement Units] tab, [Cell Units] tab or [External Accessories] tab.

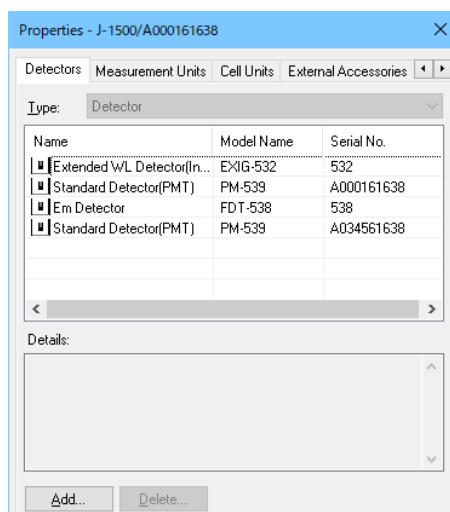


Figure 3.7 [Detectors] tab in [Properties] dialog box

- (4) Select the type from the [Type] drop-down list.

Table 3.1 Detectors

Type
Standard WL Detector (PMT): PM-539
Extended WL Detector (ExPMT): EXPM-531
Extended WL Detector (InGaAs): InGaAs-532
Extended WL Detector (APD) : EXAP-533
FDCD Detector (PMT): PMJ-534
FDCD Detector (ExPMT): EXPML-535
FDCD Detector (InGaAs): EXLGL-536
FDCD Detector (APD) : EXAPL-537
FL Intensity Detector : FDT-538
Em Detector: FDT-455
ORD-M Detector (PMT): PM-540
Extended WL Detector (InSb) : INSB-541 (only for J-1700)

Note: For J-1100, the [Detectors] tab is not displayed.

Table 3.2 Measurement Units

Type
ORD Attachment
Fluorescence Monochromator

Note: For J-1100, the [Measurement Units] tab is not displayed.

Table 3.3 Cell Units


Type
Cell Holder
Temperature Controller
Flow Cell
Integration Sphere
FDCD Attachment
Double Beam UV Attachment
MCD Controller
Stopped Flow Controller
Cryostat Holder
High Temp. Cell Unit

Note: For J-1100, only the cell holder and temperature controller can be selected.

Table 3.4 Accessories

Type
Automatic Titrator
Circulator
Autosampler
PAL Unit

- (5) Click the <Add...> button to display the dialog box in Figure 3.8.

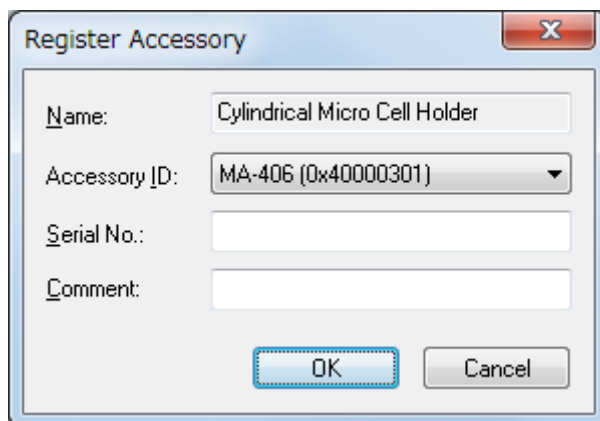

The image shows a Windows-style dialog box titled "Register Accessory". It has a standard title bar with a close button (X). The dialog contains four input fields: "Name:" with the text "Cylindrical Micro Cell Holder", "Accessory ID:" with a dropdown menu showing "MA-406 (0x40000301)", "Serial No.:" with an empty text box, and "Comment:" with an empty text box. At the bottom right, there are two buttons: "OK" and "Cancel".

Figure 3.8 [Register Accessory] dialog box

- (6) [Name] is automatically entered by selecting the accessory in the [Accessory ID] drop-down list. Type the serial number and a comment as required, and then click the <OK> button to register the accessory and display it in the list of registered accessories.

3.4 Actions of [Spectra Manager] When Recognizing an Accessory

This section describes the actions of [Spectra Manager] running a measurement program when an automatically-recognized accessory with registered programs is attached to or removed from the CD spectrometer.

Note: Do not attach/remove an automatically-recognized accessory while the program is displaying a dialog box or message.

3.4.1 When an Automatically-Recognized Accessory is Attached

When an automatically-recognized accessory is attached, a dialog box stating “An accessory has been attached” is displayed, and information on that accessory is displayed on the accessory bar.

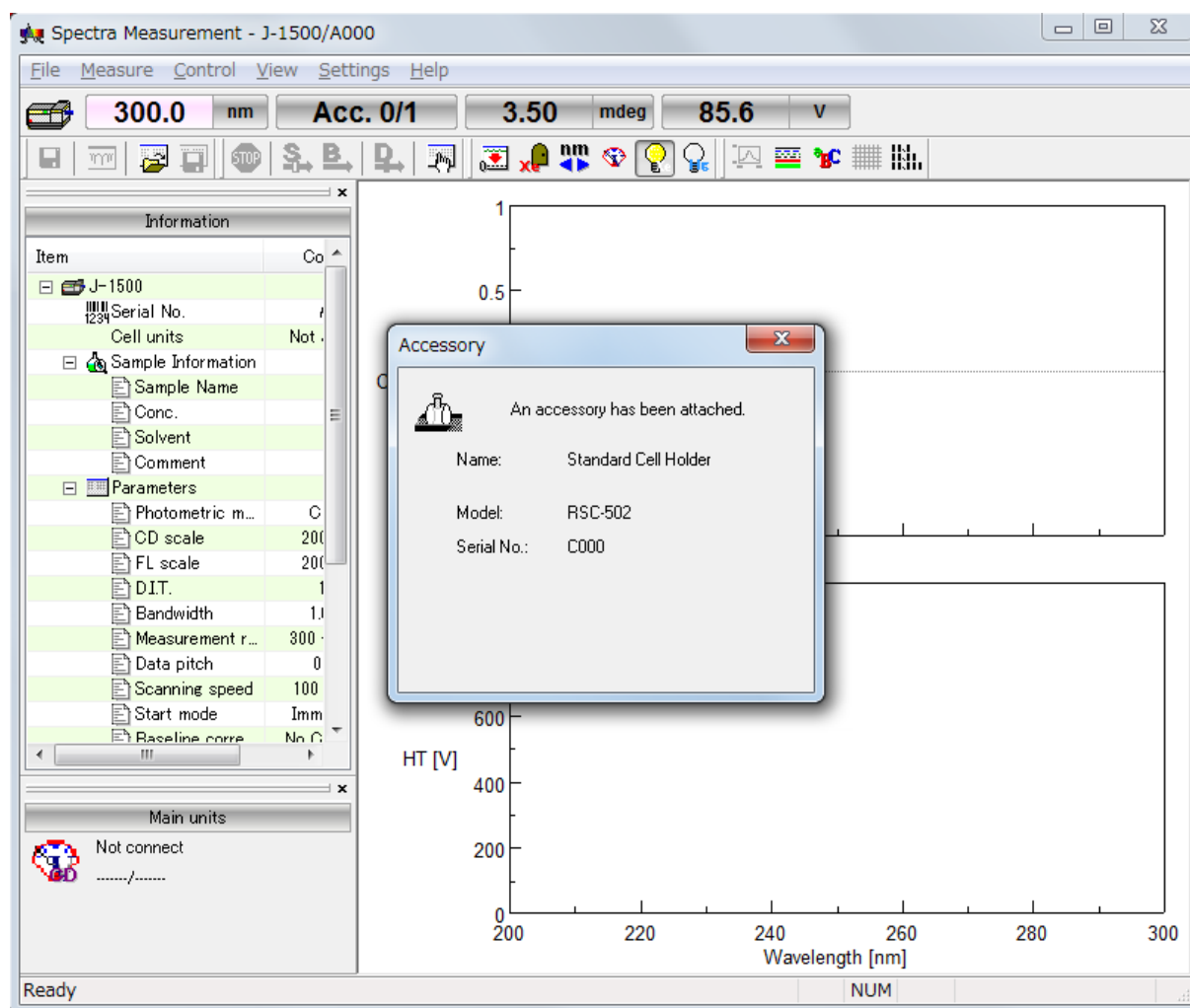


Figure 3.9 Message when an automatically-recognized accessory is being attached while [Spectra Manager] and a measurement program are running

3.4.2 When the Automatically-Recognized Accessory is Removed

When an automatically-recognized accessory is removed, a dialog box stating “An accessory was removed” is displayed, and information on that accessory disappears from the accessory bar.

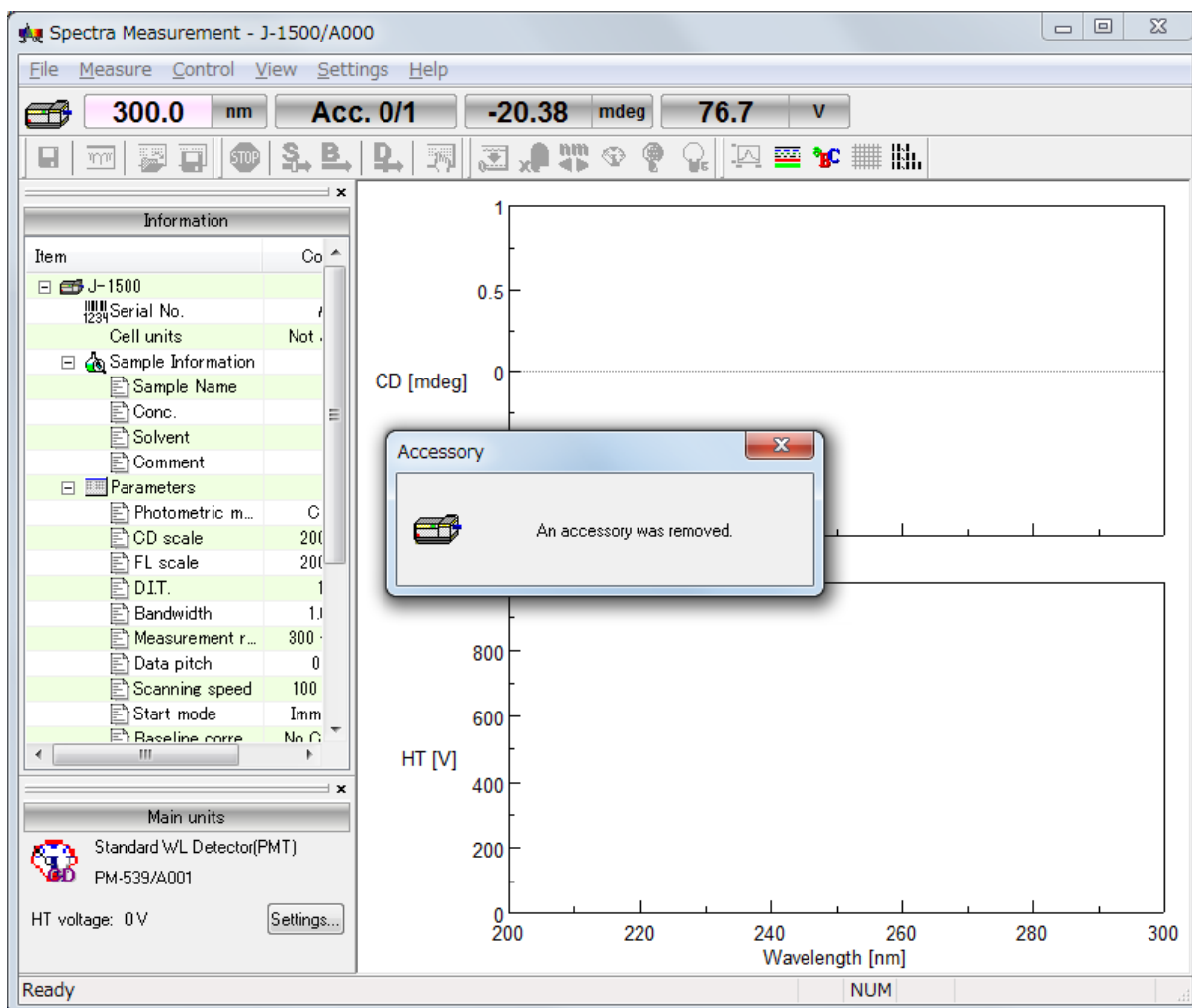


Figure 3.10 Message when an automatically-recognized accessory is removed while [Spectra Manager] and a measurement program are running

3.5 Status Display for the CD Spectrometer

This section describes the display of the CD spectrometer status in a measurement program.

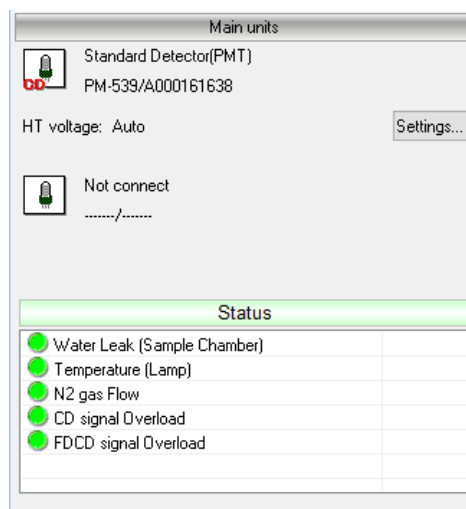


Figure3.11 Status of CD spectrometer

Water Leak (Sample Chamber)	When water leakage occurs in the sample compartment, the green indicator in the status box turns red. A stop valve, an optional accessory, automatically prevents water from leaking.
Temperature (Lamp)	When the light source temperature exceeds 80 °C, the green indicator turns red and the light source is turned off automatically.
N ₂ gas Flow	<p>When the nitrogen flow rate is less than the specified value, the green indicator turns red. This item is displayed only when the [Manually with sensor] or [Automatic (PC control)] flow meter is used.</p> <div style="border: 1px solid black; padding: 5px;"> <p><i>Note 1: For the [Manually with sensor] type, the sensor position can be changed and the flow rate can be specified.</i></p> <p><i>Note 2: For the [Automatic (PC control)] type, when the flow rate is less than the idling flow rate specified in the [Administrative Tools] program, the green indicator turns red.</i></p> </div>
CD signal Overload	When the CD signal exceeds the upper limit of the measurement range specified by [CD scale] in the [Parameters] dialog box, the green indicator turns red. Specify a larger CD scale.
FDCD signal Overload	When the FDCD signal exceeds the upper limit of the measurement range specified by [FL scale] in the [Parameters] dialog box, the green indicator turns red. Specify a larger FL scale.

4. Introduction to the [Spectra Measurement] Program

This chapter describes how to use the [Spectra Measurement] program, including procedures for starting the [Spectra Measurement] program, measuring the spectrum of a sample, and saving and printing data, for practice in operating a JASCO CD spectrometer.

Explanations of the parameters have been kept to a minimum, with a focus on the flow of operations. Follow the measurement procedures outlined below for practice in operating the CD spectrometer. Refer to Chapter 6 “Measurement Program Reference” for details about toolbar icons and menus, and to Chapter 7 “[Spectra Measurement] Program Menu Reference” for details about the [Parameters...] command.

4.1 Procedures for the [Spectra Measurement] Program

This section describes procedures ranging from measuring a CD spectrum of an ammonium d-10-camphorsulfonate solution to saving and printing the measurement data.

Starting the [Spectra Measurement] Program	Refer to Section 4.2
↓	
Specifying and Saving Measurement Parameters	Refer to Section 4.3
↓	
Reviewing and Approving the Parameter File*	Refer to Section 4.4
↓	
Measuring the Sample	Refer to Section 4.5
↓	
Saving Spectra*	Refer to Section 4.6
↓	
Printing Results	Refer to Section 4.7
↓	
Exiting the Program	Refer to Section 4.8

*The step in section 4.4 is only for the CFR version and 4.6 is only for the non CFR version

Note 1: In the CFR version, Managers or Analysts authority is required to specify measurement parameters.

Note 2: In the CFR version, measurement can basically be performed using only reviewed and approved measurement parameters. However, only the person who created parameter file can perform measurement without review and approval.

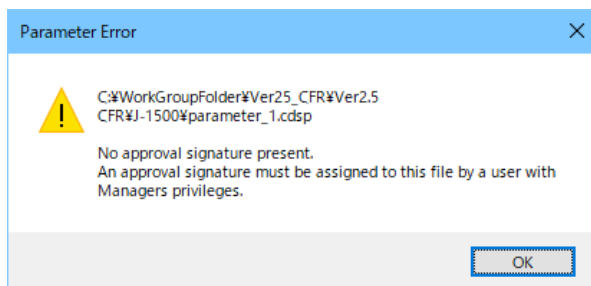


Figure 4.1 Message when measurement is attempted using unapproved and unsigned measurement parameters

4.2 Starting the [Spectra Measurement] Program

To start the program, double-click [Spectra Measurement] from the [Instrument] list of [Spectra Manager CFR]. The window in Figure 4.2 is displayed.

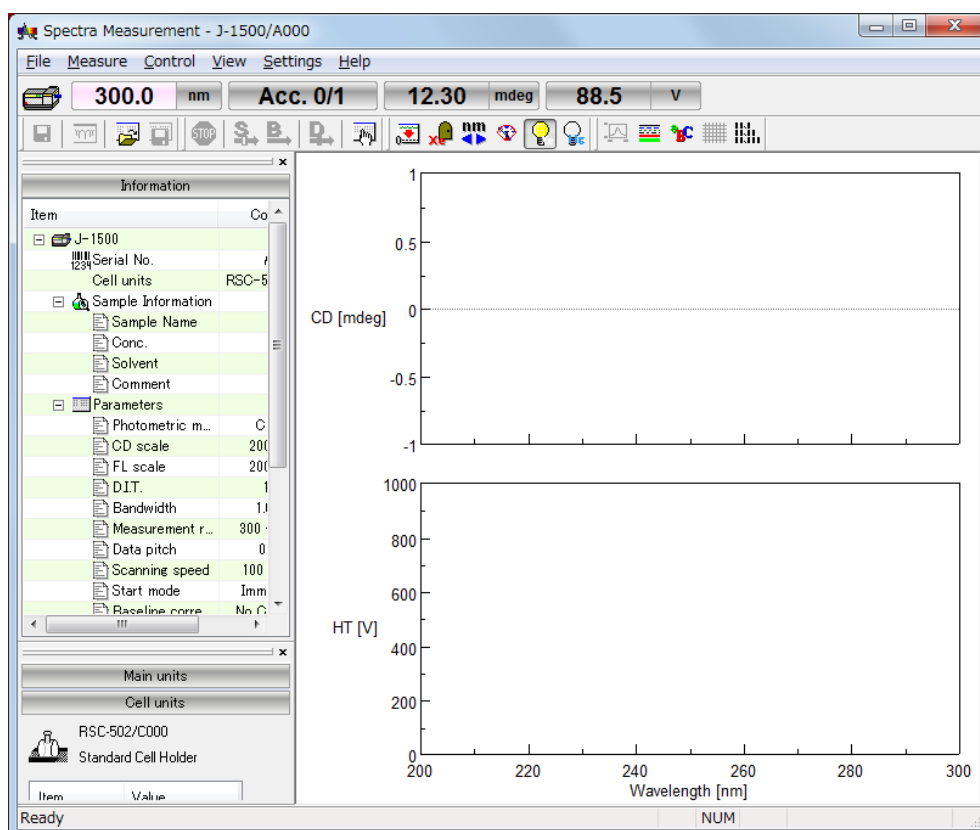


Figure 4.2 [Spectra Measurement] window

Note: In the CFR version, if [Monitor] is deselected in the [System Policy Settings] dialog box in the [Administrative Tools] program, "Off" is displayed in the monitor bar of the program main window. This setting can be selected only by users with [Spectra Manager CFR] Administrator privileges. For details about this setting, refer to the [System Policy Settings...] section in the instruction manual for the [Spectra Manager Ver.2.5 CFR Administrative Tools] program.

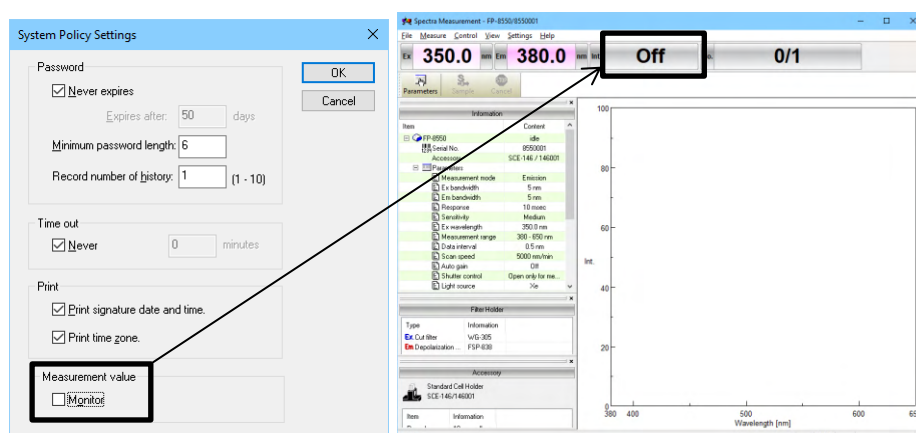



Figure 4.3 Clearing [Monitor] in [System Policy Settings] dialog box and displaying Off in the monitor bar in the CFR version

4.3 Specifying and Saving Measurement Parameters

- (1) Select [Parameters...] from the [Measure] menu (or click the  button) to display the dialog box in Figure 4.4. The [Parameters] dialog box contains the [General], [Control], [Information], and [Data] tabs. The [General] tab is open by default in the [Parameters] dialog box. The page can be changed by clicking each tab at the top of the dialog box.

Notes: In the CFR version,

1. *Managers or Analysts authority is required to specify parameters.*
2. *In order to perform a measurement by users with Operators authority, a measurement parameter file with an approval signature is required.*
3. Table 4.1 lists the privileges of different users with regard to inspection parameter files.

Table 4.1 Parameter file signature status and user privileges to perform measurements

Inspection parameter file			Measurement		
Creation	Review	Approval	Managers	Analysts	Operators
Managers Analysts	Managers Analysts	Managers			
■	□	□	Possible (Only with creator's own parameter file)	Possible (Only with creator's own parameter file)	Impossible
■	■	□	Impossible	Impossible	Impossible
■	■	■	Possible	Possible	Possible

■: Signed, □: Not signed

The parameters can be switched between two modes: Basic mode and Advanced mode. In Basic mode, the choice of parameters is partially limited. In Advanced mode, each parameter is specified individually. The parameters described in this section are specified in Basic mode. To switch between Basic and Advanced modes, click the <Basic Mode> or <Advanced Mode> button.

(2) Specify the measurement parameters in the [General] tab as follows.

Parameters Basic

General Control Information Data

Photometric mode

Channels Num: 2

Channel 1: CD

Channel 2: HT

Channel 3: HT

Channel 4: UV single(Abs)

Start: 350 nm

End: 220 nm

Data pitch: 0.1 nm

Start mode: Immediately

Scanning mode: Continuous

Scanning speed: 100 nm/min

CD scale: 200 mdeg/0.1 dOD

FL scale: 200 mdeg/0.1 dOD

D.I.T.: 2 sec

Bandwidth: 1.00 nm

Accumulation/cycle

☐ Accumulation

No. of cycles: 1

Advanced Mode Open... Save... Default OK Cancel

Figure 4.4 [General] tab in [Parameters] dialog box

Photometric mode	
[Channels Num]	2
[Channel 1]	[CD]
[Channel 2]	[HT]
[CD scale]	200 mdeg/1.0 dOD
[FL scale]	200 mdeg/1.0 dOD
[D. I. T.]	2 sec
[Bandwidth]	1 nm
[Start]	350 nm
[End]	220 nm
[Data pitch]	0.1 nm
[Start mode]	[Immediately]
[Scanning mode]	[Continuous]
[Scanning speed]	100 nm/min
Accumulation/cycle	
[Accumulation]	Deselected
[No. of accumulations]	1

Note: Select the appropriate [CD scale] for the sample to be measured.

Parameters can be set by using two methods: selecting from available options and specifying numeric values and characters. An example of setting parameters by using each method is given below.

1) Selecting from available options

The following example describes how to select the CD scale. Select the CD scale from the available options in the [CD scale] drop-down list. Click the arrow to the right of the box to display the full list of [CD scale] options (see Figure 4.5). Select the desired CD scale (in this example, 200 mdeg/1.0 dOD).



Figure 4.5 [CD scale] drop-down list

2) Specifying numeric values and characters

The following example describes how to specify the start wavelength. Click the [Start] text box to edit the start wavelength. When the cursor appears in the text box, values can be entered. Type the desired wavelength (in the example in Figure 4.4, 350 nm) in the text box to set the start wavelength.

(3) Specify the measurement parameters in the [Control] tab as follows.

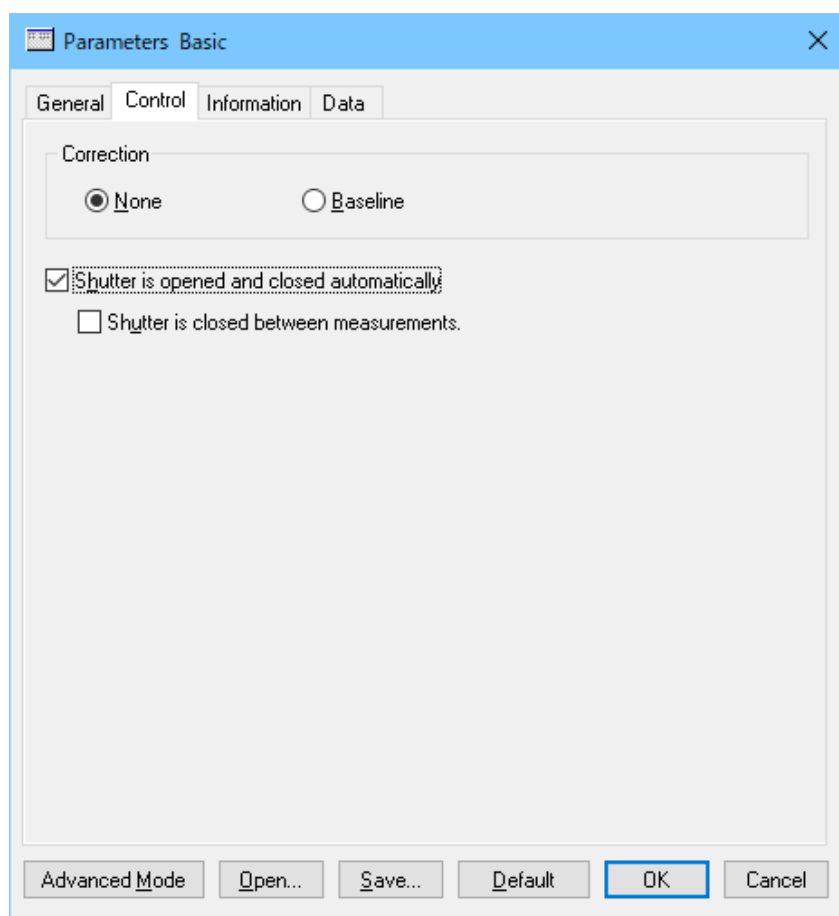


Figure 4.6 [Control] tab in [Parameters] dialog box

Correction	[None]
[Shutter is opened and closed automatically]	Selected
[Shutter is closed between measurements.]	Deselected

- (4) Click the [Information] tab. Type information in the [Sample name], [Username], [Division], [Comment], [Concentration] and [Solvent] text boxes as desired. The information entered here is saved as comment information together with the spectral data when the spectra are saved.

Figure 4.7 [Information] tab in [Parameters] dialog box

[Sample name]	Enter as desired (maximum 63 characters).
[Username]	Enter a username (maximum 63 characters).
[Division]	Enter a division name for the operator (maximum 127 characters).
[Comment]	Enter a comment (maximum 127 characters).
[Concentration]	Enter a sample concentration.
[Solvent]	Enter a solvent name (maximum 127 characters).
[Display the [Comments] dialog box before measurement]	Deselected

Note: In the CFR version, the [Username] and [Division] text boxes cannot be changed by the user.

- (5) Click the [Data] tab. Specify the [Save data], [Send data], and [Print] options for the measured spectrum. In this example, select the [To Spectra Analysis] check box.

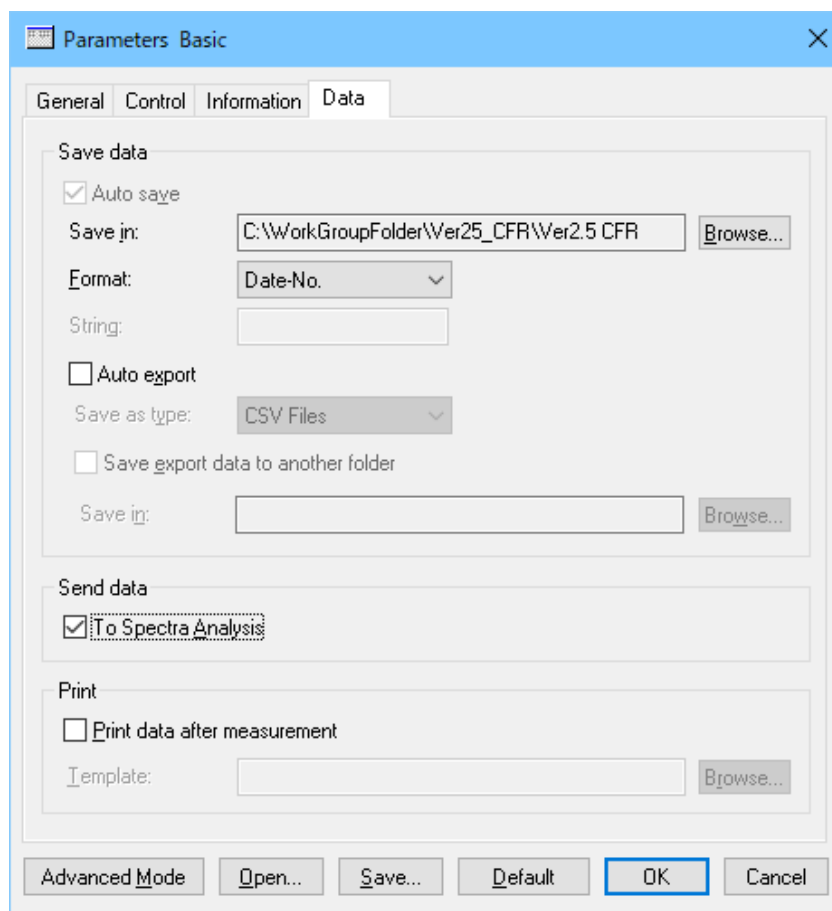


Figure 4.8 [Data] tab in [Parameters] dialog box

[To Spectra Analysis]	Selected
[Print data after measurement]	Deselected

Note : In the CFR version, all measurement data must be saved. Select the folder specified for the workgroup that the user belongs to, or a subfolder.

- (6) In the CFR version, always save changes to the parameters. Click the <Save...> button to display the dialog box in Figure 4.9.

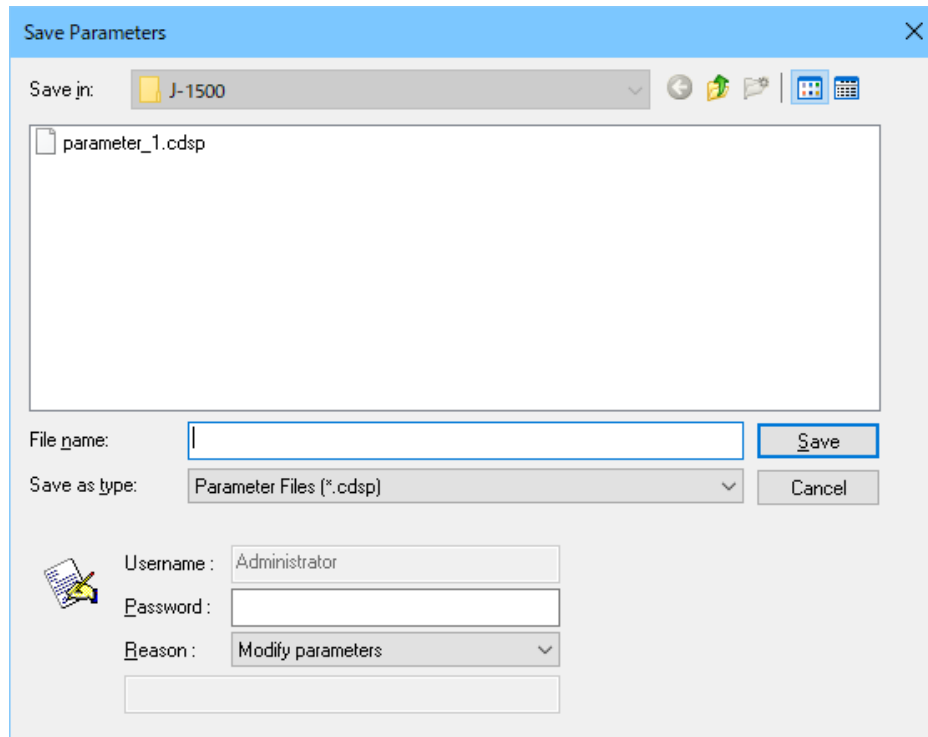


Figure 4.9 [Save Parameters] dialog box

- (7) After entering the file name, type a password and then click the <Save> button.

Note: In the CFR version, when the parameters have not been saved, the following message is displayed.

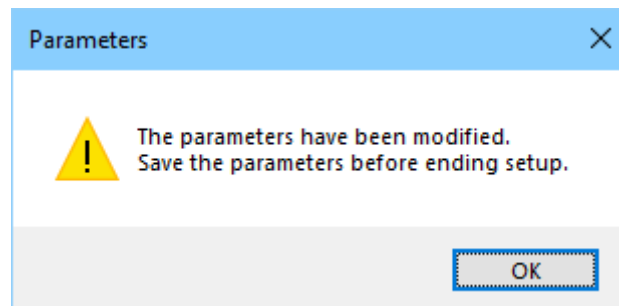


Figure 4.10 Message when an attempt was made to close the parameter file without saving it

4.4 Reviewing and Approving the Parameter File

Note: In the CFR version, Managers or Analysts authority is required for the review signature, and Managers authority is required for the approval signature. Log on to [Spectra Manager] as a user who has authority to provide a signature.

- (1) Click [Spectra Measurement] in the [Instrument] list of [Spectra Manager CFR].
- (2) Select [Parameter Files (*.cdspp)] in the [Files of type] drop-down list to display the file name of a saved calibration curve. Clicking the calibration file name displays the [Information], [History], and [Signature] tabs. Click the [Signature] tab.

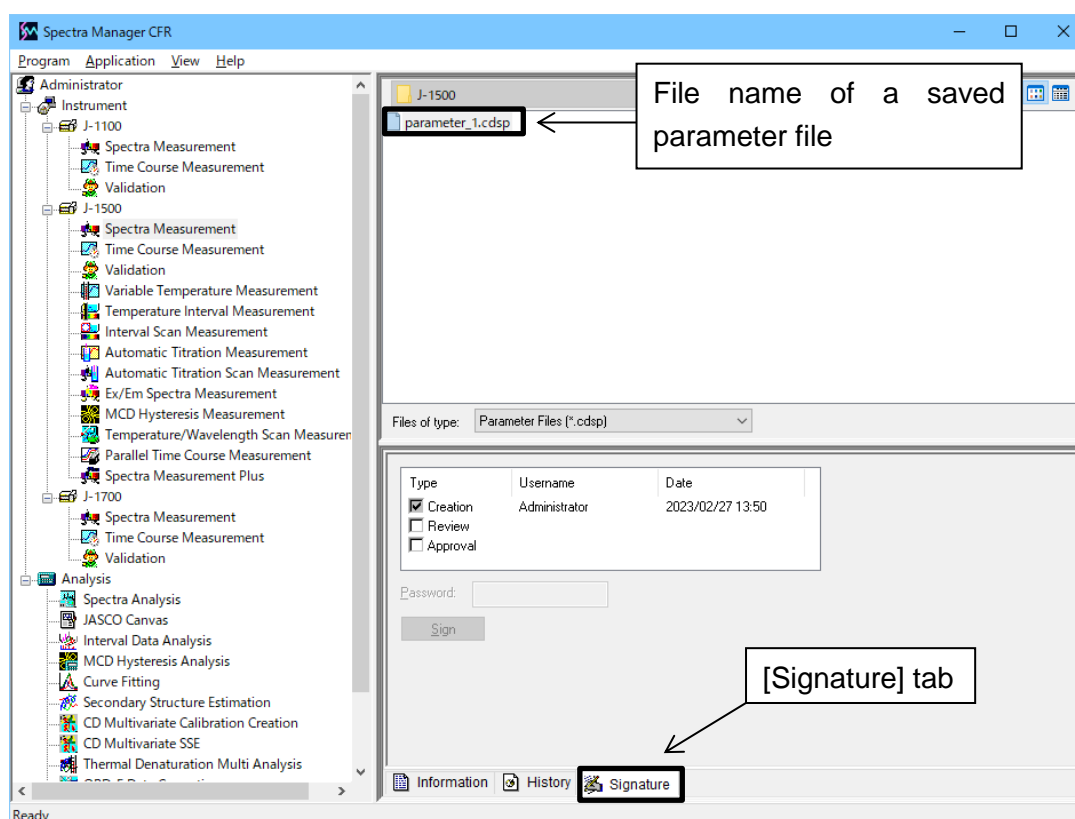


Figure 4.11 Assigning signatures to parameter file

- (3) Select the [Review] and/or [Approval] check boxes, type the password, and then click the <Sign> button.

- (4) Start the [Spectra Measurement] program, set the data mode and select [Open Parameters...] from the [File] menu. The dialog box in Figure 4.12 is displayed.

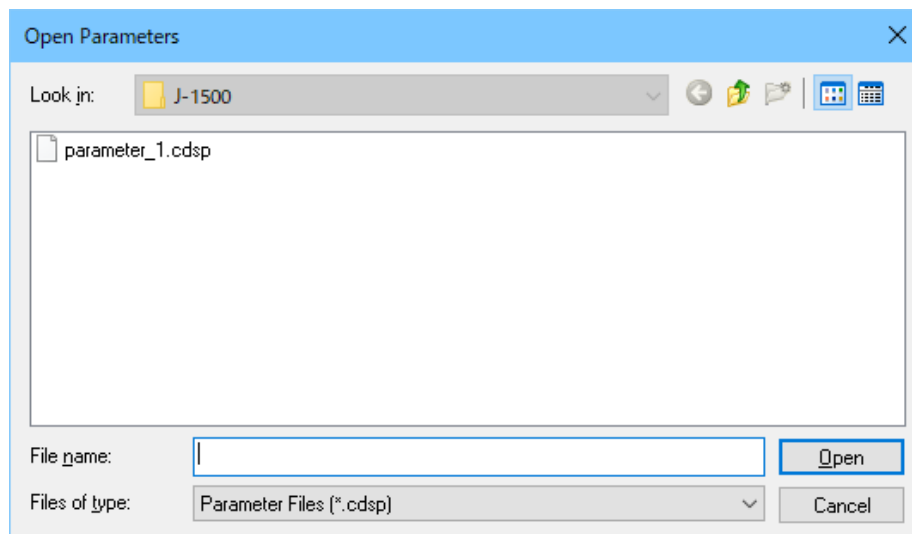


Figure 4.12 [Open Parameters] dialog box

Note: In the CFR version, a folder higher than a workgroup folder cannot be used. However, a subfolder within a prespecified workgroup folder can be selected.


- (5) Select a parameter file and click the <Open> button. The parameter file opens in the [Spectra Measurement] window.

4.5 Measuring the Sample

(1) Inserting the sample

Fill the cell with the sample, place it in the cell holder, and then close the lid of the sample compartment.

(2) Measuring the sample

Select [Sample Measurement] from the [Measure] menu (or click the  button) to measure the spectrum for the sample. In the CFR version, the dialog box shown in Figure 4.13 is displayed at the start of the measurement. Type a password, and then click the <OK> button. The progress of the measurement is displayed in the window (see Figure 4.14).

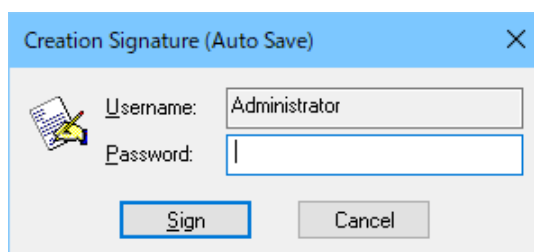


Figure 4.13 [Creation Signature (Auto Save)] dialog box

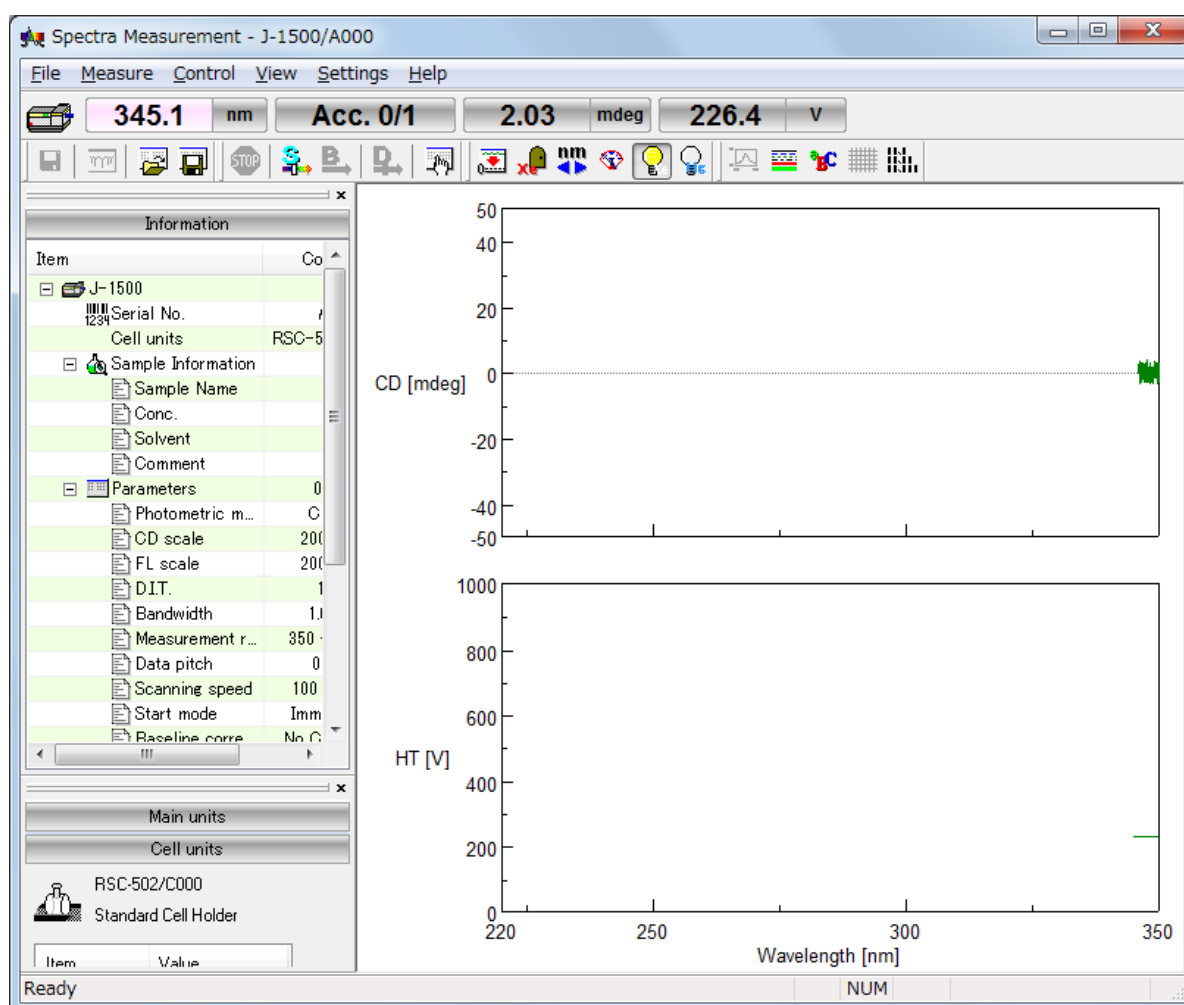



Figure 4.14 [Spectra Measurement] window

After measurement, the [Spectra Analysis] program starts and the measured data is displayed in a window. This window is called a "View".

Note 1: When the [Spectra Analysis] program is already running, it is not automatically displayed as the active window after measurement. Click [Spectra Analysis] on the Windows task bar to display the program as the active window.

Note 2: Data is not automatically transferred unless the [To Spectra Analysis] check box is selected in the [Data] tab of the [Parameters] dialog box. To transfer data, select [Send to Analysis] from the [File] menu (or click the  button).

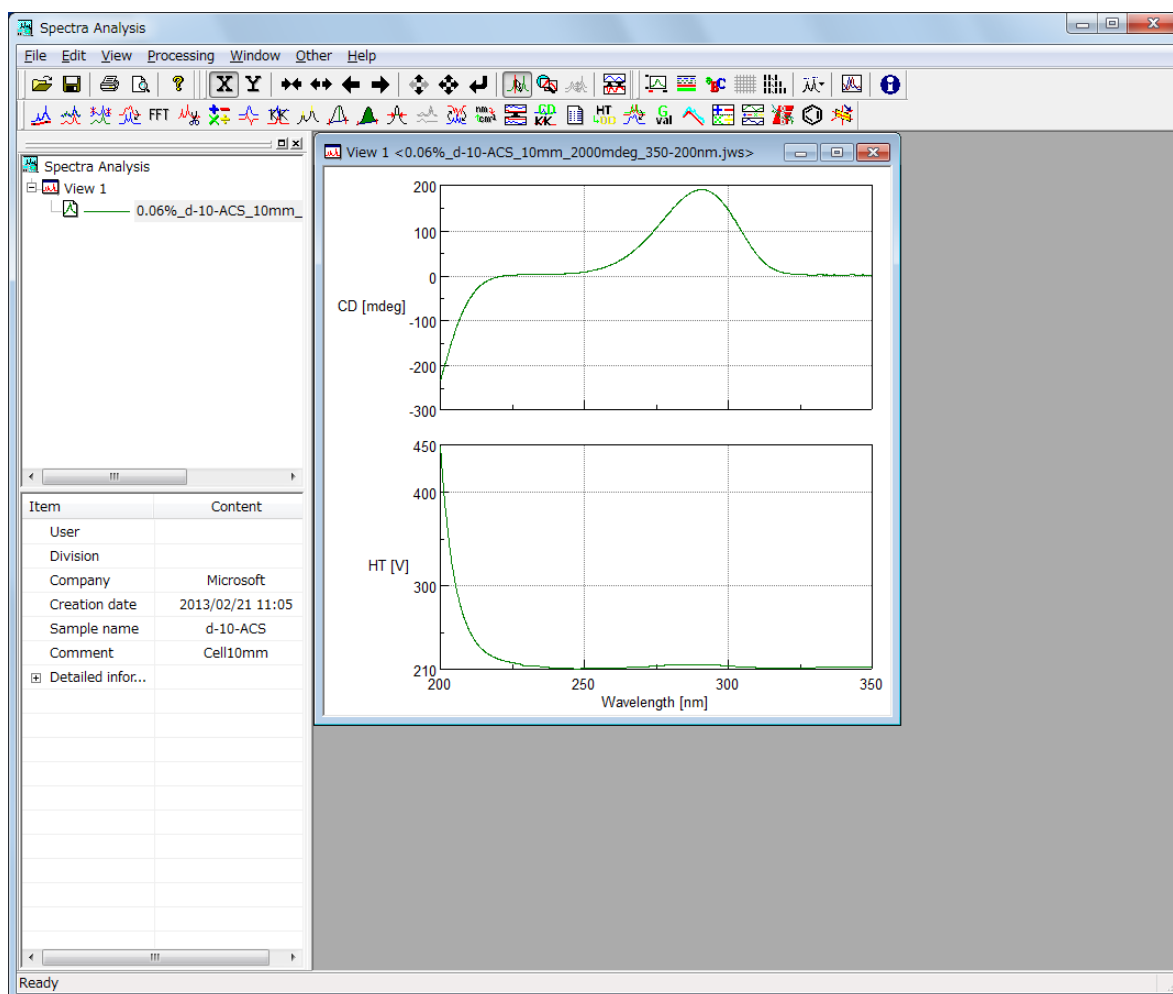



Figure 4.15 Spectrum View

4.6 Saving Spectra

This section describes the procedure for saving spectra using the [Spectra Analysis] program. In the CFR version, since all measurement data are always saved, the steps in this section are not required.

Note: Spectra can also be saved by selecting [Save Data...] from the [File] menu (or by clicking the  button) in the [Spectra Measurement] program.

- (1) Select [Save As...] from the [File] menu in the [Spectra Analysis] program to display the dialog box in Figure 4.16.

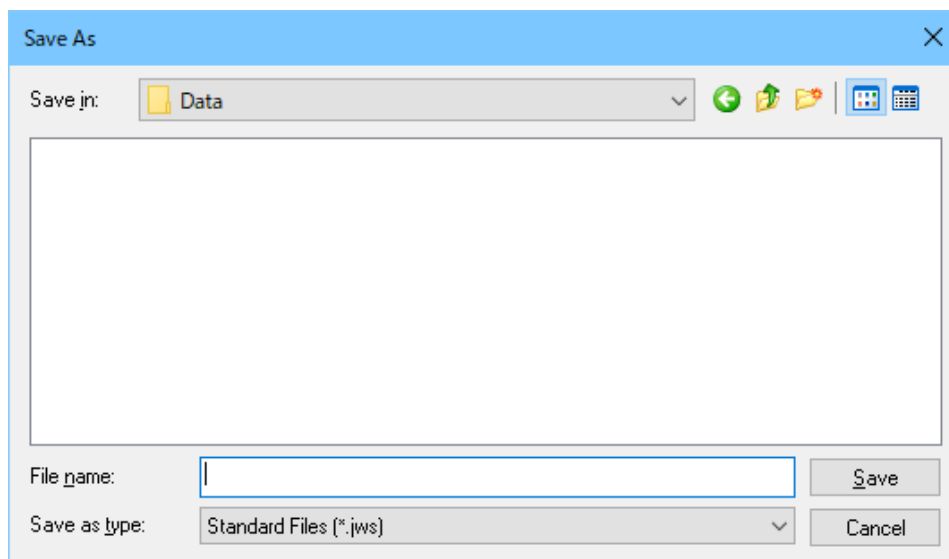


Figure 4.16 [Save As] dialog box

- (2) From the [Save in] drop-down list, select the folder in which to save the spectrum.
- (3) Type the file name in the [File name] text box. In this example, type “d-10-ACS” as the file name.
- (4) After entering the file name, type a password and then click the <Save> button.
- (5) The file name is appended with the extension “.jws” (“d-10-ACS.jws”), which is the standard file type.
- (6) After saving the data, the title bar of the View changes to “View (d-10-ACS.jws)”.

4.7 Printing Results

Print the measured spectra using a connected printer as follows.

- (1) Select [Print Setup...] from the [File] menu to display the dialog box in Figure 4.17. The content of this dialog box varies depending on the connected printer. Configure the printer settings if necessary.

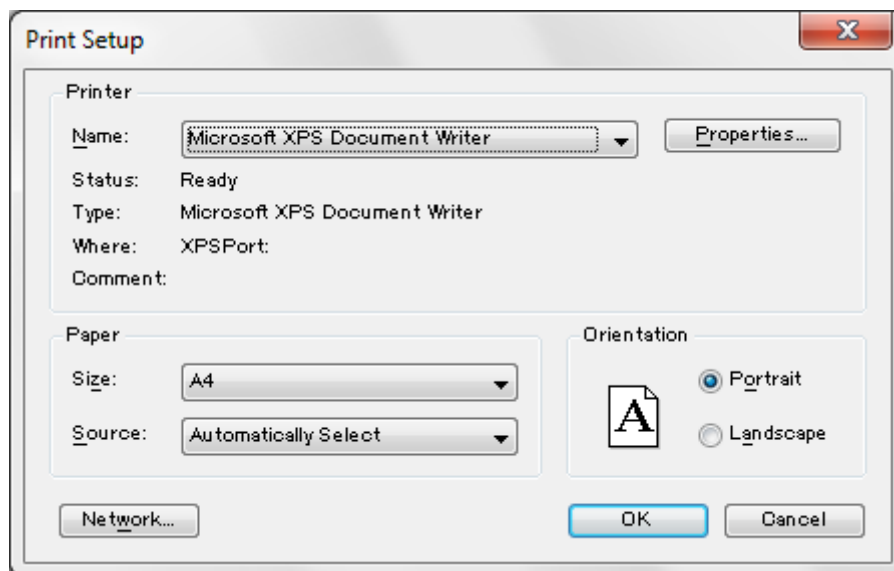


Figure 4.17 Example of [Print Setup] dialog box

- (2) Select [Print...] from the [File] menu to print the data.

Note: Refer to the [Spectra Analysis] program manual for details about printing a spectrum.

4.8 Exiting the Program

(1) Exiting the [Spectra Analysis] program

Select [Exit] from the [File] menu. The [Spectra Analysis] window closes, leaving the [Spectra Measurement] window displayed.

Note: When there are unsaved spectra in the window, a warning message is displayed. Perform the action recommended by the message. A message is displayed for each unsaved spectrum.

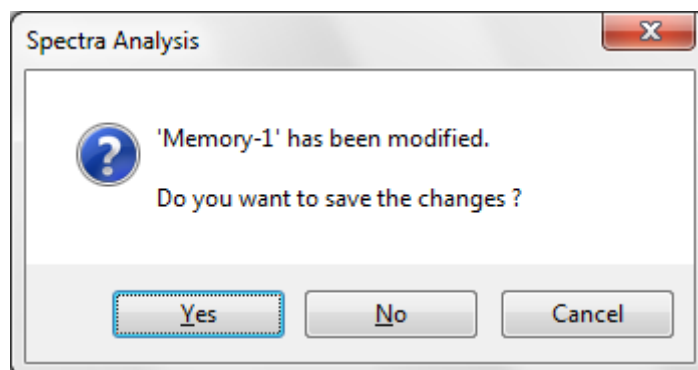


Figure 4.18 Message when data has not been saved

(2) Exiting the [Spectra Measurement] program

Select [Exit] from the [File] menu to close the [Spectra Measurement] program, leaving the [Spectra Manager] window displayed.

(3) Exiting the [Spectra Manager] program

Select [Exit] from the [Program] menu.

(4) Exiting Windows

Exit Windows in accordance with standard Windows operation procedures.

(5) Turning off the PC and CD spectrometer

First, turn off the PC and monitor. Ensure that the sample compartment is empty, and then turn off the CD spectrometer.

5. Introduction to the [Time Course Measurement] Program

This chapter describes how to use the [Time Course Measurement] program, including procedures for starting the [Time Course Measurement] program, measuring the spectrum of a sample, and saving and printing data, for practice in operating a JASCO CD spectrometer. Explanations of the parameters have been kept to a minimum, with a focus on the flow of operations. Follow the measurement procedures outlined below for practice in operating the CD spectrometer. Refer to Chapter 6 “Measurement Program Reference” for details about toolbar icons and menus, and refer to Chapter 8 “[Time Course Measurement] Program Menu Reference” for details about the [Parameters...] command.

5.1 Procedures for the [Time Course Measurement] Program

This section describes procedures ranging from measuring CD time-course data for an ammonium d-10-camphorsulfonate solution to saving and printing the measurement data.

Starting the [Time Course Measurement] Program	Refer to Section 5.2
↓	
Specifying and Saving Measurement Parameters	Refer to Section 5.3
↓	
Reviewing and Approving the Parameter File*	Refer to Section 5.4
↓	
Measuring the Sample	Refer to Section 5.5
↓	
Saving Time-Course Data	Refer to Section 5.6
↓	
Printing Results	Refer to Section 5.7
↓	
Exiting the Program	Refer to Section 5.7

*The step in section 5.4 is only for the CFR version and 5.6 is only for the non CFR version

Note 1: In the CFR version, Managers or Analysts authority is required to specify measurement parameters.

Note 2: In the CFR version, measurement can basically be performed using only reviewed and approved measurement parameters. However, only the person who created parameter file can perform measurement without review and approval.

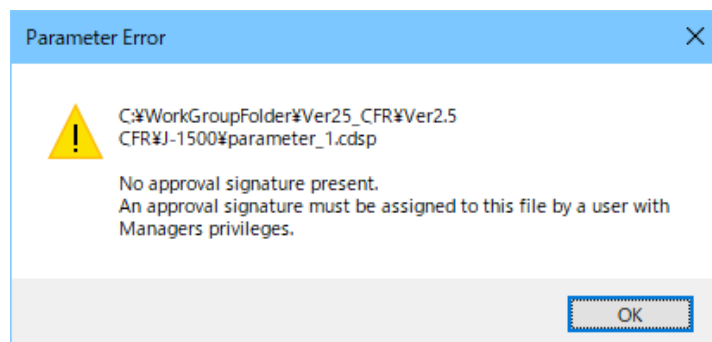


Figure 5.1 Message when measurement is attempted using unapproved and unsigned measurement parameters

5.2 Starting the [Time Course Measurement] Program

To start the program, double-click [Time Course Measurement] from the [Instrument] list of [Spectra Manager CFR]. The window in Figure 5.2 is displayed.

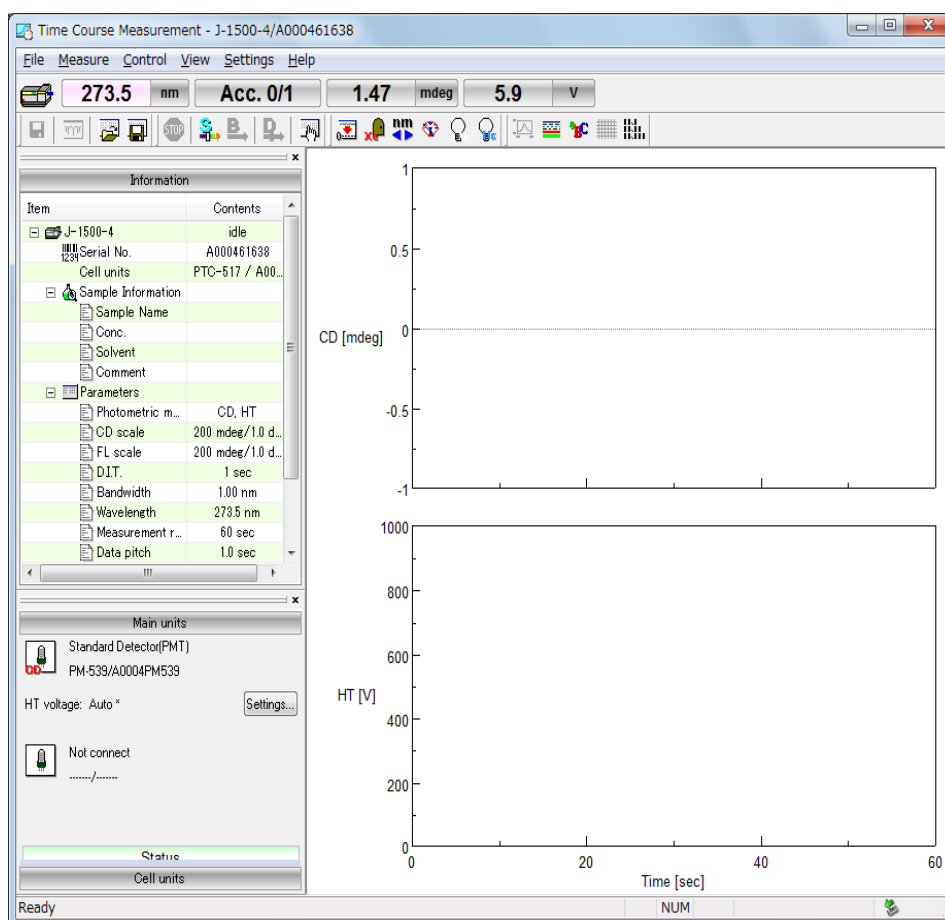


Figure 5.2 [Time Course Measurement] window

Note: In the CFR version, if [Monitor] is deselected in the [System Policy Settings] dialog box in the [Administrative Tools] program, “Off” is displayed in the monitor bar of the program main window. This setting can be selected only by users with [Spectra Manager CFR] Administrator privileges. For details about this setting, refer to the [System Policy Settings...] section in the instruction manual for the [Spectra Manager Ver.2.5 CFR Administrative Tools] program.

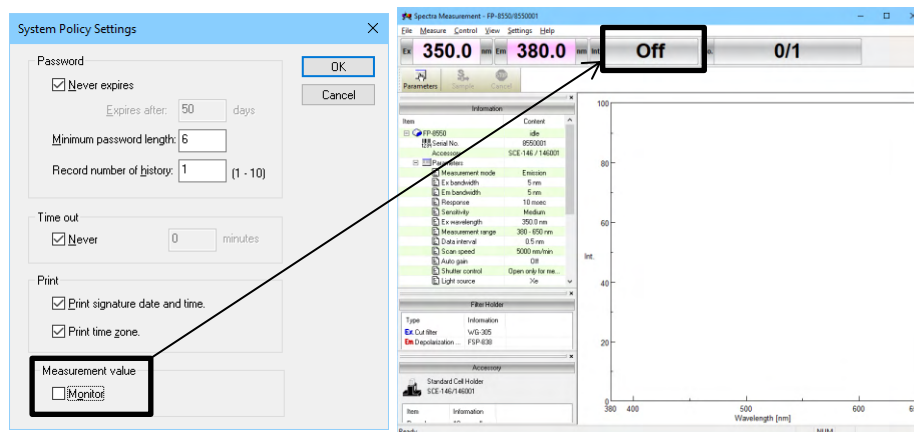



Figure 5.3 Clearing [Monitor] in [System Policy Settings] dialog box and displaying Off in the monitor bar in the CFR version

5.3 Specifying and Saving Measurement Parameters

- (1) Select [Parameters...] from the [Measure] menu (or click the  button) to display the dialog box in Figure 5.4. The [Parameters] dialog box contains the [General], [Control], [Information], and [Data] tabs. The [General] tab is open by default in the [Parameters] dialog box. The page can be changed by clicking each tab at the top of the dialog box.

Notes: In the CFR version,

1. Managers or Analysts authority is required to specify parameters.
2. In order to perform a measurement by users with Operators authority, a measurement parameter file with an approval signature is required.
3. Table 4.1 lists the privileges of different users with regard to inspection parameter files.

The parameters can be switched between two modes: Basic mode and Advanced mode. In Basic mode, the choice of parameters is partially limited. In Advanced mode, each parameter is specified individually. The parameters described in this section are specified in Basic mode. To switch between Basic and Advanced modes, click the <Basic Mode> or <Advanced Mode> button.

(2) Specify the measurement parameters in the [General] tab as follows.



Figure 5.4 [General] tab in [Parameters] dialog box

Photometric mode	
[Channels Num]	2
[Channel 1]	[CD]
[Channel 2]	[HT]
[CD scale]	200 mdeg/1.0 dOD
[FL scale]	200 mdeg/1.0 dOD
[D. I. T.]	1 sec
[Bandwidth]	1 nm
[Wavelength]	291 nm
[Measuring time]	60 sec
[Data pitch]	1 sec
[Start mode]	[Immediately]

Note: Select the appropriate [CD scale] for the sample to be measured.

Parameters can be set by using two methods: selecting from available options and specifying numeric values and characters. An example of setting parameters using each method is given below.

1) Selecting from available options

The following example describes how to select the CD scale. Select the CD scale from the available options in the [CD scale] drop-down list. Click the arrow to the right of the box to display the full list of [CD scale] options (see Figure 5.5). Select the desired CD scale (in this example, 200 mdeg/1.0 dOD).

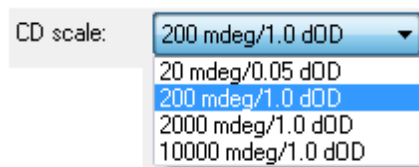


Figure 5.5 [CD scale] drop-down list

2) Specifying numeric values and characters

The following example describes how to specify the wavelength. Click the [Wavelength] text box to edit the wavelength. When the cursor appears in the text box, values can be entered. Type the desired wavelength (in the example in Figure 5.4, 291 nm) in the text box to set the wavelength.

(3) Specify the measurement parameters in the [Control] tab as follows.

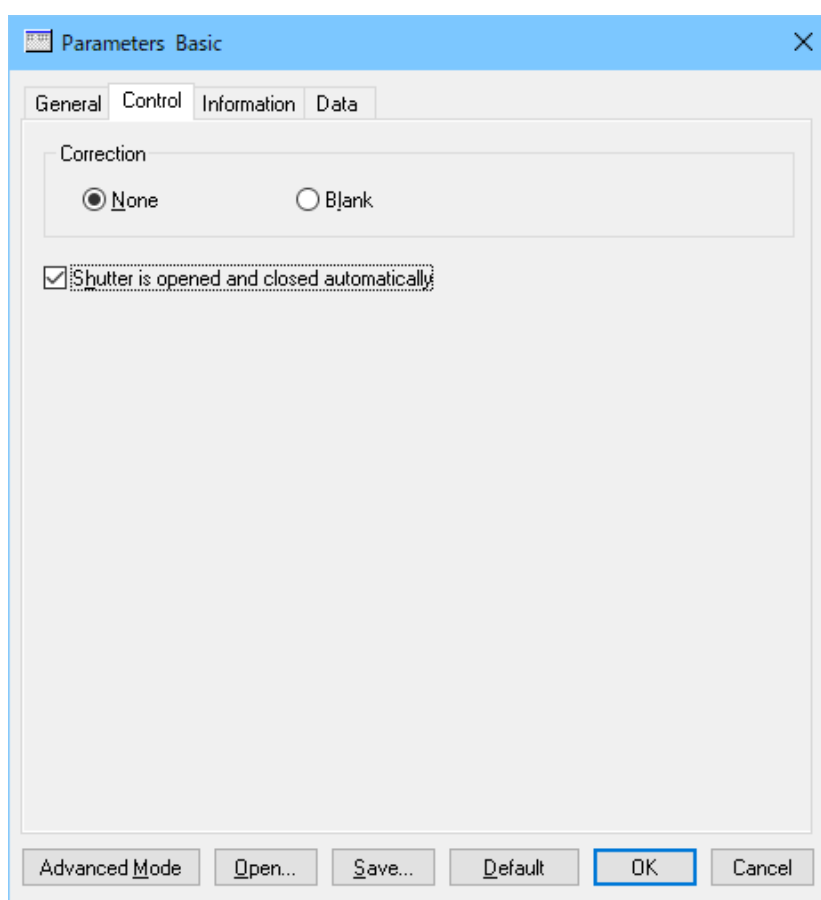


Figure 5.6 [Control] tab in [Parameters] dialog box

Correction	[None]
[Shutter is opened and closed automatically]	Selected

- (4) Click the [Information] tab. Type information in the [Sample name], [Username], [Division], [Comment], [Concentration] and [Solvent] text boxes as desired. The information entered here is saved as comment information together with the time-course data when the data are saved.

Figure 5.7 [Information] tab in [Parameters] dialog box

[Sample name]	Enter as desired (maximum 63 characters).
[Username]	Enter a username (maximum 63 characters).
[Division]	Enter a division name for the operator (maximum 127 characters).
[Comment]	Enter as desired (maximum 127 characters).
[Concentration]	Enter as desired.
[Solvent]	Enter as desired (maximum 127 characters).
[Display the [Comments] dialog box before measurement]	Deselected

Note: In the CFR version, the [Username] and [Division] text boxes cannot be changed by the user.

- (5) Click the [Data] tab. Specify the [Save data], [Send data], and [Print] options for the measured data. In this example, select the [To Spectra Analysis] check box.

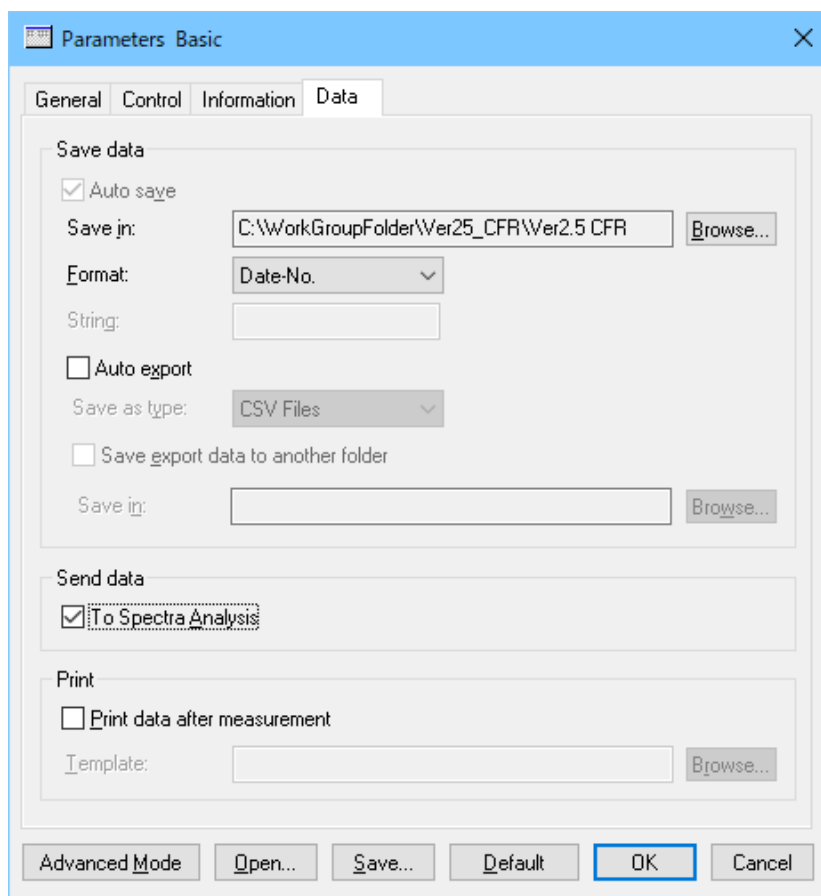


Figure 5.8 [Data] tab in [Parameters] dialog box

[Auto save]	Deselected
[To Spectra Analysis]	Selected
[Print data after measurement]	Deselected

Note: In the CFR version, all measurement data must be saved. Select the folder specified for the workgroup that the user belongs to, or a subfolder.

- (6) In the CFR version, always save changes to the parameters. Click the <Save...> button to display the dialog box in Figure 5.9.

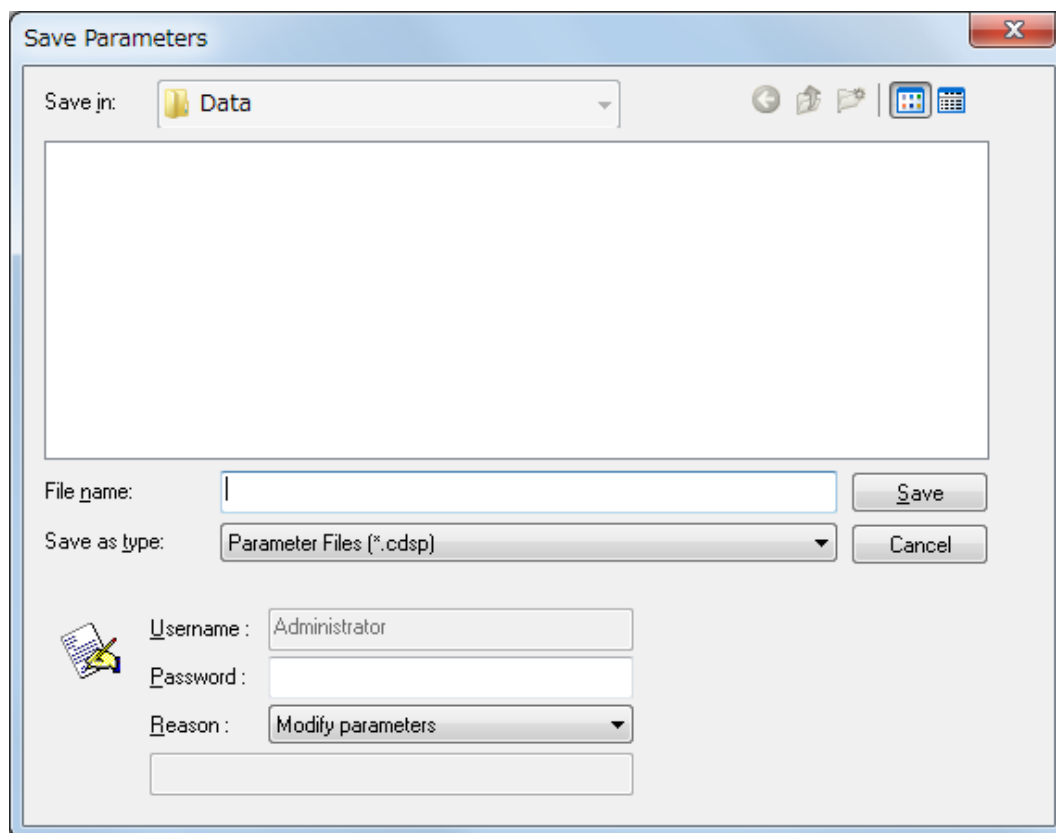


Figure 5.9 [Save Parameters] dialog box

- (7) After entering the file name, type a password and then click the <Save> button.

Note: When the parameters have not been saved, the following message is displayed.

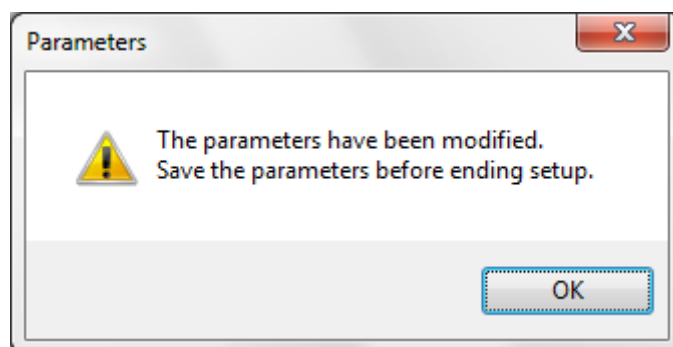


Figure 5.10 Message when the parameters have not been saved

5.4 Reviewing and Approving the Parameter File

Note : In the CFR version, Managers or Analysts authority is required for the review signature, and Managers authority is required for the approval signature. Log on to [Spectra Manager] as a user who has authority to provide a signature.

- (1) Click [Time Course Measurement] in the [Instrument] list of [Spectra Manager CFR].
- (2) Select [Parameter Files (*.cdtm)] in the [Files of type] drop-down list to display the file name of a saved calibration curve. Clicking the calibration file name displays the [Information], [History], and [Signature] tabs. Click the [Signature] tab.

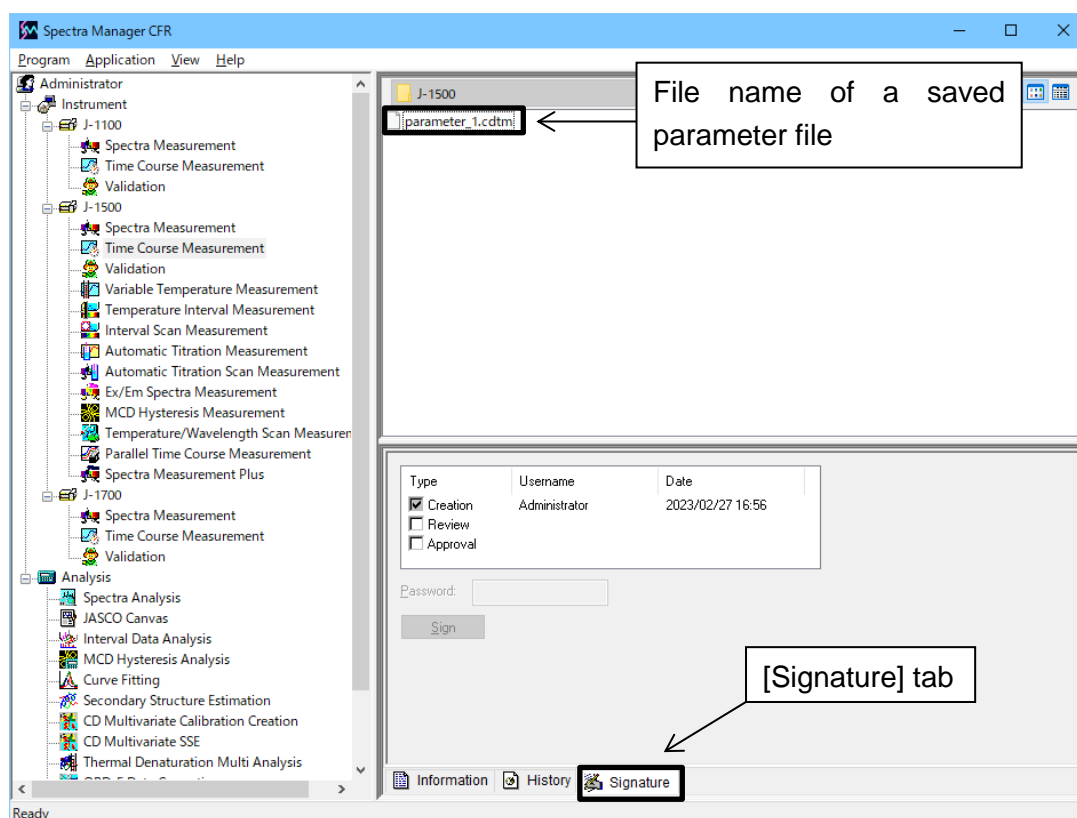


Figure 5.11 Assigning signatures to parameter file

- (3) Select the [Review] and/or [Approval] check boxes, type the password, and then click the <Sign> button.

- (4) Start the [Time Course Measurement] program, set the data mode and select [Open Parameters...] from the [File] menu. The dialog box in Figure 5.12Figure 4.12 is displayed.

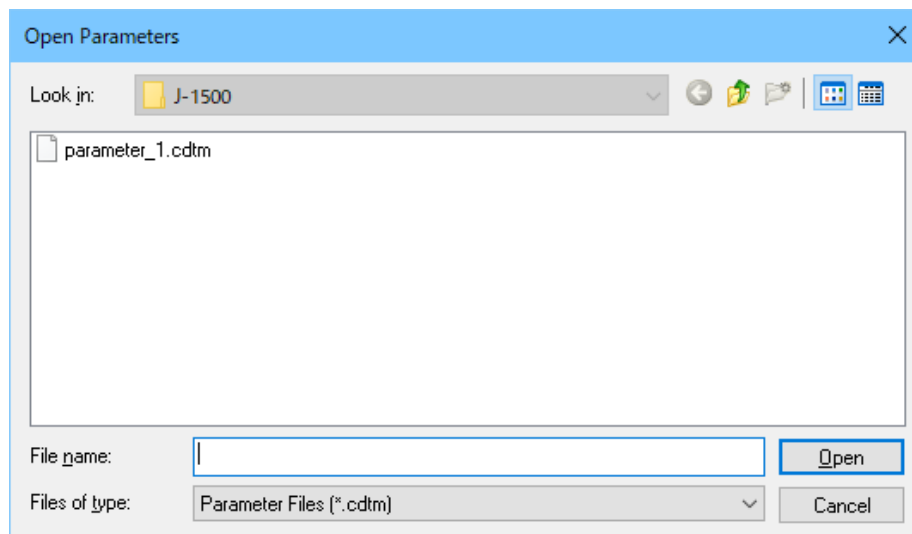


Figure 5.12 [Open Parameters] dialog box

Note: In the CFR version, a folder higher than a workgroup folder cannot be used. However, a subfolder within a prespecified workgroup folder can be selected.


- (5) Select a parameter file and click the <Open> button. The parameter file opens in the [Spectra Measurement] window.

5.5 Measuring the Sample

(1) Inserting the sample

Fill the cell with the sample, place it in the cell holder, and then close the lid of the sample compartment.

(2) Measuring the sample

Select [Sample Measurement] from the [Measure] menu (or click the  button) to measure the time-course data for the sample. In the CFR version, the dialog box shown in Figure 5.13 is displayed at the start of the measurement. Type a password, and then click the <OK> button. The progress of the measurement is displayed in the window (see Figure 5.14).

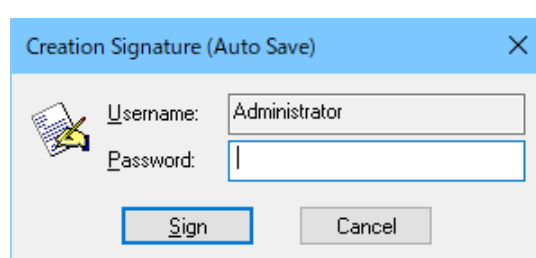


Figure 5.13 [Creation Signature (Auto Save)] dialog box

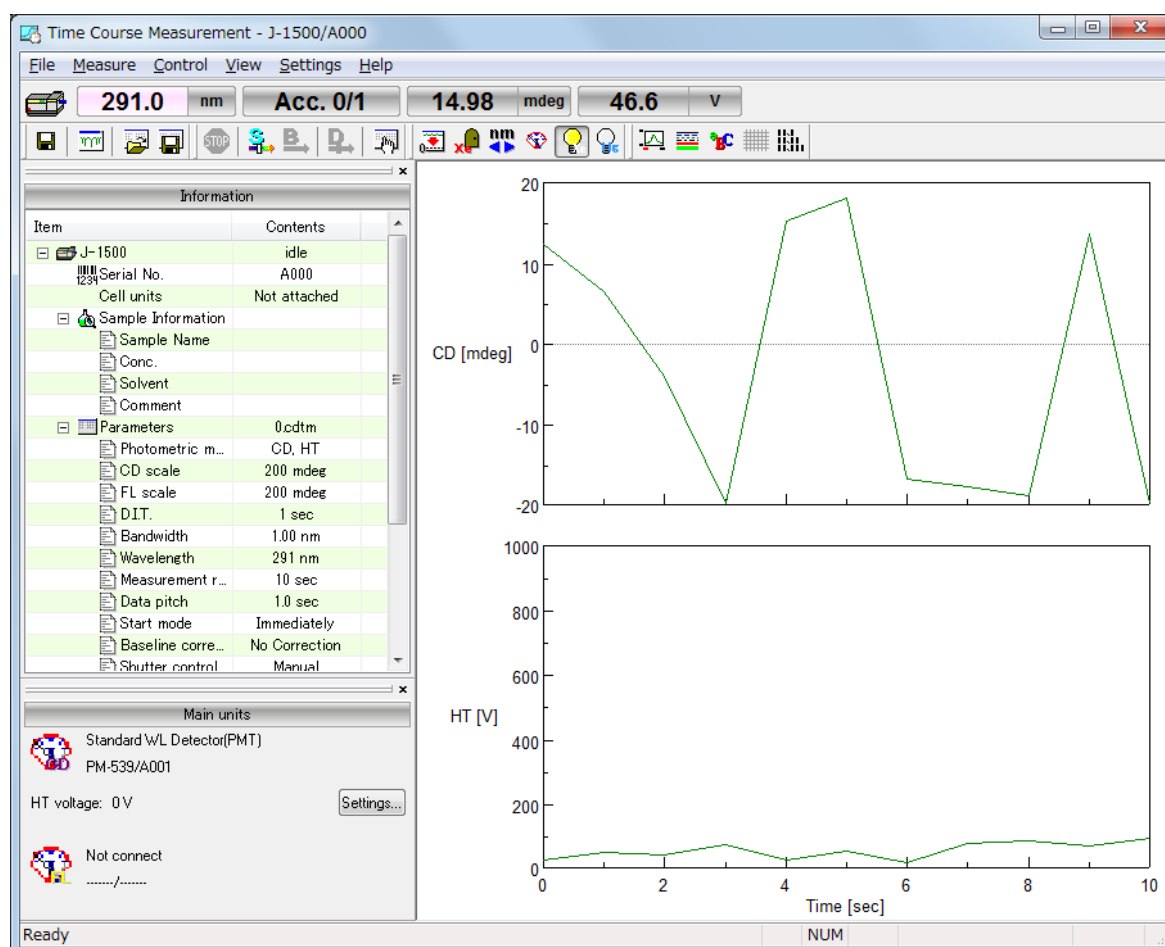



Figure 5.14 [Spectra Measurement] window

After measurement, the [Spectra Analysis] program starts and the measured data is displayed in the window. This window is called a "View".

Note 1: When the [Spectra Analysis] program is already running, it is not automatically displayed as the active window after measurement. Click [Spectra Analysis] on the Windows task bar to display the program as the active window.

Note 2: Data is not automatically transferred unless the [To Spectra Analysis] check box is selected in the [Data] tab of the [Parameters] dialog box. To transfer data, select [Send to Analysis] from the [File] menu (or click the  button).

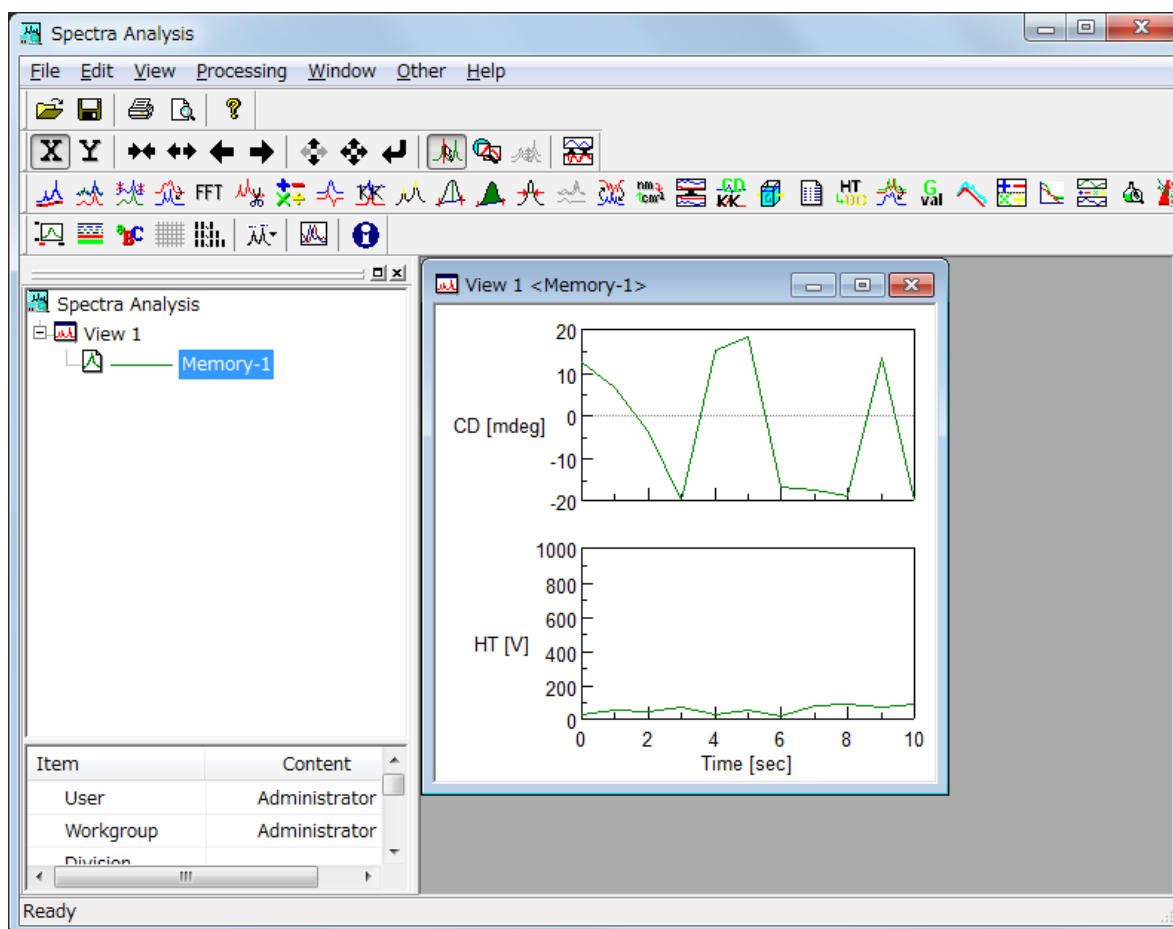



Figure 5.15 Spectrum View

5.6 Saving Time-Course Data

This section describes the procedure for saving time-course data using the [Spectra Analysis] program. In the CFR version, since all measurement data are always saved, the steps in this section are not required.

Note: Data can also be saved by selecting [Save Data...] from the [File] menu (or by clicking the  button) in the [Spectra Measurement] program.

- (1) Select [Save As...] from the [File] menu in the [Spectra Analysis] program to display the dialog box in Figure 5.16.

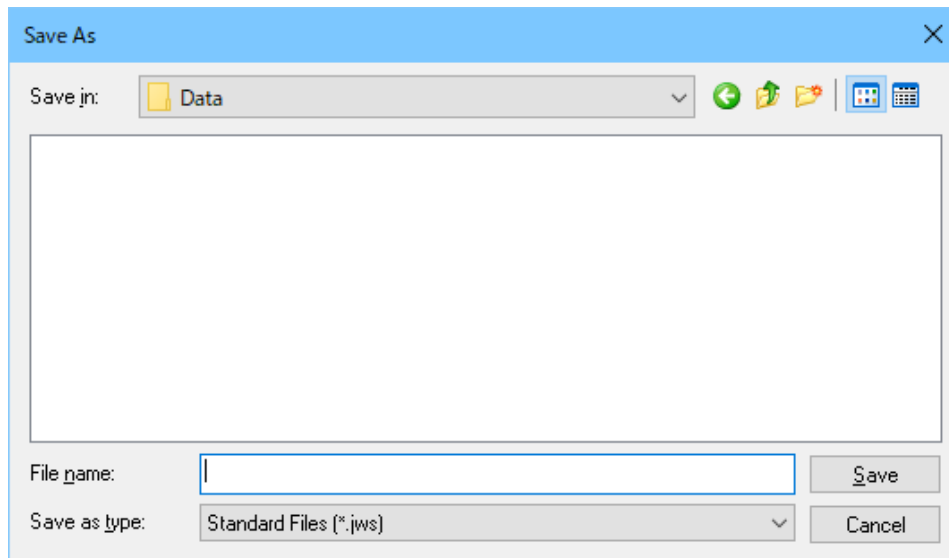


Figure 5.16 [Save As] dialog box

- (2) From the [Save in] drop-down list, select the folder in which to save the data.
- (3) Type the file name in the [File name] text box. In this example, type “d-10-ACS” as the file name.
- (4) After entering the file name, type a password and then click the <Save> button.
- (5) The file name is appended with the extension “.jws” (“d-10-ACS.jws”), which is the standard file type.
- (6) After saving the data, the title bar of the View changes to “View (d-10-ACS.jws)”.

5.7 Printing Results

Print the measured spectra using a connected printer as follows.

- (1) Select [Print Setup...] from the [File] menu to display the dialog box in Figure 5.17. The content of this dialog box varies depending on the connected printer. Configure the printer settings if necessary.

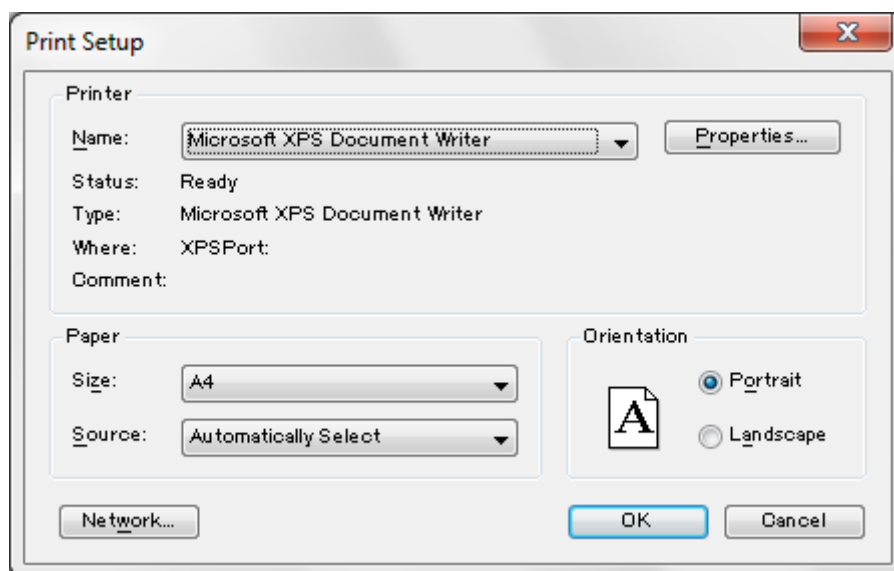


Figure 5.17 Example of [Print Setup] dialog box

- (2) Select [Print...] from the [File] menu to print the data.

Note: Refer to the [Spectra Analysis] program manual for details about printing a spectrum.

5.8 Exiting the Program

(1) Exiting the [Spectra Analysis] program

Select [Exit] from the [File] menu. The [Spectra Analysis] window closes, leaving the [Spectra Measurement] window displayed.

Note: When there are unsaved data in the window, a warning message is displayed. Perform the action recommended by the message. A message is displayed for each unsaved set of data.

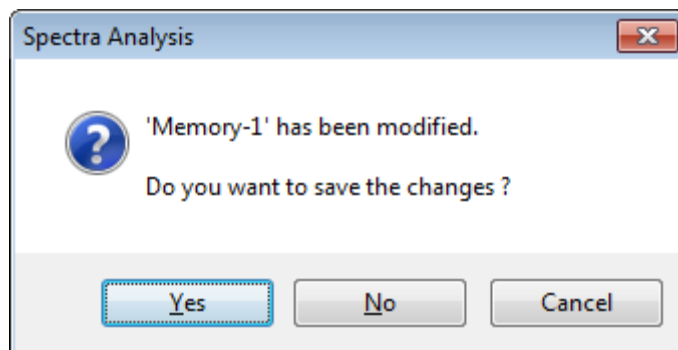


Figure 5.18 Message when data have not been saved

(2) Exiting the [Time Course Measurement] program

Select [Exit] from the [File] menu to close [Time Course Measurement], leaving the [Spectra Manager] window displayed.

(3) Exiting the [Spectra Manager] program

Select [Exit] from the [Program] menu.

(4) Exiting *Windows*

Exit *Windows* in accordance with standard *Windows* operation procedures.

(5) Turning off the PC and CD spectrometer

First, turn off the PC and monitor. Ensure that the sample compartment is empty, and then turn off the CD spectrometer.

6. Measurement Program Reference

This chapter describes common toolbar and menu commands for the [Spectra measurement], [Time Course Measurement] and [Data Monitor] programs. Refer to Chapters 7 to 9 or the [Help] menu for details about specific parameters for each measurement program.

Note 1: Refer to Table 6.1 for information about the availability of toolbar and menu commands in the three programs.

Note 2: All toolbar and menu commands are not necessarily available for every program.

Note 3: In the CFR version, the [Data Monitor] program is not available.

The [Spectra Measurement] program reference describes the features of the [Spectra Measurement] program for measuring a sample spectrum.

6.1 Overview of Measurement Programs

Starting the measurement program displays the window in Figure 6.1.

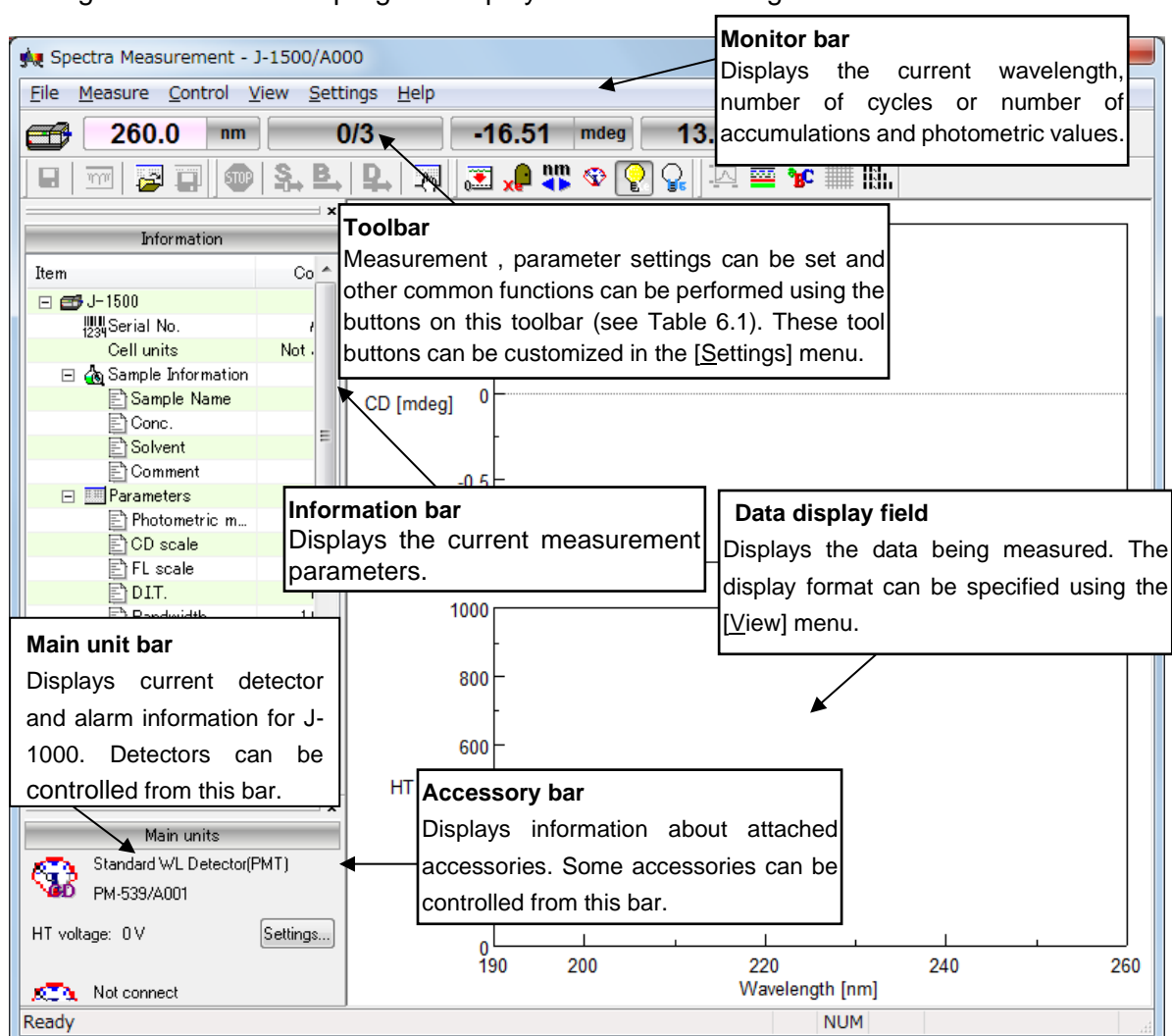





Figure 6.1 Measurement program window ([Spectra Measurement])

Table 6.1 Toolbar buttons and names

Button	Name	Reference Section	Spectra Measurement	Time Course Measurement	Data Monitor
	[Save]	6.2.1 [Save Data...] <div>Note: In the CFR version, this function is not available.</div>	✓	✓	
	[Send to Analysis]	6.2.2 [Send to Analysis]	✓	✓	
	[Open Parameters]	6.2.3 [Open Parameters...]	✓	✓	
	[Save Parameters]	6.2.4 [Save Parameters...] <div>Note: In the CFR version, this function is not available.</div>	✓	✓	
	[Cancel]	6.3.1 [Cancel]	✓	✓	
	[Sequence Measurement]	6.3.2 [Sequence Measurement]	✓	✓	
	[Sample Measurement]	6.3.3 [Sample Measurement]	✓	✓	
	[Blank/Baseline Measurement]	6.3.4 [Baseline Measurement] 6.3.5 [Blank Measurement]	✓	✓	
	[Dark Measurement]	6.3.6 [Dark Measurement]	✓	✓	
	[Parameters]	6.3.7 [Parameters...]	✓	✓	✓
	[Zero Clear]	6.4.1 [Zero Clear]	✓	✓	✓
	[Shutter]	6.4.2 [Shutter...]	✓	✓	✓
	[Move Wavelength]	6.4.3 [Move Wavelength...]	✓	✓	✓
	[Detector]	6.4.4 [Detector...]	✓	✓	✓
 	[Xe Lamp] [WI Lamp]	6.4.5 [Light Source...] <div>Note: Turns the xenon lamp on or off.</div>	✓	✓	✓
	[Hg Lamp]	6.4.5 [Light Source...] <div>Note: Turn off the Hg lamp.</div>	✓	✓	✓
	[Scales]	6.5.1 [Scales...]	✓	✓	
	[Patterns]	6.5.2 [Patterns...]	✓	✓	✓
	[Fonts]	6.5.3 [Fonts ...]	✓	✓	✓
	[Gridlines]	6.5.4 [Gridlines...]	✓	✓	✓

	[Styles]	6.5.5 [Styles...]	✓	✓	✓
	[Digital]	6.5.12 [View Mode]			✓
	[Graph]	6.5.12 [View Mode]			✓

Window

Title bar	Displays the program name.
Menu bar	Displays the menus for operation.
Monitor bar	Displays from left to right the current wavelength, number of cycles or number of accumulations and photometric values.
Toolbar	Displays available tool buttons. These tool buttons can be customized in the [Settings] menu.
Information bar	Displays the current measurement parameters.
Main unit bar	Displays current detector information and alarm information for the J-1000 CD spectrometer.
Accessory bar	Displays information about the attached accessory.
Spectrum display field	Displays the spectrum or time-course data being measured.
Status bar	Displays the current state of the instrument and an explanation of the selected menu.

Menu

[File] menu		Spectra Measurement	Time Course Measurement	Data Monitor
[Save Data...]	Saves the spectrum with a new file name. <i>Note: In the CFR version, this function is not available.</i>	✓	✓	
[Send to Analysis]	Transfers the measured data to the [Spectra Analysis] program.	✓	✓	
[Open Parameters...]	Selects a parameter file and opens those parameters.	✓	✓	
[Save Parameters...]	Saves the currently specified measurement parameters with the entered file name. <i>Note: In the CFR version, this function is not available.</i>	✓	✓	
[Open Baseline...]	Selects a baseline file and opens that baseline spectrum. <i>Note: In the CFR version, this function is not available.</i>	✓		

[Save Baseline...]	Saves the currently measured baseline spectrum with the entered file name. <i>Note: In the CFR version, this function is not available.</i>	✓		
[Exit]	Exits the measurement program and returns to [Spectra Manager]. <i>Note: This command is in the [Measure] menu in the [Data Monitor] program.</i>	✓	✓	✓
[Measure] menu				
[Cancel]	Cancels the measurement.	✓	✓	
[Sequence Measurement]	Executes the sequence.	✓		
[Sample Measurement]	Measures the spectrum or time-course data for the sample.	✓	✓	
[Baseline Measurement]	Measures the baseline spectrum for baseline correction.	✓		
[Blank Measurement]	Measures the blank for blank correction.		✓	
[Dark Measurement]	Measures the dark spectrum for dark correction.	✓	✓	
[Parameters...]	Specifies parameters and saves or opens the measurement parameters.	✓	✓	✓
[Control] menu				
[Zero Clear]	Sets the photometric value of ORD-M or fluorescence intensity at the current wavelength to zero.	✓	✓	✓
[Shutter...]	Opens/closes the shutter.	✓	✓	✓
[Move Wavelength...]	Moves the CD spectrometer wavelength to the desired wavelength.	✓	✓	✓
[Detector...]	Sets the detector sensitivity.	✓	✓	✓
[Light Source]	Turns on or off the xenon lamp.	✓	✓	✓
[Nitrogen Flowrate...]	Specifies the flow rate of nitrogen gas. This command is only displayed when [Automatic (PC control)] is selected for the nitrogen flow meter.	✓	✓	✓
[Select Detector...]	Selects a detector.	✓	✓	✓
[Select Accessory...]	Selects an accessory.	✓	✓	✓


[View] menu				
[Scales...]	Specifies the display scale for the spectrum or time-course data.	✓	✓	
[Patterns...]	Selects the display color and line style for the spectrum or time-course data.	✓	✓	
[Fonts...]	Specifies the display font for the spectrum or time-course data.	✓	✓	
[Gridlines...]	Shows/hides the gridlines for the spectrum or time-course data.	✓	✓	
[Styles...]	Specifies the display style for the spectrum or time-course data.	✓	✓	
[Overlay]	Overlays spectra or time-course data when cycle measurement is selected.	✓	✓	
[Select Graph...]	Selects the graph to be displayed on the CD spectrum display field.	✓	✓	
[Information Bar]	Shows/hides the information bar.	✓	✓	✓
[Accessory Bar]	Shows/hides the accessory bar.	✓	✓	✓
[Toolbars]	Shows/hides toolbars.	✓	✓	✓
[Status Bar]	Shows/hides the status bar.	✓	✓	✓
[View Mode]	Displays the photometric value in digital or graph mode.			✓
[Graph View]	Selects the graph to be displayed when graph mode is selected for the photometric value display.			✓
[Settings] menu				
[Default Parameters...]	Specifies measurement parameters for running the [Spectra Measurement] and [Time Course Measurement] programs based on the currently recognized accessory. <div style="border: 1px solid black; padding: 5px; margin-top: 10px;"><i>Note: In the CFR version, this function is not available.</i></div>	✓	✓	
[Customize Toolbar...]	Customizes and shows/hides toolbars.	✓	✓	✓
[Monitor Bar...]	Customizes the style of the monitor bar.	✓	✓	✓
[Help] menu				
[About...]	Displays version information for the program.	✓	✓	✓

6.2 [File] Menu

Saves or opens a measured data or parameter file.

6.2.1 [Save Data...]

Saves the measured data with a new file name.

Note 1: This operation can also be performed by clicking the tool button  [Save].

Note 2: In the CFR version, since all measurement data are always saved, this function is not available.

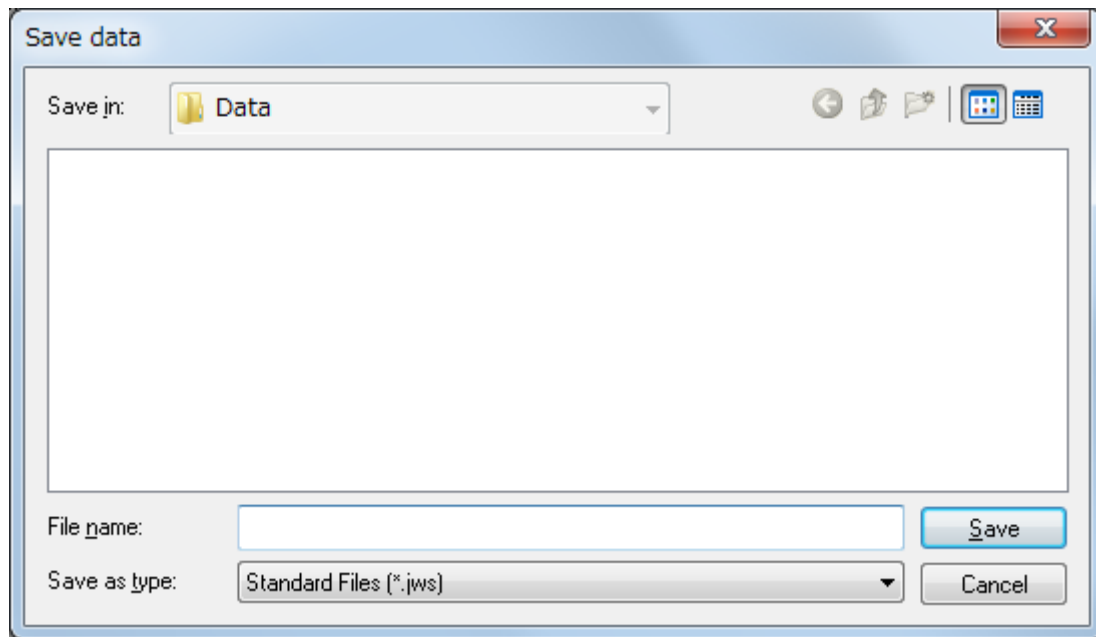


Figure 6.2 [Save data] dialog box

[Save in] list	Selects a drive or folder to browse using the drop-down list. Files saved in the currently opened folder are displayed in the file name list.
File name list	Lists the files in the currently opened folder. Use as a reference when naming a file.
[File name]	Enters a file name for saving the measured data. An extension may be omitted.
[Save as type]	Selects the files (type of extension) to display in the file name list. Files other than Standard Files (*.jws) cannot be selected.
<Save>	Saves the measured data and closes the dialog box.
<Cancel>	Closes the dialog box without saving the measured data.

6.2.2 [Send to Analysis]

Transfers the measured data to the [Spectra Analysis] program. Selecting the [To Spectra Analysis] check box in the [Data] tab of the [Parameters] dialog box from the [Measure] menu automatically transfers the measured spectrum to the [Spectra Analysis] program after measurement.

When a value of 2 or more is specified for the number of cycles in the [Parameters] dialog box, the dialog box in Figure 6.3 is displayed.

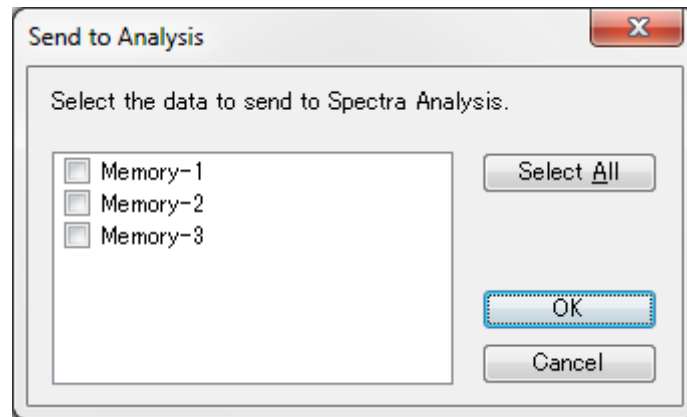




Figure 6.3 [Send to Analysis] dialog box

Select the check boxes for the data to send to the [Spectra Analysis] program (or click the <Select All> button to send all data), and then click the <OK> button.

Note: This operation can also be performed by clicking the tool button  [Send to Analysis].

6.2.3 [Open Parameters...]

Selects a parameter file and opens those parameters.

Note 1: This operation can also be performed by clicking the tool button  [Open Parameters].

Note 2: The dialog box in Figure 6.4 displays only the parameter files for the currently attached accessory.

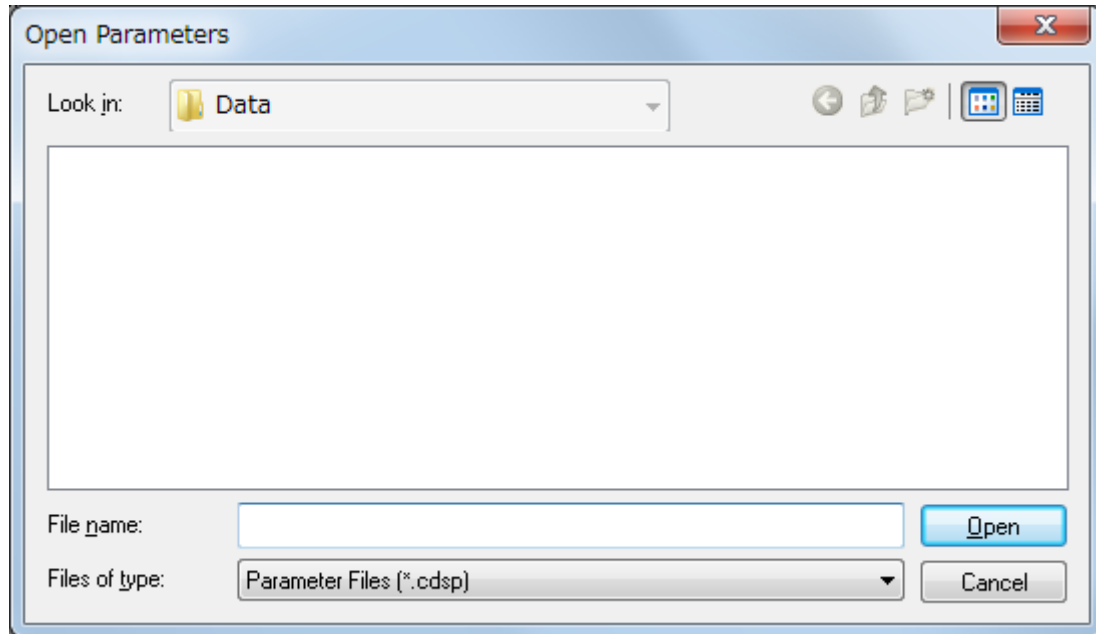



Figure 6.4 [Open Parameters] dialog box

[Look in] list	Selects a drive and folder to browse using the drop-down list. Files saved in the currently opened folder are displayed in the file name list.
File name list	Lists the files in the currently opened folder. Selects the file name of the parameters to open.
[File name]	Enters the file name for the parameters to open. An extension may be omitted since the appropriate extension for the measurement program is automatically appended.
[Files of type]	Selects the files (type of extension) to display in the file name list. Files other than Parameter Files (having the appropriate extension for the measurement program) cannot be selected.
<Open>	Opens a parameter file and closes the dialog box.
<Cancel>	Closes the dialog box without opening a parameter file.

Note: In the CFR version, the [Look in] drop-down list cannot be used to select a folder. However, a subfolder within the prespecified folder can be selected.

6.2.4 [Save Parameters...]

Saves the currently specified measurement parameters with the entered file name.

Note1: This operation can also be performed by clicking the tool button  [Save Parameters].

Note2: In the CFR version, since all measurement parameters are always saved, this function is not available.

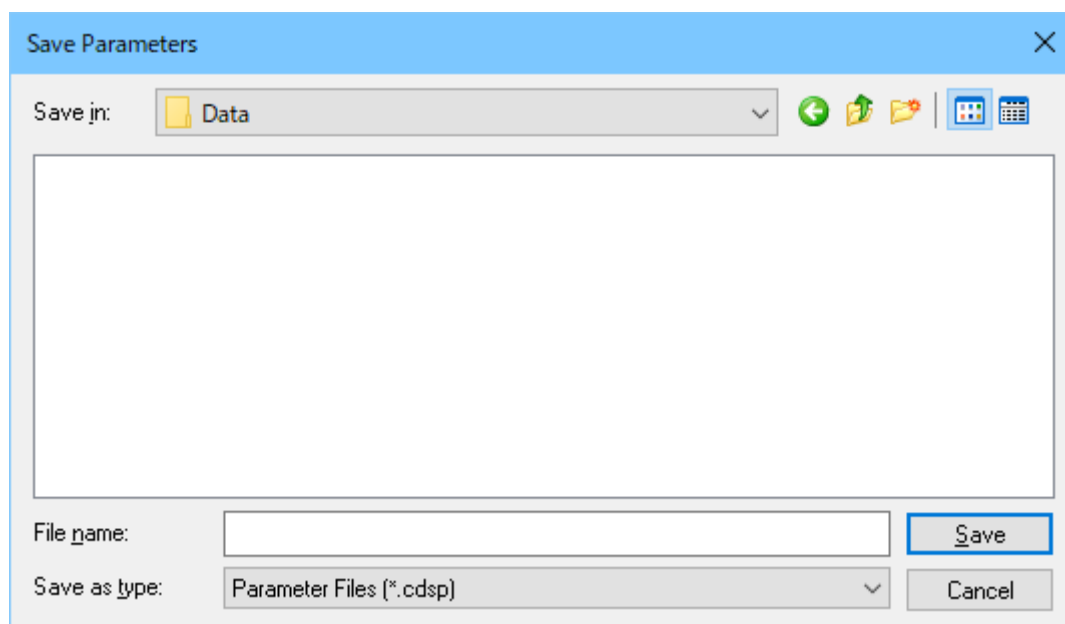


Figure 6.5 [Save Parameters] dialog box

[Save in] list	Selects a drive or folder to browse using the drop-down list. Files saved in the currently opened folder are displayed in the file name list.
File name list	Lists the files in the currently opened folder. Use as a reference when naming a file.
[File name]	Enters a file name for saving parameters. An extension may be omitted since the appropriate extension for the measurement program is automatically appended.
[Save as type]	Selects the files (type of extension) to display in the file name list. Files other than Parameter Files (having the appropriate extension for the measurement program) cannot be selected.
<Save>	Saves the parameter file and closes the dialog box.
<Cancel>	Closes the dialog box without saving the parameter file.

6.2.5 [Open Baseline...]

Selects a baseline file and opens that baseline. When a baseline file is opened, the sample spectrum is corrected by using the opened baseline spectrum.

Note 1: In the CFR version, this function is not available.

Note 2: The following dialog box displays only the baseline files for the currently attached accessory.

Note 3: This function is available if [Baseline] is selected in the [Control] tab of the [Parameters] dialog box.

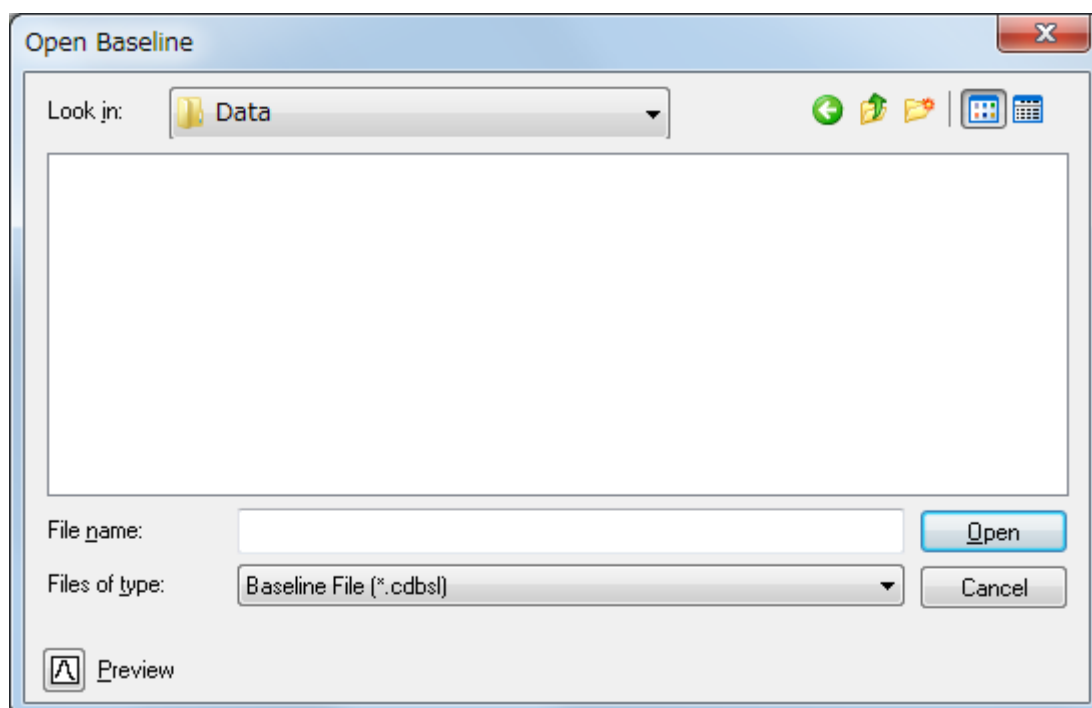



Figure 6.6 [Open Baseline] dialog box

Click the  <Preview> button in Figure 6.6 to display the spectrum of the selected baseline file (see Figure 6.7).

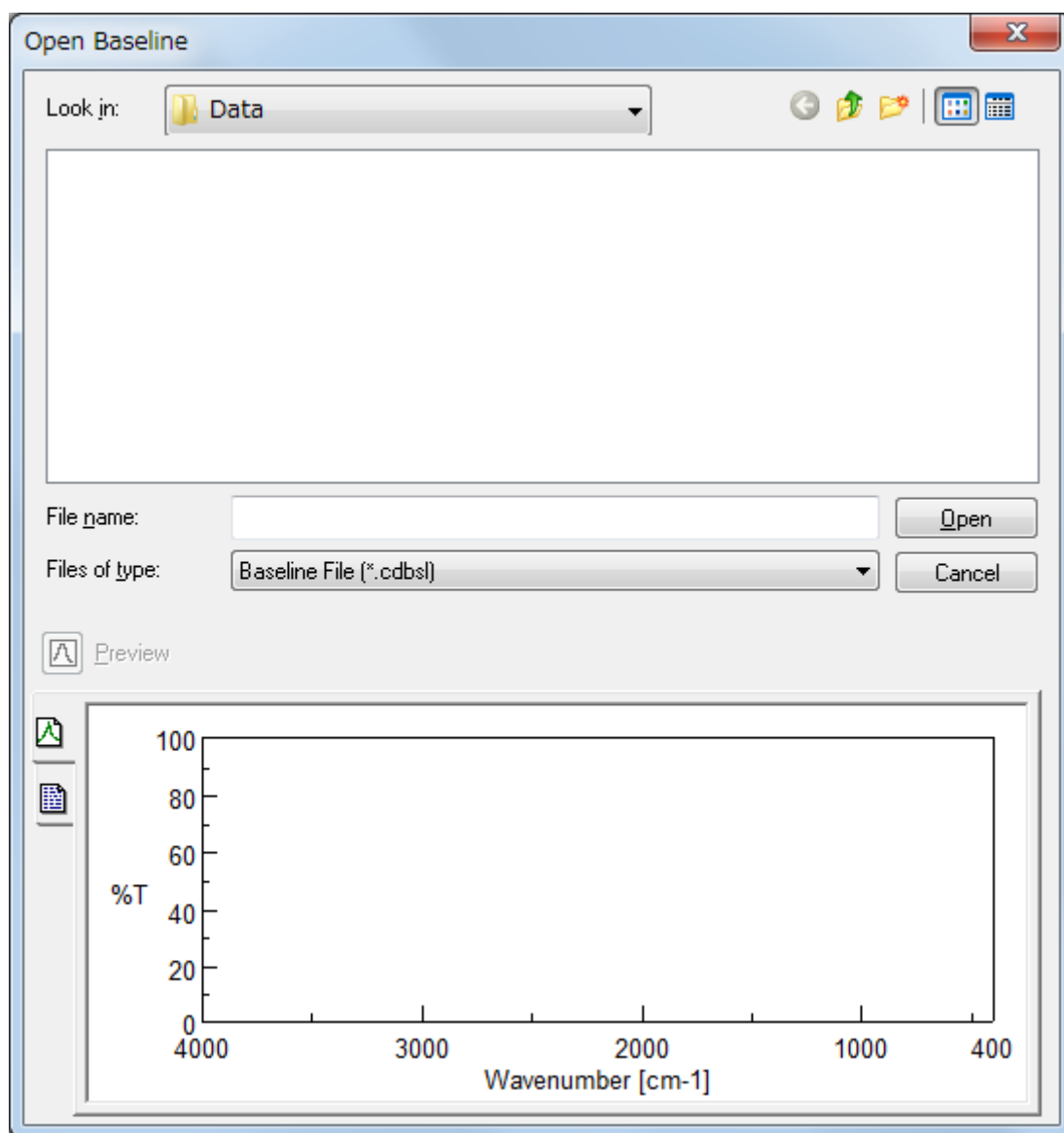


Figure 6.7 [Open Baseline] dialog box

[Look in] list	Selects a drive or folder to browse using the drop-down list. Files saved in the currently opened folder are displayed in the file name list.
File name list	Lists the files in the currently opened folder. Selects the file name of the baseline spectrum to open.
[File name]	Enters the file name of the baseline spectrum to open. An extension may be omitted since the extension ".cdbsl" is automatically appended.
[Files of type]	Selects the files (type of extension) to display in the file name list. Files other than Baseline Files (*.cdbsl) cannot be selected.
<Preview>	Displays the baseline spectrum and information about the selected file (see Figure 6.7).
<Open>	Opens the baseline file and closes the dialog box.
<Cancel>	Closes the dialog box without opening a baseline file.

6.2.6 [Save Baseline...]

Saves the currently measured baseline spectrum with the entered file name.

Note: In the CFR version, this function is not available.

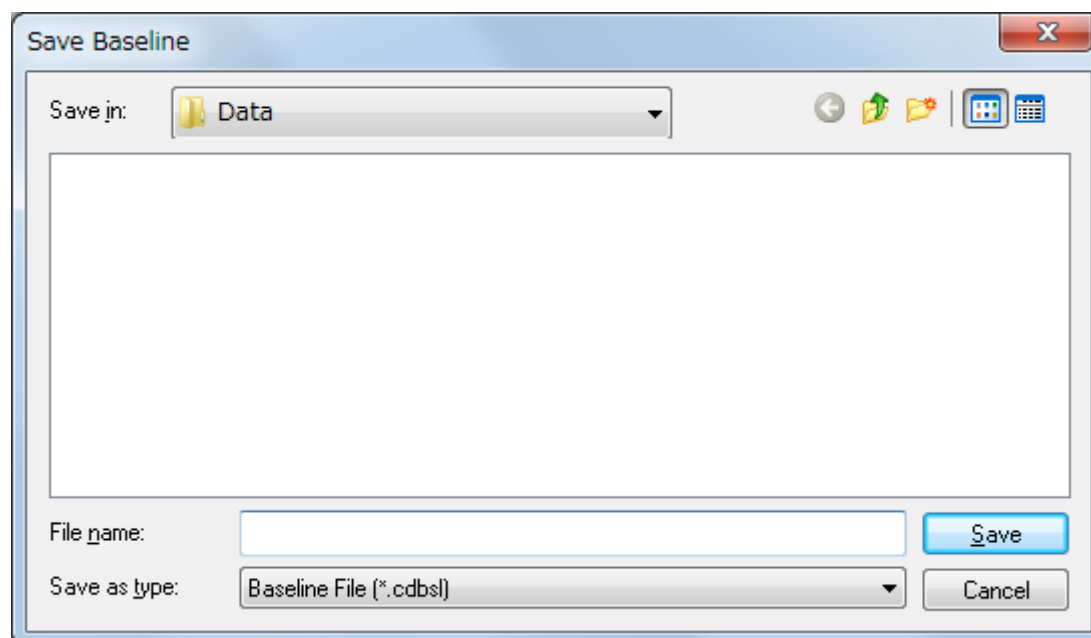


Figure 6.8 [Save Baseline] dialog box

[Save in] list	Selects a drive or folder to browse using the drop-down list. Files saved in the currently opened folder are displayed in the file name list.
File name list	Lists the files in the currently opened folder. Use as a reference when naming a file.
[File name]	Enters a file name for saving the baseline spectrum. An extension may be omitted since the extension ".cdbsl" is automatically appended.
[Save as type]	Selects the files (type of extension) to display in the file name list. Files other than Baseline Files (*.cdbsl) cannot be selected.
<Save>	Saves the measured baseline spectrum and closes the dialog box.
<Cancel>	Closes the dialog box without saving the measured baseline spectrum.

6.2.7 [Print Setup...]

Specifies the target printer and print parameters. The content of this dialog box varies depending on the connected printer. Configure the printer settings if necessary.

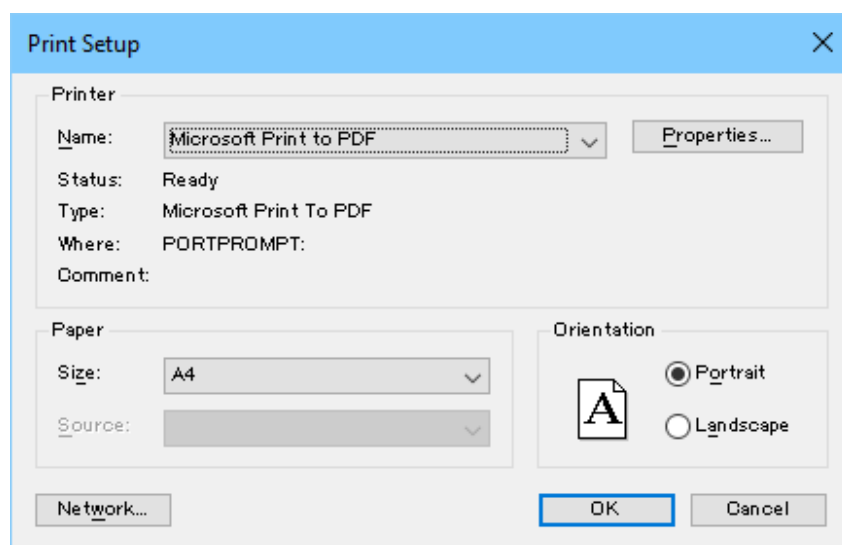


Figure 6.9 [Print Setup] dialog box

Printer	Displays the available printers. To add a new printer, select [Control Panel] - [Printers] from the <i>Windows</i> task bar and select [Add Printer...].
Paper	Selects the paper size and paper source tray.
Orientation	Selects the orientation of the paper for printing.
<OK>	Applies the changes to the [Print Setup] settings and closes the dialog box.
<Cancel>	Closes the dialog box without applying any changes to the [Print Setup] settings.

6.2.8 [Exit]

Exits the [Spectra Measurement] program.

6.3 [Measure] Menu

Starts or cancels the measurement, and specifies the measurement parameters.

6.3.1 [Cancel]

Cancels the measurement. A dialog box to confirm whether to keep the measurement data is displayed.

Note: This operation can also be performed by clicking the tool button  [Cancel].

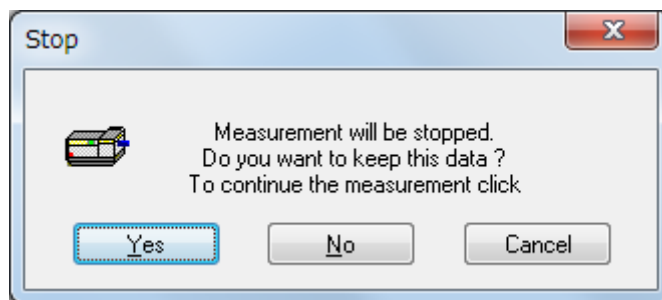



Figure 6.10 [Cancel] dialog box

<Yes>	Keeps the measured spectrum.
<No>	Discards the measured spectrum.
<Cancel>	Continues the measurement.


6.3.2 [Sequence Measurement]

Executes the sequence measurement. This function is enabled when autosampler is connected.

Note: This operation can also be performed by clicking the tool button  [Sequence Measurement].

6.3.3 [Sample Measurement]

Measures the CD spectrum of the sample. Selecting the [To Spectra Analysis] check box in the [Data] tab of the [Parameters] dialog box from the [Measure] menu automatically transfers the measured data to the [Spectra Analysis] program after measurement. The measured data is displayed in a new View.


Note: This operation can also be performed by clicking the tool button  [Sample Measurement].

6.3.4 [Baseline Measurement]

Measures the baseline spectrum for baseline correction. Available if [Baseline] is selected in the [Control] tab of the [Parameters] dialog box.

Baseline correction removes the effects of solvents and additional agents other than the target sample. The true sample spectrum is obtained by subtracting the baseline spectrum from the measured spectrum.

Confirm that the sample for baseline measurement has been inserted, and then click the <Measure> button.

Note: This operation can also be performed by clicking the tool button  [Baseline Measurement].

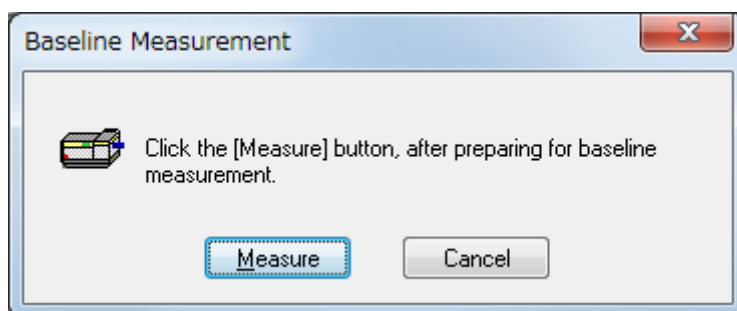


Figure 6.11 [Baseline Measurement] dialog box


<Measure>	Measures the baseline spectrum for baseline correction.
<Cancel>	Closes the [Baseline Measurement] dialog box without measuring the baseline

6.3.5 [Blank Measurement]

Measures the blank data for blank correction. Available if [Blank] is selected in the [Control] tab of the [Parameters] dialog box.

Blank correction removes the effects of solvents and additional agents other than the target sample. The sample signal is obtained by subtracting the blank data from the data obtained during measurement of the sample.

Confirm that the blank has been inserted, and then click the <Measure> button.

Note: This operation can also be performed by clicking the tool button  [Blank Measurement].

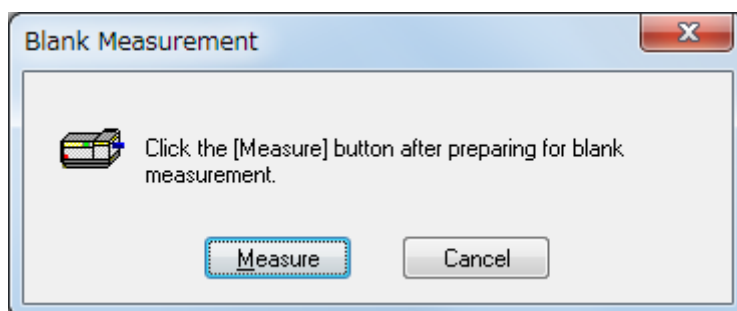



Figure 6.12 [Blank Measurement] dialog box

<Measure>	Measures the blank spectrum for blank correction.
<Cancel>	Closes the [Blank Measurement] dialog box without measuring the blank.

6.3.6 [Dark Measurement]

Measures the photometric value for dark correction. Available if the InGaAs or APD detector is used.

Note: This operation can also be performed by clicking the tool button  [Dark Measurement].

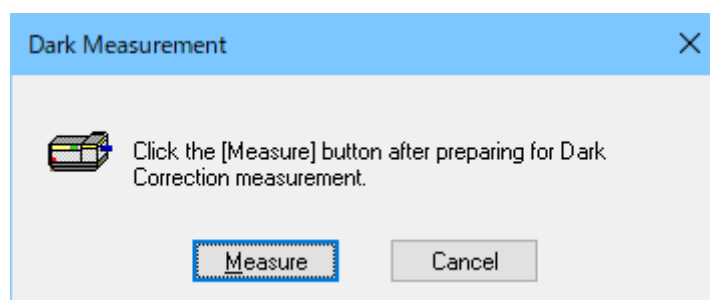


Figure 6.13 [Dark Measurement] dialog box

<Measure>	Measures the photometric value for dark correction.
<Cancel>	Closes the [Dark Measurement] dialog box without measuring the dark current.


6.3.7 [Parameters...]

Specifies the measurement parameters. Refer to each measurement program reference.

6.4 [Control] menu


6.4.1 [Zero Clear]

Sets the photometric value of the ORD-M or fluorescence intensity to zero. Available if an ORD-M or fluorescence detector is connected.

Note: This operation can also be performed by clicking the tool button  [Zero Clear].

6.4.2 [Shutter]

Opens/closes the shutter.

Note: This operation can also be performed by clicking the tool button  [Shutter].

6.4.3 [Move Wavelength...]

Moves the wavelength of the CD spectrometer to the desired wavelength. When the measurement is started, data acquisition begins once the wavelength returns to the value specified in [Parameters].

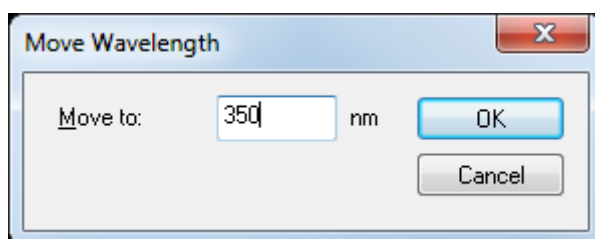



Figure 6.14 [Move Wavelength] dialog box

[Move to]	Specifies the wavelength. J-1100: 180 to 600 nm J-1500: 163 to 1600 nm J-1700: 163 to 2500 nm
<OK>	Moves the wavelength of the CD spectrometer to the specified wavelength and closes the dialog box.
<Cancel>	Closes the dialog box without applying any changes to the [Move Wavelength] settings.

Note: This operation can also be performed by clicking the tool button  [Move Wavelength].

6.4.4 [Detector ...]

Specifies the sensitivity for monitoring the photometric value.

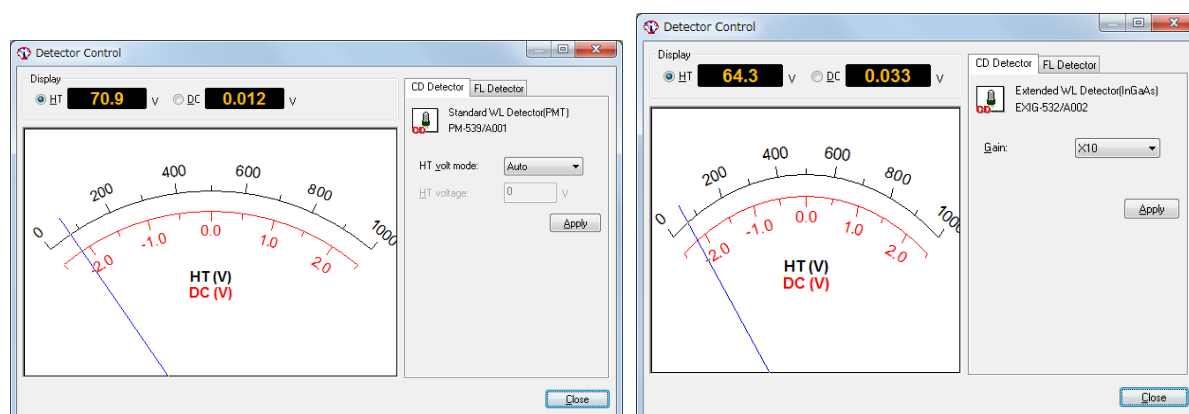


Figure 6.15 [Detector Control] dialog box (left: PMT, APD) (right: InGaAs, InSb)

Display	
[HT]	Displays the HT voltage for the detector.
[DC]	Displays the DC voltage for the detector.
[HT volt mode]	
[Auto]	Automatically sets the applied voltage for the photomultiplier tube.
[Manual]	Manually sets the applied voltage for the photomultiplier tube.
[Off]	Sets the detector voltage off.
[HT voltage]	Specifies the photomultiplier tube voltage when [Manual] is selected for [HT volt mode]. Setting range: 0 to 1000 V
[Gain]	Selects the voltage gain for the InGaAs or InSb detector. Options: x1, x10, x100
<Apply>	Sets the sensitivity of the CD spectrometer to the specified sensitivity.
<Cancel>	Closes the dialog box without applying any changes to the [Detector Control] settings.

Note: This operation can also be performed by clicking the tool button



6.4.5 [Light Source...]

Turn on/off the light source.

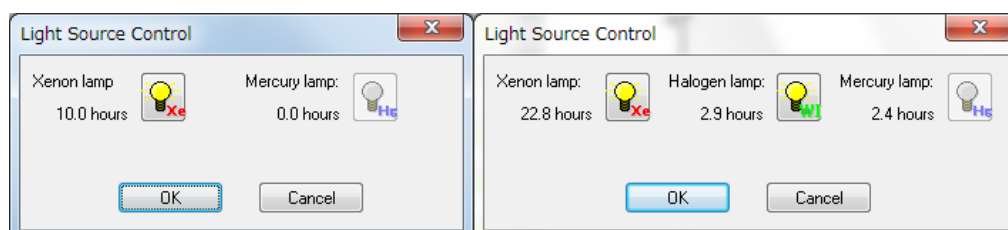






Figure 6.16 [Light Source Control] dialog box J-1100/1500 (left) J-1700 (right)

Note 1: This operation can also be performed by clicking the tool button    [Light Source].

Note 2: The tool button  [Mercury lamp] is available to turn off only.

6.4.6 [Nitrogen Flowrate...]

Controls the nitrogen gas flow rate. Available if a PC-controlled nitrogen-gas auto flow meter is selected.

Note: Refer to the software manual “J-1000 series Nitrogen Flowmeter” for detail.

6.4.7 [Select Detector...]

Recognizes a detector manually. Select the detector to use and click the <OK> button. A dialog box stating “Connecting to the detector...” is displayed.

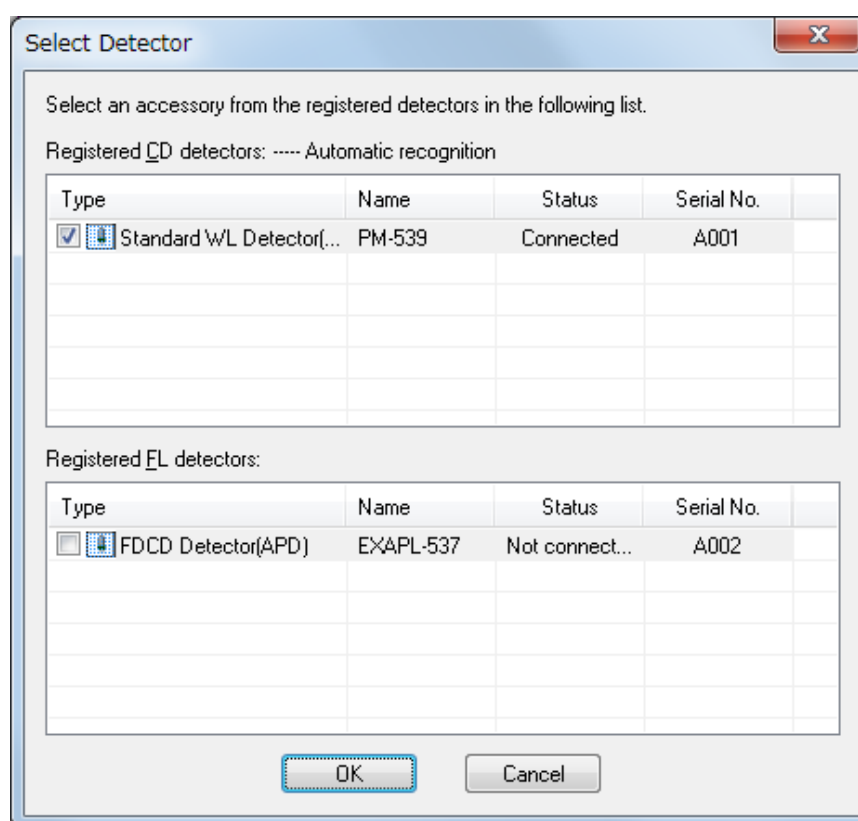


Figure 6.17 [Select Detector] dialog box

6.4.8 [Select Accessory...]

Selects an accessory (cell unit, external accessories or measurement units) to be used. After clicking the <OK> button, the message "The accessory was attached successfully." is displayed.

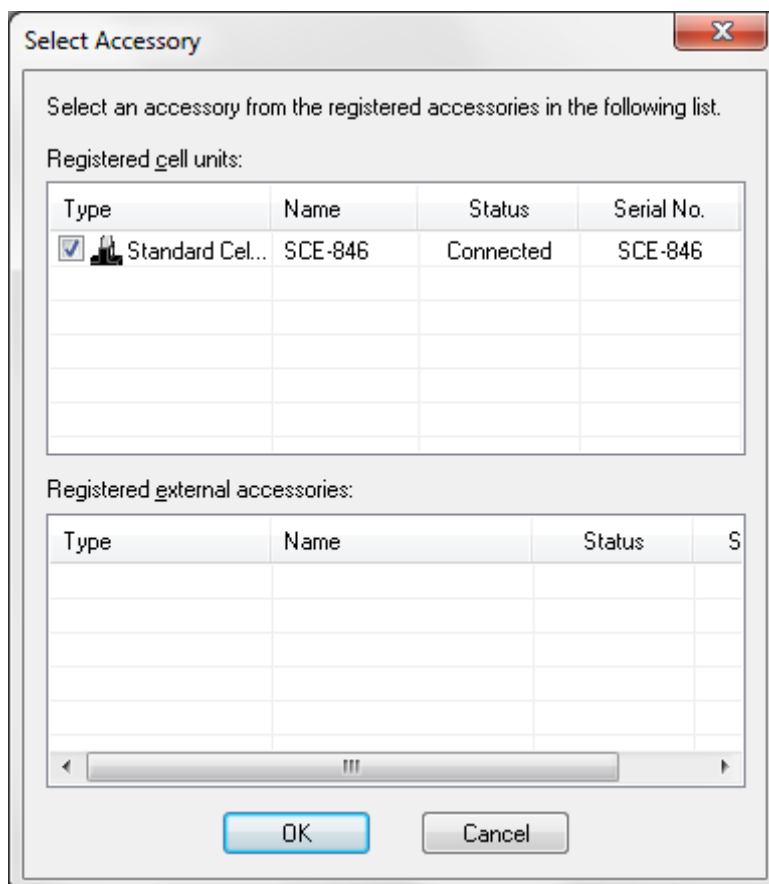


Figure 6.18 [Select Accessory] dialog box

Note: In the CFR version, the function for automatically starting registered programs for accessories is not provided.

6.4.9 [Accessory Model Number]

When an accessory that has been registered is attached, the accessory name is displayed as a command on the [Control] menu (see Figure 6.19). Select the command to control the attached accessory.

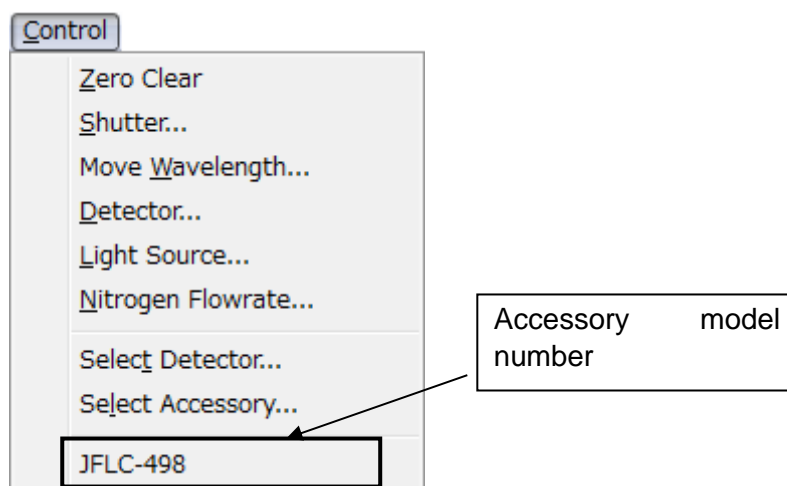


Figure 6.19 [Control] dialog box

Note 1: Use this function to return the temperature of the Peltier device to room temperature after measuring data at high or low temperatures to set a target temperature after inserting a sample.

Note 2: This operation can also be performed by clicking the <Setting...> button in the accessory bar.

6.5 [View] Menu

Specifies the view options.

6.5.1 [Scales...]

Specifies the display scale for the spectrum.

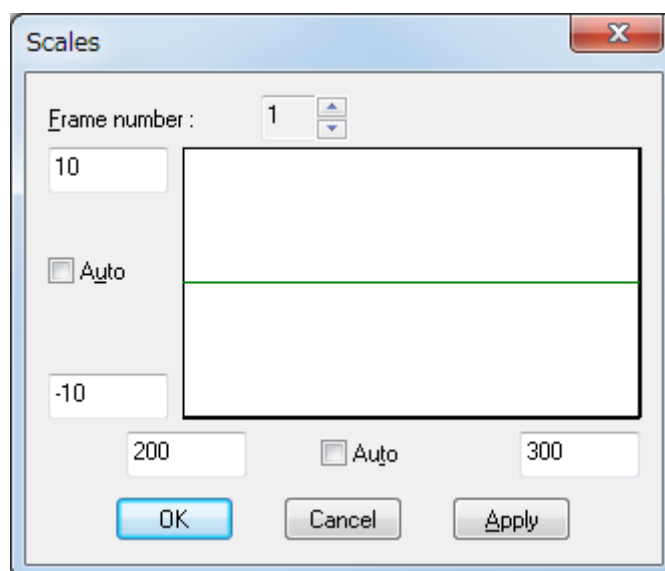


Figure 6.20 [Scales] dialog box

[Frame number]	Selects the graph for which to change the scales.
Horizontal axis	Enters the scale for the horizontal axis. Selecting the [Auto] check box displays the entire range irrespective of the entered values.
Vertical axis	Enters the scale for the vertical axis. Selecting the [Auto] check box displays the graph at the optimal scale for the specified horizontal axis range.
<OK>	Applies the changes to the [Scales] settings and closes the dialog box.
<Cancel>	Closes the dialog box without applying any changes to the [Scales] settings.
<Apply>	Applies the specified scales to the graph selected in [Frame number].

Note: This operation can also be performed by clicking the tool button  [Scales].

6.5.2 [Patterns...]

Selects the display color and line style for the spectrum.

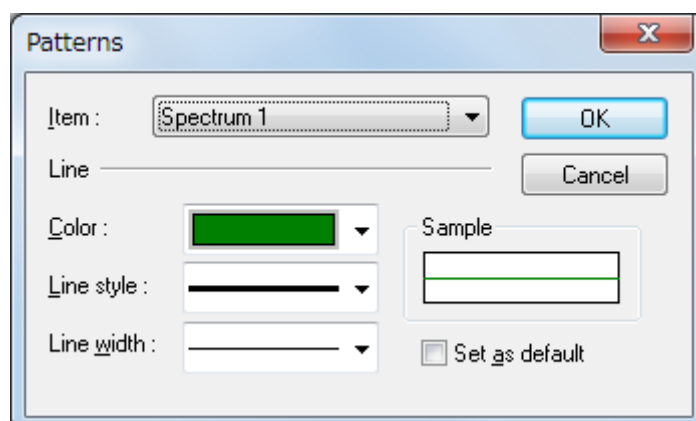



Figure 6.21 [Patterns] dialog box

[Item]	Selects an item from spectra 1 to 16, frame, gridlines, and auxiliary gridlines to change its color and line style.
[Color]	Shows the available colors for the selected item in the drop-down list.
[Line style]	Shows the available line styles. The currently selected item in the [Item] drop-down list is displayed with the selected line style.
[Line width]	Shows the available line widths. The currently selected item in the [Item] drop-down list is displayed with the selected line width.
Sample	Displays a sample of the selected pattern.
[Set as default]	Applies the selected patterns to all subsequently opened Views.
<OK>	Applies the changes to the [Patterns] settings and closes the dialog box.
<Cancel>	Closes the dialog box without applying any changes to the [Patterns] settings.

Note: This operation can also be performed by clicking the tool button  [Patterns].

6.5.3 [Fonts...]

Specifies the display font for the data.

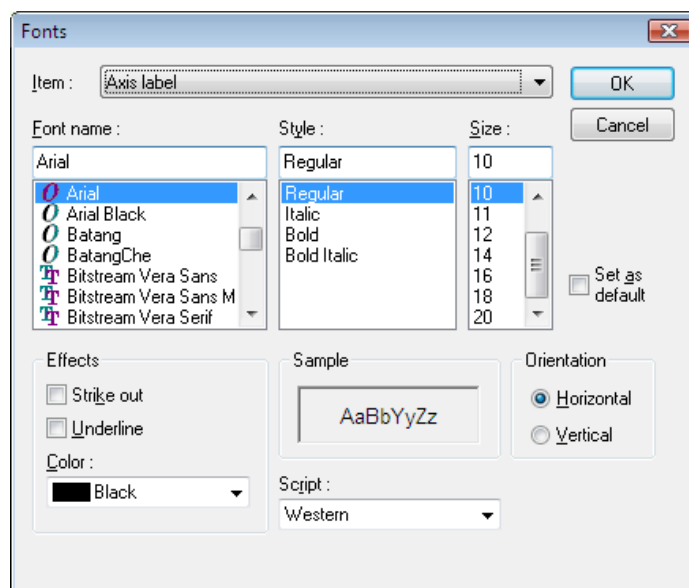



Figure 6.22 [Fonts] dialog box

[Item]	<p>Selects an item to change its font settings.</p> <p>[Axis label]: Character ([CD (mdeg)], [nm], or other)</p> <p>[Scale label]: Numeric value</p>
[Font name]	Lists available font names.
[Style]	Lists available styles.
[Size]	Lists available sizes.
Effects	<p>Selects strikeout and underline effects.</p> <p>[Strike out]: Draws a line through the middle of the text.</p> <p>[Underline]: Underlines the text.</p>
Sample	Displays a sample of the specified font.
Orientation	<p>Selects the orientation for displaying the axis label.</p> <p>[Horizontal]: Displays the y-axis label horizontally</p> <p>[Vertical]: Displays the y-axis label vertically</p>
[Set as default]	Applies the specified fonts to all subsequently opened Views.
[Color]	Shows the available colors for the selected item in the drop-down list.
[Script]	Selects the language for the specified font.
<OK>	Applies the changes to the [Fonts] settings and closes the dialog box.
<Cancel>	Closes the dialog box without applying any changes to the [Fonts] settings.

Note: This operation can also be performed by clicking the tool button  [Fonts].

6.5.4 [Gridlines...]

Shows/hides the gridlines for the data.

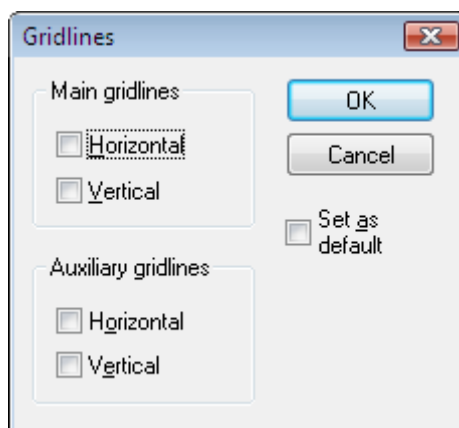



Figure 6.23 [Gridlines] dialog box

Main gridlines	
[Horizontal]	Displays major horizontal axis gridlines.
[Vertical]	Displays major vertical axis gridlines.
Auxiliary gridlines	
[Horizontal]	Displays minor horizontal axis gridlines.
[Vertical]	Displays minor vertical axis gridlines.
[Set as default]	Applies the selected gridlines to all subsequently opened Views.
<OK>	Applies the changes to the [Gridlines] settings and closes the dialog box.
<Cancel>	Closes the dialog box without applying any changes to the [Gridlines] settings.

Note: This operation can also be performed by clicking the tool button  [Gridlines].

6.5.5 [Styles...]

Specifies the display style for the spectrum.

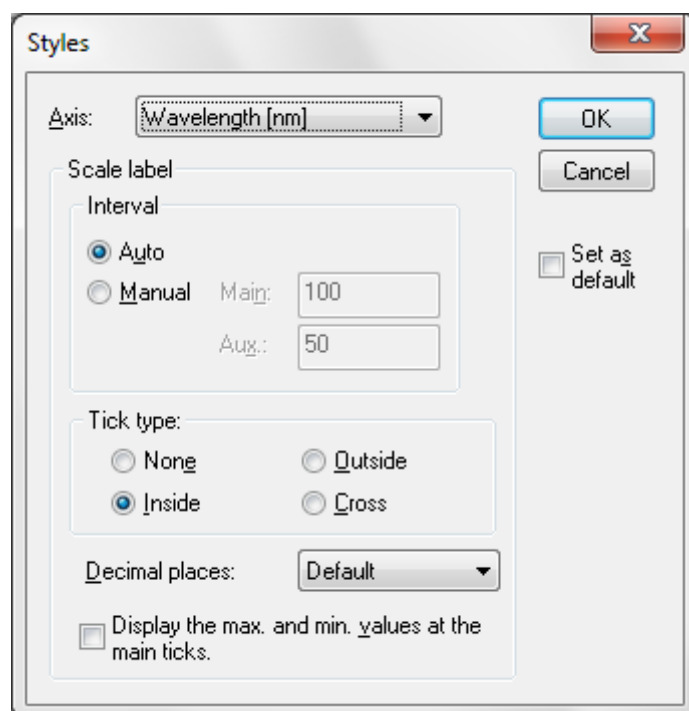



Figure 6.24 [Styles] dialog box

[Axis]	Selects an item to change its display style.
Interval	
[Auto]	Automatically selects the gridline format.
[Manual]	Manually specifies the gridline format for the main and auxiliary scales.
Tick type	
[None]	Displays no ticks on the axis
[Outside]	Displays ticks on the outside of the axis.
[Inside]	Displays ticks on the inside of the axis.
[Cross]	Displays ticks on both sides of the axis.
[Decimal places]	Selects the number of decimal places for the display values on the horizontal and vertical axes.
[Display the max. and min. values at the main ticks]	Displays the maximum and minimum values at the main ticks. Values at the auxiliary ticks are not displayed.
[Set as default]	Applies the specified styles to all subsequently opened Views.
<OK>	Applies the changes to the [Styles] settings and closes the dialog box.
<Cancel>	Closes the dialog box without applying any changes to the [Styles] settings.

Note: This operation can also be performed by clicking the tool button  [Styles].

6.5.6 [Overlay]

Overlays the spectra in the spectrum measurement display field when cycle measurement is selected.

6.5.7 [Select Graph...]

Selects the graphs to display in the data display field.

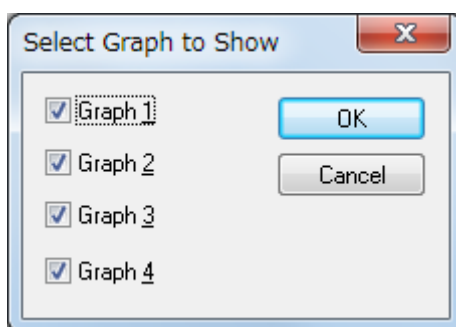


Figure 6.25 [Select Graph to Show] dialog box

[Graph 1]	Displays the graph for photometric mode 1 as graph 1 in the data display field.
[Graph 2]	Displays the graph for photometric mode 2 as graph 2 in the data display field.
[Graph 3]	Displays the graph for photometric mode 3 as graph 3 in the data display field.
[Graph 4]	Displays the graph for photometric mode 4 as graph 4 in the data display field.
<OK>	Selects the graphs to display and closes the dialog box.
<Cancel>	Closes the dialog box without selecting a graphs to display.

6.5.8 [Information Bar]

Shows/hides the information bar.

6.5.9 [Accessory Bar]

Shows/hides the accessory bar.

6.5.10 [Toolbars]

Shows/hides toolbars.

[File]	Shows/hides the toolbar corresponding to the [File] menu.
[Measure]	Shows/hides the toolbar corresponding to the [Measure] menu.
[Control]	Shows/hides the toolbar corresponding to the [Control] menu.
[View]	Shows/hides the toolbar corresponding to the [View] menu.

6.5.11 [Status Bar]

Shows/hides the status bar.

6.5.12 [View Mode]

Displays the photometric value in digital or graph mode in the photometric value monitor.

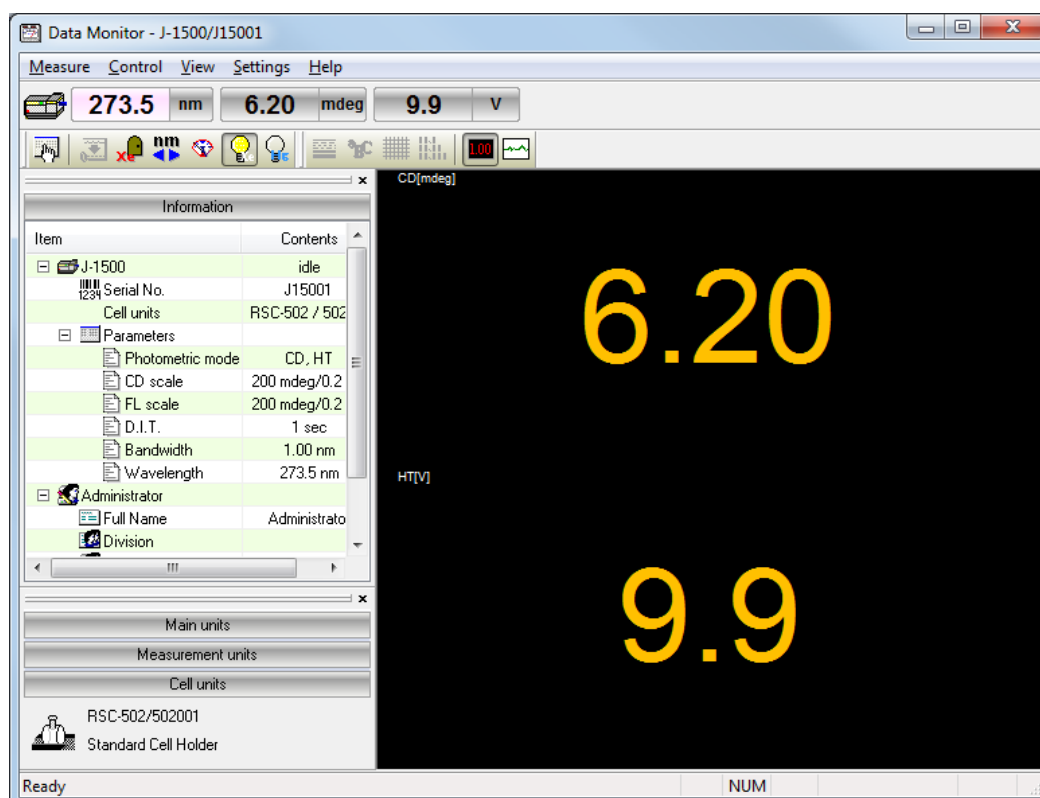


Figure 6.26 Digital mode of the photometric value monitor

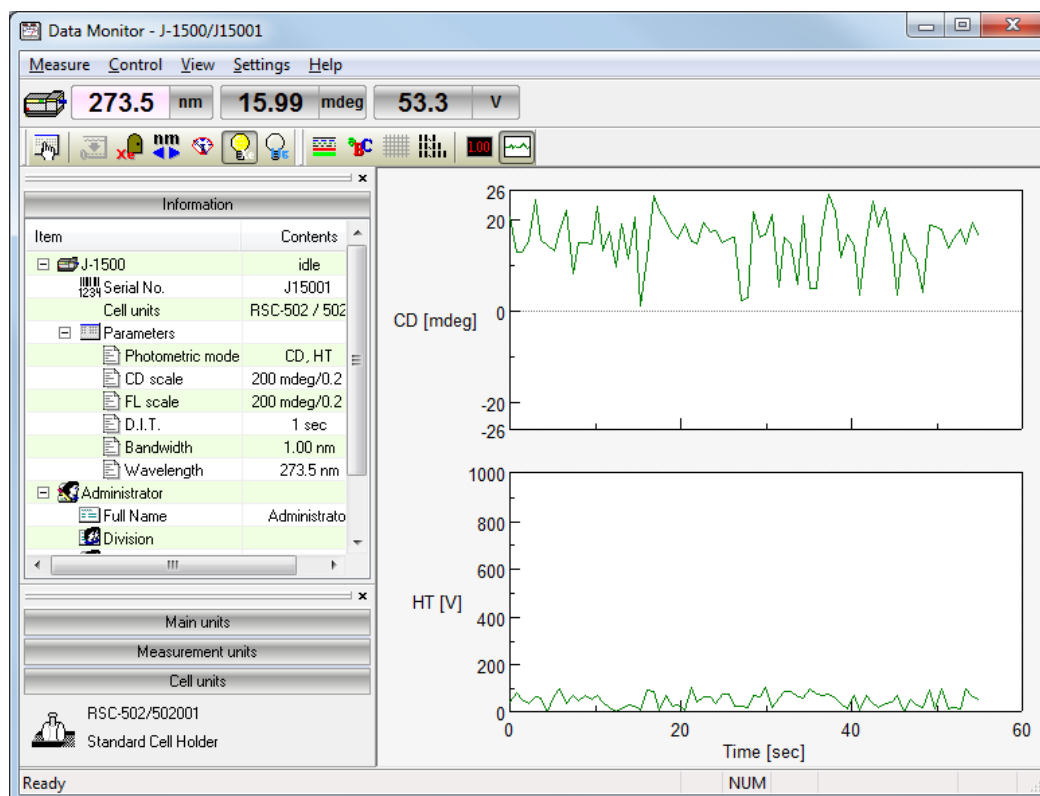


Figure 6.27 Graph mode of the photometric value monitor

6.5.13 [Graph View]

Specifies [Patterns], [Fonts], [Gridlines] and [Styles] when graph mode is selected for the photometric value monitor. Refer to Sections 6.5.2 [Patterns...], 6.5.3 [Fonts...], 6.5.4 [Gridlines...] and 6.5.5 [Styles...] for details about setting each command.

6.6 [Settings] Menu

6.6.1 [Default Parameters...]

Note: In the CFR version, this function is not available.

Specifies measurement parameters for running the measurement program based on the currently recognized accessory. Default parameters can be specified for each accessory. Attach the accessory to be used and specify its default parameters.

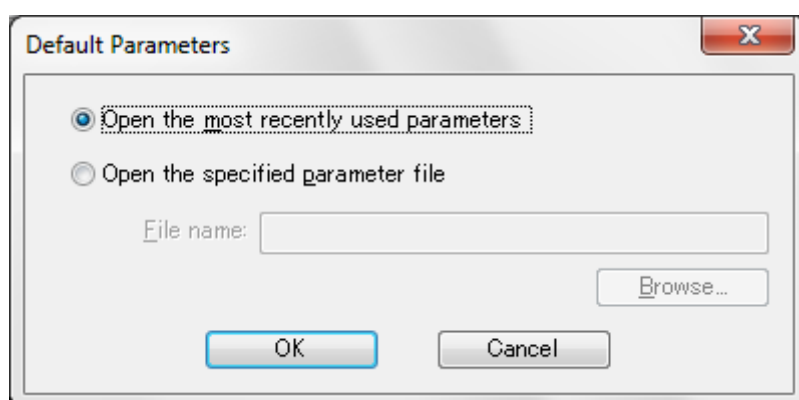


Figure 6.28 [Default Parameters] dialog box

[Open the most recently used parameters]	Starts with the parameters previously specified before exiting the measurement program.
[Open the specified parameter file]	Starts with a specified parameter file.
[File name]	Displays the name of path to the parameter file when the [Open the specified parameter file] option button is selected.
<Browse...>	Selects the drive and file name for a saved parameter file (see Figure 6.29).
<OK>	Applies the changes to the [Default Parameters] settings and closes the dialog box.
<Cancel>	Closes the dialog box without applying any changes to the [Default Parameters] settings.

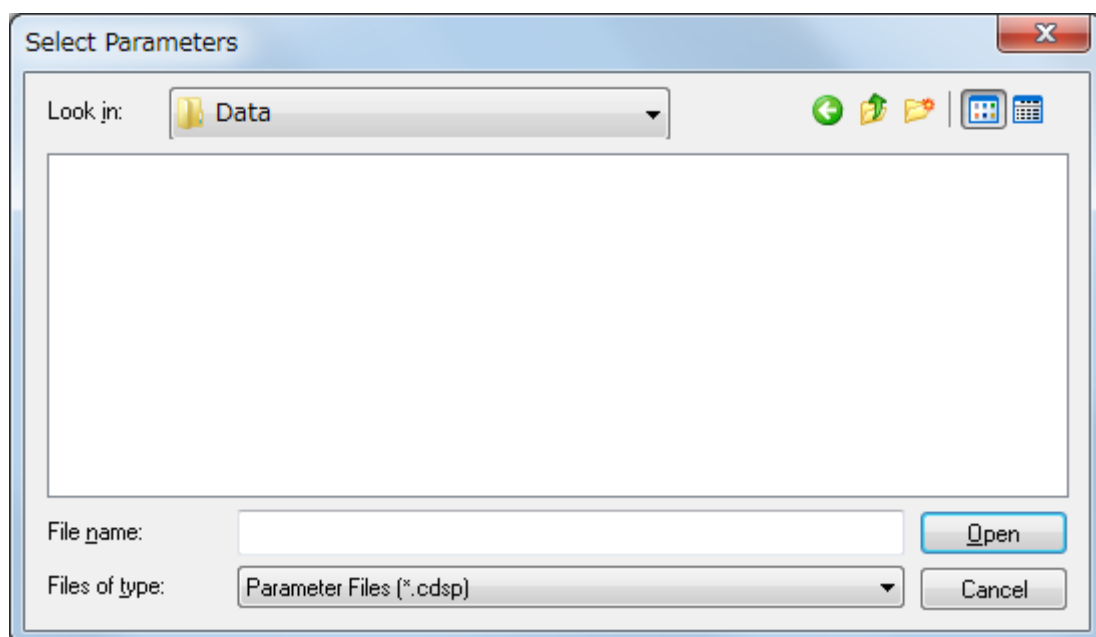


Figure 6.29 [Select Parameters] dialog box (when <Browse...> is clicked in [Default Parameters])

6.6.2 [Customize Toolbar...]

Customizes and shows/hides toolbars.

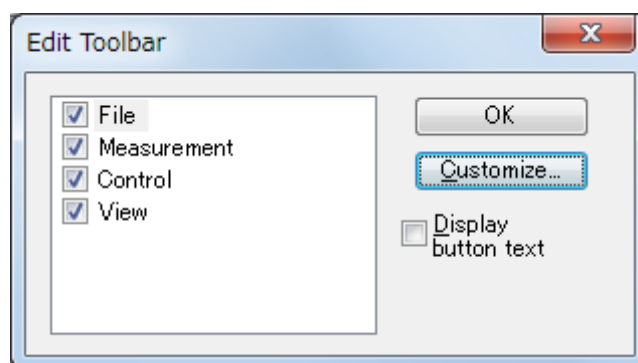




































































Figure 6.30 [Edit Toolbar] dialog box

[File]	<p>Shows/hides the toolbar corresponding to the [File] menu and adds/removes the following tool buttons.</p> <table border="1"> <thead> <tr> <th>Button</th><th>Name</th></tr> </thead> <tbody> <tr> <td></td><td>[Save]</td></tr> <tr> <td></td><td>[Send to Analysis]</td></tr> <tr> <td></td><td>[Open Parameters]</td></tr> <tr> <td></td><td>[Save Parameters]</td></tr> </tbody> </table>	Button	Name		[Save]		[Send to Analysis]		[Open Parameters]		[Save Parameters]
Button	Name										
	[Save]										
	[Send to Analysis]										
	[Open Parameters]										
	[Save Parameters]										

[Measurement]	<p>Shows/hides the toolbar corresponding to the [Measure] menu and adds/removes the following tool buttons.</p> <table border="1"> <thead> <tr> <th>Button</th><th>Name</th></tr> </thead> <tbody> <tr> <td></td><td>[Cancel]</td></tr> <tr> <td></td><td>[Sequence Measurement]</td></tr> <tr> <td></td><td>[Sample Measurement]</td></tr> <tr> <td></td><td>[Baseline Measurement]/ [Blank Measurement]</td></tr> <tr> <td></td><td>[Dark Measurement]</td></tr> <tr> <td></td><td>[Parameters]</td></tr> </tbody> </table>	Button	Name		[Cancel]		[Sequence Measurement]		[Sample Measurement]		[Baseline Measurement]/ [Blank Measurement]		[Dark Measurement]		[Parameters]		
Button	Name																
	[Cancel]																
	[Sequence Measurement]																
	[Sample Measurement]																
	[Baseline Measurement]/ [Blank Measurement]																
	[Dark Measurement]																
	[Parameters]																
[Control]	<p>Shows/hides the toolbar corresponding to the [Control] menu and adds/removes the following tool buttons.</p> <table border="1"> <thead> <tr> <th>Button</th><th>Name</th></tr> </thead> <tbody> <tr> <td></td><td>[Auto Zero]</td></tr> <tr> <td></td><td>[Shutter]</td></tr> <tr> <td></td><td>[Move Wavelength]</td></tr> <tr> <td></td><td>[Sensitivity]</td></tr> <tr> <td></td><td>[Light Source][Hg Lamp]</td></tr> </tbody> </table>	Button	Name		[Auto Zero]		[Shutter]		[Move Wavelength]		[Sensitivity]		[Light Source][Hg Lamp]				
Button	Name																
	[Auto Zero]																
	[Shutter]																
	[Move Wavelength]																
	[Sensitivity]																
	[Light Source][Hg Lamp]																
[View]	<p>Shows/hides the toolbar corresponding to the [View] menu and adds/removes the following tool buttons.</p> <table border="1"> <thead> <tr> <th>Button</th><th>Name</th></tr> </thead> <tbody> <tr> <td></td><td>[Scales]</td></tr> <tr> <td></td><td>[Patterns]</td></tr> <tr> <td></td><td>[Fonts]</td></tr> <tr> <td></td><td>[Gridlines]</td></tr> <tr> <td></td><td>[Styles]</td></tr> <tr> <td></td><td>[Digital]</td></tr> <tr> <td></td><td>[Graph]</td></tr> </tbody> </table>	Button	Name		[Scales]		[Patterns]		[Fonts]		[Gridlines]		[Styles]		[Digital]		[Graph]
Button	Name																
	[Scales]																
	[Patterns]																
	[Fonts]																
	[Gridlines]																
	[Styles]																
	[Digital]																
	[Graph]																
[Display button text]	Displays the button name under each tool button.																
<Customize...>	Displays the dialog box for customizing toolbars (see Figure 6.31).																

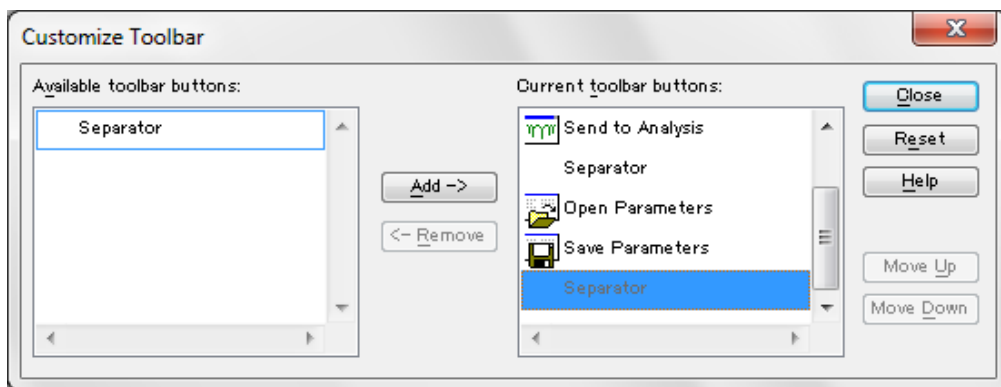


Figure 6.31 [Customize Toolbar] dialog box

6.6.3 [Monitor Bar...]

Customizes the style for the monitor bar.

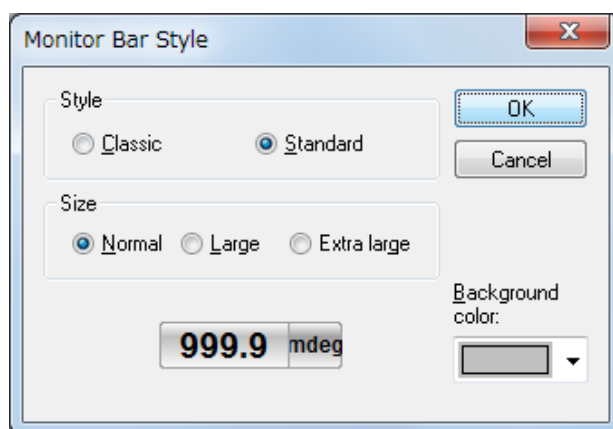
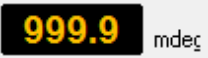
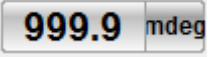


Figure 6.32 [Monitor Bar Style] dialog box

[Style]	<p>Selects the style for the monitor bar.</p> <p>[Classic]: </p> <p>[Standard]: </p>
[Size]	<p>Selects the size for the monitor bar.</p> <p>Size: [Normal], [Large], [Extra large]</p>
[Background color]	<p>Selects the background color for the monitor bar when [Standard] is selected for [Style].</p>

6.7 [Help] Menu

6.7.1 [Contents]

Displays the help window with the contents tab selected.

6.7.2 [Search Topic]

Displays the help window with the index tab selected.

6.7.3 [About...]

Displays version information for the program.

7. [Spectra Measurement] Program Reference


The [Spectra Measurement] program measures the change in the sample signal intensity while scanning the wavelength. This chapter describes specific measurement parameters for the [Spectra Measurement] program.

7.1 [Parameters...]

Specifies measurement parameters. The dialog box has four tabs for parameter settings, as shown in Figure 7.2 [General], Figure 7.3 [Control], Figure 7.4 [Information], and Figure 7.5 [Data]. The [General] tab is open by default in the [Parameters] dialog box. The page can be changed by clicking each tab at the top of the dialog box.

The parameters can be switched between two modes: Basic mode and Advanced mode. In Basic mode, the choice of parameters is partially limited. In Advanced mode, each parameter is specified individually.

<Basic Mode/Advanced Mode>	Switches between Basic and Advanced modes. Advanced mode can only be used for J-1500.
<Open...>	Opens a measurement parameter file.
<Save...>	Saves the measurement parameters. In the CFR version, an electronic signature is required for saving parameters.
<Default>	Restores the parameters of the [General] and [Control] tabs to their default values.
<OK>	Applies the changes to the [Parameters] settings and closes the dialog box.
<Cancel>	Closes the dialog box without applying any changes to the [Parameters] settings.

Note 1: This operation can also be performed by clicking the tool button  [Parameters].

Note 2: In the CFR version, Managers or Analysts authority is required to specify the parameters.

Note 3: In the CFR version, when the parameters are specified or changed, a signature is required. Clicking the <OK> button when the parameters have not been saved displays the following message.

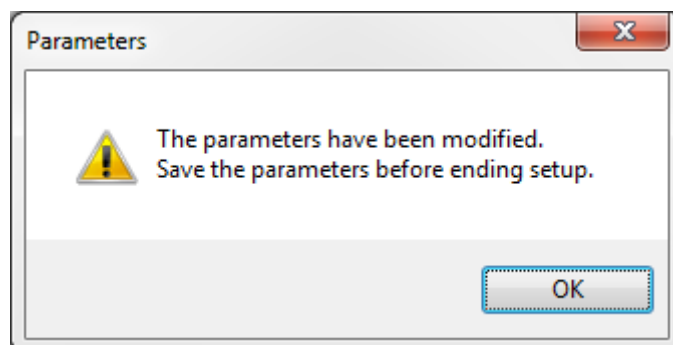


Figure 7.1 Message when the signature for the parameters is missing

7.1.1 [General] tab

The 'Parameters Advanced' dialog box, General tab, for J-1100/1500. The 'Basic Mode' button is highlighted with a blue dashed border.

Photometric mode

Channels Num: 2

Channel 1: CD

Channel 2: HT

Channel 3: CD

Channel 4: CD

Start: 300 nm

End: 200 nm

Data pitch: 0.1 nm

Start mode: Immediately

Scanning mode: Continuous

Scanning speed: 100 nm/min

CD scale: 200 mdeg/0.1 dOD

FL scale: 200 mdeg/0.1 dOD

D.I.T.: 1 sec

Bandwidth: 1.00 nm

☐ Slit width: 100 μ m

Accumulation/cycle

☒ Accumulation:

No. of accumulations: 1

Buttons: Basic Mode, Open..., Save..., Default, OK, Cancel

Figure 7.2 (a) J-1100/1500 [General] tab in [Parameters] dialog box

The 'Parameters Advanced' dialog box, General tab, for J-1700. The 'Basic Mode' button is highlighted with a blue dashed border.

Photometric mode

Channels Num: 2

Channel 1: CD

Channel 2: HT

Channel 3: CD

Channel 4: CD

Start: 300 nm

End: 200 nm

Data pitch: 0.1 nm

Start mode: Immediately

Scanning mode: Continuous

Scanning speed: 100 nm/min

CD scale: 200 mdeg/0.1 dOD

FL scale: 200 mdeg/0.1 dOD

D.I.T.: 1 sec

NIR D.I.T.: 1 sec

Bandwidth: 1.00 nm

NIR bandwidth: 20 nm

☐ Slit width: 100 μ m

☐ NIR slit width: 100 μ m

Accumulation/cycle

☒ Accumulation:

No. of accumulations: 1

Buttons: Basic Mode, Open..., Save..., Default, OK, Cancel

Figure 7.2 (b) J-1700 [General] tab in [Parameters] dialog box

Photometric mode	Selects the desired measurement parameters for the photometric mode.
[Channels Num]	Selects the number of channels to display in the spectrum display field. Click the arrow (▼ or ▲) to select the number of channels. Options: 1 to 4
[Channel <i>n</i>]	Selects the photometric mode for each channel. Click the drop-down list to select the photometric mode.
[CD scale]	<p>Selects the CD/LD range for the detector in the transmission configuration.</p> <p>[20 mdeg/0.05 dOD]: High sensitivity mode. Available when the CD range is within ± 20 mdeg and the LD range is within ± 0.05 dOD.</p> <p>[200 mdeg/1.0 dOD]: Standard sensitivity mode. Available when the CD range is within ± 200 mdeg and the LD range is within ± 1.0 dOD.</p> <p>[2000 mdeg/1.0 dOD]: Wide range mode. Available when the CD range is within ± 2000 mdeg and the LD range is within ± 1.0 dOD.</p> <p>[10000 mdeg/1.0 dOD]: Maximum range mode. Available when the CD range is within ± 10000 mdeg and the LD range is within ± 1.0 dOD.</p>
[FL scale]	<p>Selects the FD CD/FD LD range for the C detector in the fluorescence configuration.</p> <p>[20 mdeg/0.05 dOD]: High sensitivity mode. Available when the FD CD range is within ± 20 mdeg and the FD LD range is within ± 0.05 dOD.</p> <p>[200 mdeg/1.0 dOD]: Standard sensitivity mode. Available when the FD CD range is within ± 200 mdeg and the FD LD range is within ± 1.0 dOD.</p> <p>[2000 mdeg/1.0 dOD]: Wide range mode. Available when the FD CD range is within ± 2000 mdeg and the FD LD range is within ± 1.0 dOD.</p> <p>[10000 mdeg/1.0 dOD]: Maximum range mode. Available when the FD CD range is within ± 10000 mdeg and the FD LD range is within ± 1.0 dOD.</p> <p><i>Note: [FL scale] is not displayed on the J-1100.</i></p>
[D.I.T.]	<p>Selects the data acquisition time (integration time) per point.</p> <p>J-1100: 5 msec to 8 sec</p> <p>J-1500/1700: 0.1 msec to 30 sec</p>
[Bandwidth]	<p>Selects the spectral bandwidth.</p> <p>J-1100: 1 nm (fixed)</p> <p>J-1500/1700: 0.01 nm to 16.00 nm</p>

[Slit width]	<p>Specifies the slit width within the range from 1 to 4000 μm. Selecting [Slit width] activates the text box for entering the slit width.</p> <p><i>Note: [Slit width] is not displayed on the Basic Mode.</i></p>						
[Start]	<p>Specifies the measurement start wavelength.</p> <p>J-1100: 180 to 600 nm J-1500: 163 to 1600 nm J-1700: 163 to 2500 nm</p>						
[End]	<p>Specifies the measurement end wavelength.</p> <p>J-1100: 180 to 600 nm J-1500: 163 to 1600 nm J-1700: 163 to 2500 nm</p>						
[Data pitch]	<p>Selects the wavelength interval for data acquisition.</p> <p><i>Note: For J-1700, when the measurement start wavelength is 950 nm or more, set the data pitch to 0.1 nm or more.</i></p>						
[Start mode]	<p>Selects the measurement start condition.</p> <p>Options: [Immediately], [Wait for ext trigger] Use [Wait for ext. trigger] when an optional accessory such as a stopped flow system is used.</p>						
[Scanning mode]	<p>Selects the scanning mode.</p> <p>Options:</p> <table border="0"> <tr> <td data-bbox="550 1108 726 1153">[Continuous]:</td><td data-bbox="863 1108 1428 1187">Continuous scan. Scans the wavelength at the speed specified in [Scanning speed].</td></tr> <tr> <td data-bbox="550 1198 638 1243">[Step]:</td><td data-bbox="863 1198 1428 1568">Step scan. Increments the wavelength in steps corresponding to the data wavelength interval, and performs measurements at each wavelength for a time specified by [D.I.T.] while the drive is stopped. The scanning speed depends on [D.I.T.]. In this case, the D.I.T. value is chosen automatically and cannot be selected.</td></tr> <tr> <td data-bbox="550 1579 726 1624">[Step (Auto)]:</td><td data-bbox="863 1579 1428 1960">Performs a step scan while automatically adjusting D.I.T. to obtain a constant signal-to-noise ratio. Selecting [Step (Auto)] displays [D.I.T. Upper] and [D.I.T. Lower] instead of [D.I.T.]. Click the arrows to display the drop-down lists and select the desired values. In the low signal-to-noise ratio region, the D.I.T. is automatically set between [D.I.T. Upper] and [D.I.T. Lower].</td></tr> </table> <p><i>Note: [Step (Auto)] is not displayed on the Basic Mode.</i></p>	[Continuous]:	Continuous scan. Scans the wavelength at the speed specified in [Scanning speed].	[Step]:	Step scan. Increments the wavelength in steps corresponding to the data wavelength interval, and performs measurements at each wavelength for a time specified by [D.I.T.] while the drive is stopped. The scanning speed depends on [D.I.T.]. In this case, the D.I.T. value is chosen automatically and cannot be selected.	[Step (Auto)]:	Performs a step scan while automatically adjusting D.I.T. to obtain a constant signal-to-noise ratio. Selecting [Step (Auto)] displays [D.I.T. Upper] and [D.I.T. Lower] instead of [D.I.T.]. Click the arrows to display the drop-down lists and select the desired values. In the low signal-to-noise ratio region, the D.I.T. is automatically set between [D.I.T. Upper] and [D.I.T. Lower].
[Continuous]:	Continuous scan. Scans the wavelength at the speed specified in [Scanning speed].						
[Step]:	Step scan. Increments the wavelength in steps corresponding to the data wavelength interval, and performs measurements at each wavelength for a time specified by [D.I.T.] while the drive is stopped. The scanning speed depends on [D.I.T.]. In this case, the D.I.T. value is chosen automatically and cannot be selected.						
[Step (Auto)]:	Performs a step scan while automatically adjusting D.I.T. to obtain a constant signal-to-noise ratio. Selecting [Step (Auto)] displays [D.I.T. Upper] and [D.I.T. Lower] instead of [D.I.T.]. Click the arrows to display the drop-down lists and select the desired values. In the low signal-to-noise ratio region, the D.I.T. is automatically set between [D.I.T. Upper] and [D.I.T. Lower].						

[Scanning speed]	Specifies the scanning speed. Available if [Continuous] is selected for [Scanning mode]. J-1100: 1 nm/min to 5000 nm/min J-1500/1700: 1 nm/min to 10000 nm/min
[NIR D. I. T.]	Selects the data acquisition time (integration time) per point for NIR region. J-1700: 16 msec to 30 sec
[NIR bandwidth]	Selects the spectral bandwidth for NIR region. J-1700: 5nm to 50 nm
[NIR slit width]	Specifies the slit width within the range from 1 to 4000 μm . Selecting [NIR slit width] activates the text box for entering the slit width. (only for J-1700)
Accumulation/cycle	
[No. of cycles]	Specifies the number of measurements for each sample. When 2 or more measurements are specified, the [Cycle interval] text box is displayed. Input range: 1 to 999 <div style="border: 1px solid black; padding: 5px; margin-top: 10px;"><i>Note: When the [Accumulation] check box is selected, [No. of cycles] is not displayed.</i></div>
[Accumulation]	Specifies that cumulative averaging should be used to measure the spectrum. Selecting [Accumulation] displays [No. of accumulations].
[No. of accumulations]	Specifies the number of measurements used to average the spectrum. If measurement is stopped, all completed accumulations are averaged.

7.1.2 [Control] tab

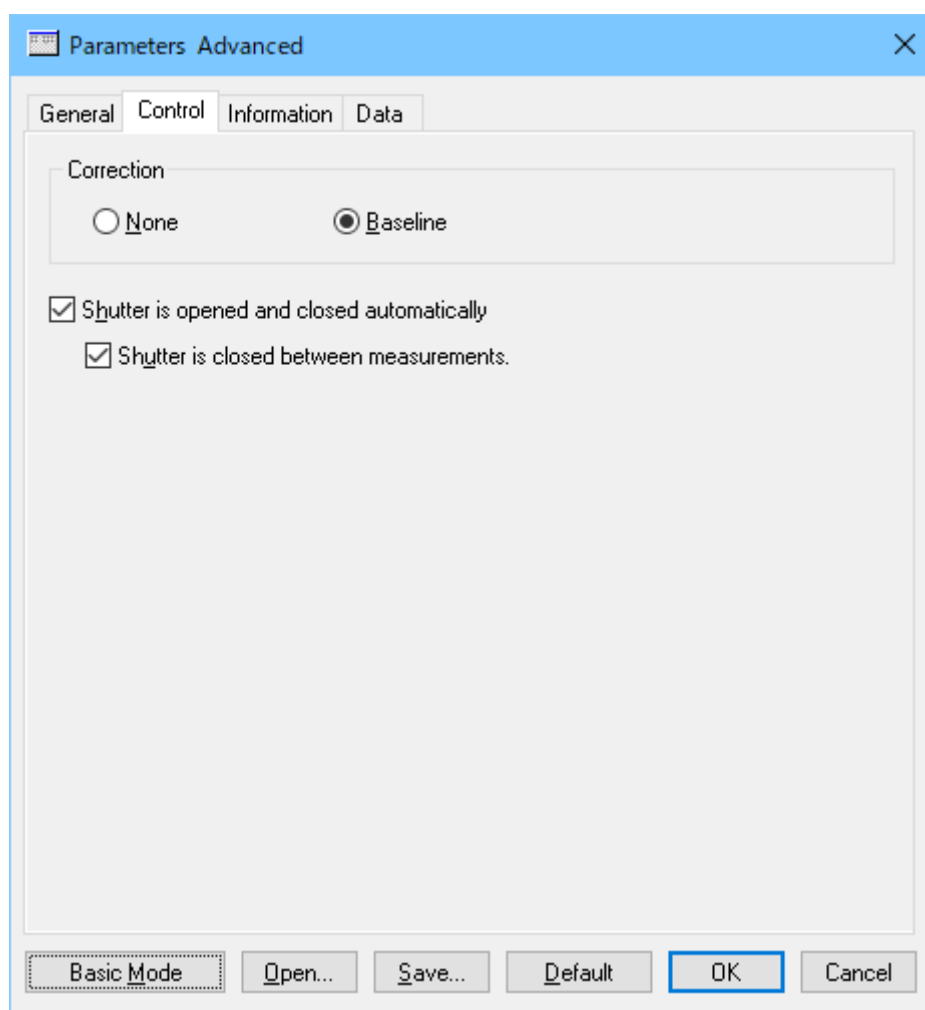


Figure 7.3 [Control] tab in [Parameters] dialog box

Correction	
[None]	Selects the use of no correction.
[Baseline]	Selects baseline correction using a measured baseline.
[Shutter is opened and closed automatically]	Automatically opens the shutter just before the measurement starts and closes it after the measurement finishes, regardless of the current shutter status. Even when this option is selected, if cycle measurement is selected, the shutter is not closed between measurements.
[Shutter is closed between measurements]	Automatically closes the shutter between measurements, if cycle measurement is selected.

7.1.3 [Information] tab

Parameters Advanced

General Control **Information** Data

Sample name:

Username:

Division:

Comment:

Concentration: mol/L

Solvent:

☐ Display the [Comments] dialog box before measurement

Basic Mode Open... Save... Default **OK** Cancel

Figure 7.4 [Information] tab in [Parameters] dialog box

[Sample name]	Enters a sample name (maximum 63 characters).
[Username]	Enters a username (maximum 63 characters).
[Division]	Enters a division name for the operator (maximum 127 characters).
[Comment]	Enters a comment (maximum 127 characters).
[Concentration]	Enters a sample concentration. There is no character limit.
[Solvent]	Enters a solvent name (maximum 127 single-byte characters).
[Display the [Comments] dialog box before measurement]	Displays the [Comments] dialog box for each measurement.

Note: In the CFR version, the [Username] and [Division] text boxes cannot be changed by the user.

7.1.4 [Data] tab

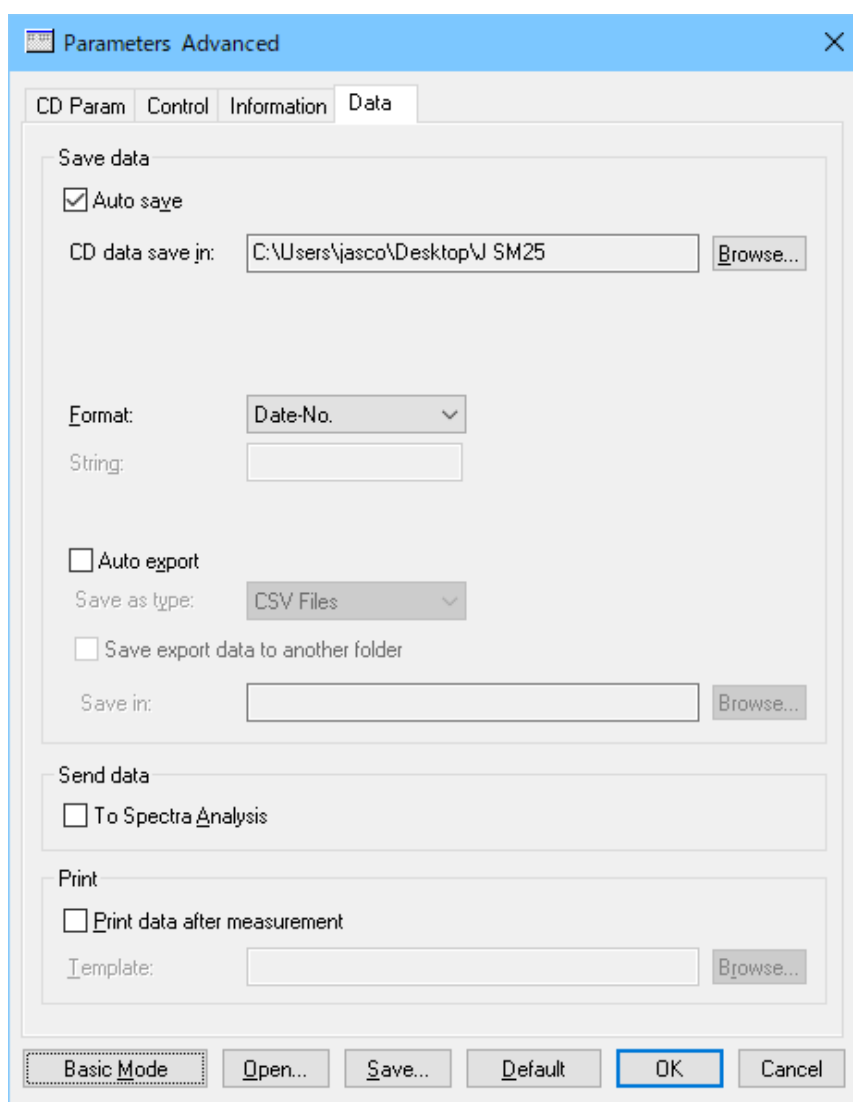


Figure 7.5 [Data] tab in [Parameters] dialog box

Save data

Selects to automatically save data. A file name is automatically allocated to the measured data and the data is saved to the specified directory.

Note: In the CFR version, when [Auto save] is selected, the following dialog box is displayed when the sample measurement starts. Start the measurement after typing a signature.

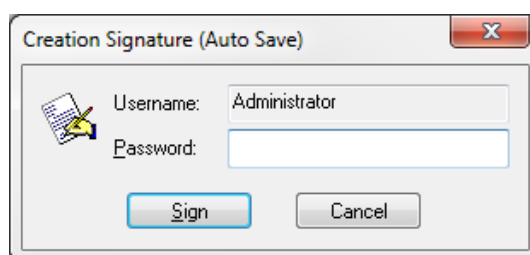


Figure 7.6 [Creation Signature (Auto Save)] dialog box

[Auto save]	Automatically saves measured data.
[CD data save in]	Displays the drive and folder name for saving data.
<Browse...>	Browses to the [CD data save in] drive and folder.
[Format]	Selects the file name to be assigned during auto saving. The file format is Standard Files (*.jws).
[Date-No.]	Sets [Date] + [3-digit serial number] as the file name. The date and serial number are connected with a hyphen "-".
[Sample-No.]	Sets [Sample name] + [3-digit serial number] as the file name. [Sample name] is the name entered in the [Information] tab.
[Comment-No.]	Sets [Comment] + [3 digit serial number] as the file name. [Comment] is the comment entered in the [Information] tab.
[Sample-Comment-No.]	Sets [Sample name] + [Comment] + [3-digit serial number] as the file name. [Sample name] and [Comment] are the name and comment entered in the [Information] tab, respectively.
[String-No.]	Sets [String-No.] + [3-digit serial number] as the file name. Type the character string in the text box below the drop-down list (maximum 63 single-byte characters).
[String]	Enters a character string when [String-No.] is selected for [Format] (maximum 63 single-byte characters).
[Auto export]	Exports the measured data.
[Save as type]	Selects the file type for the exported data. Options: CSV Files, ASCII Files
[Save export data to another folder]	Selects to save the exported data to a separate folder.
[Save in]	Displays the drive and folder name for saving exported data. <div style="border: 1px solid black; padding: 5px; margin-top: 5px;"><i>Note: In the CFR version, export data can be saved in a folder other than the workgroup folder.</i></div>
<Browse...>	Browses to the [Save in] drive and folder.
Send data	Selects to automatically send measured data to the [Spectra Analysis] program after measurement.
[To Spectra Analysis]	Automatically sends measured data to the [Spectra Analysis] program after measurement.
Print	Selects to automatically print data after measurement.
[Print data after measurement]	Automatically prints data after measurement.
[Template]	Displays the drive and file name of a saved template file.
<Browse...>	Selects the drive and file name of a saved template file.

8. [Time Course Measurement] Program Reference

The [Time Course Measurement] program measures the temporal change in the sample signal intensity at a fixed wavelength. This chapter describes specific measurement parameters for the [Time Course Measurement] program.


8.1 [Parameters...]

Specifies the measurement parameters.

The window has four tabs for parameter settings, as shown in Figure 8.2 [General], Figure 8.3 [Control], Figure 8.4 [Information], and Figure 8.5 [Data]. The [General] tab is open by default in the [Parameters] dialog box. The page can be changed by clicking each tab at the top of the dialog box.

The parameters can be switched between two modes: Basic mode and Advanced mode. In Basic mode, the choice of parameters is partially limited. In Advanced mode, each parameter is specified individually.

<Basic Mode/Advanced Mode>	Switches between Basic and Advanced modes. Advanced mode can only be used for J-1500.
<Open...>	Opens a measurement parameter file.
<Save...>	Saves the measurement parameters. In the CFR version, an electronic signature is required for saving parameters.
<Default>	Restores the parameters of the [General] and [Control] tabs to their default values.
<OK>	Applies the changes to the [Parameters] settings and closes the dialog box.
<Cancel>	Closes the dialog box without applying any changes to the [Parameters] settings.

Note 1: This operation can also be performed by clicking the tool button  [Parameters].

Note 2: In the CFR version, Managers or Analysts authority is required to specify the parameters.

Note 3: In the CFR version, when the parameters are specified or changed, a signature is required. Clicking the <OK> button when the parameters have not been saved displays the following message.

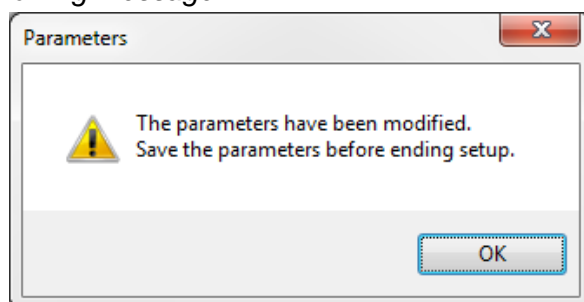


Figure 8.1 Message when the signature for the parameters is missing

8.1.1 [General] tab

The screenshot shows the 'Parameters Advanced' dialog box with the 'General' tab selected. The 'Photometric mode' section includes 'Channels Num' set to 2, and 'Channel 1' through 'Channel 4' all set to 'CD'. The 'Wavelength' is 253.7 nm, 'Measuring time' is 60 sec, 'Data pitch' is 1.0 sec, and 'Start mode' is 'Immediately'. The 'CD scale' and 'FL scale' are both set to '200 mdeg/0.1 dOD'. 'D.I.T.' is 1 sec, 'Bandwidth' is 1.00 nm, and 'Slit width' is 100 um. The 'Accumulation/cycle' section has 'Accumulation' checked and 'No. of accumulations' set to 1. At the bottom, the 'Basic Mode' button is highlighted with a blue border.

Parameters Advanced

General Control Information Data

Photometric mode

Channels Num: 2

Channel 1: CD

Channel 2: HT

Channel 3: CD

Channel 4: CD

Wavelength: 253.7 nm

Measuring time: 60 sec

Data pitch: 1.0 sec

Start mode: Immediately

CD scale: 200 mdeg/0.1 dOD

FL scale: 200 mdeg/0.1 dOD

D.I.T.: 1 sec

Bandwidth: 1.00 nm

☐ Slit width: 100 um

Accumulation/cycle

☒ Accumulation:

No. of accumulations: 1

Basic Mode Open... Save... Default OK Cancel

Figure 8.2 (a) J-1100/1500 [General] tab in [Parameters] dialog box

The screenshot shows the 'Parameters Advanced' dialog box with the 'General' tab selected. The 'Photometric mode' section includes 'Channels Num' set to 2, and 'Channel 1' through 'Channel 4' all set to 'CD'. The 'Wavelength' is 253.7 nm, 'Measuring time' is 60 sec, 'Data pitch' is 1.0 sec, and 'Start mode' is 'Immediately'. The 'CD scale' and 'FL scale' are both set to '200 mdeg/0.1 dOD'. 'D.I.T.' is 1 sec, 'Bandwidth' is 1.00 nm, and 'Slit width' is 100 um. The 'NIR D.I.T.' is 1 sec, 'NIR bandwidth' is 20 nm, and 'NIR slit width' is 100 um. The 'Accumulation/cycle' section has 'Accumulation' checked and 'No. of accumulations' set to 1. At the bottom, the 'Basic Mode' button is highlighted with a blue border.

Parameters Advanced

General Control Information Data

Photometric mode

Channels Num: 2

Channel 1: CD

Channel 2: HT

Channel 3: CD

Channel 4: CD

Wavelength: 253.7 nm

Measuring time: 60 sec

Data pitch: 1.0 sec

Start mode: Immediately

CD scale: 200 mdeg/0.1 dOD

FL scale: 200 mdeg/0.1 dOD

D.I.T.: 1 sec

Bandwidth: 1.00 nm

☐ Slit width: 100 um

NIR D.I.T.: 1 sec

NIR bandwidth: 20 nm

☐ NIR slit width: 100 um

Accumulation/cycle

☒ Accumulation:

No. of accumulations: 1

Basic Mode Open... Save... Default OK Cancel

Figure 8.2 (b) J-1700 [General] tab in [Parameters] dialog box

Photometric mode	Selects the desired measurement parameters for the photometric mode.
[Channels Num]	Click the arrow (▼ or ▲) and select [Channels Num:] to display in the time-course data display field. Options: 1 to 4
[Channel <i>n</i>]	Selects the photometric mode for each channel. Click the drop-down list to select the photometric mode.
[CD scale]	<p>Selects the CD/LD scale range for the detector in the transmission configuration.</p> <p>[20 mdeg/0.05 dOD]: High sensitivity mode. Available when the CD range is within ± 20 mdeg and the LD range is within ± 0.05 dOD.</p> <p>[200 mdeg/1.0 dOD]: Standard sensitivity mode. Available when the CD range is within ± 200 mdeg and the LD range is within ± 1.0 dOD.</p> <p>[2000 mdeg/1.0 dOD]: Wide range mode. Available when the CD range is within ± 2000 mdeg and the LD range is within ± 1.0 dOD.</p> <p>[10000 mdeg/1.0 dOD]: Maximum range mode. Available when the CD range is within ± 10000 mdeg and the LD range is within ± 1.0 dOD.</p>
[FL scale]	<p>Selects the FD CD/FD LD scale range for the C detector in the fluorescence configuration.</p> <p>[20 mdeg/0.05 dOD]: High sensitivity mode. Available when the FD CD range is within ± 10 mdeg and the FD LD range is within ± 0.05 dOD.</p> <p>[200 mdeg/1.0 dOD]: Standard sensitivity mode. Available when the FD CD range is within ± 200 mdeg and the FD LD range is within ± 1.0 dOD.</p> <p>[2000 mdeg/1.0 dOD]: Wide range mode. Available when the FD CD range is within ± 2000 mdeg and the FD LD range is within ± 1.0 dOD.</p> <p>[10000 mdeg/1.0 dOD]: Maximum range mode. Available when the FD CD range is within ± 10000 mdeg and the FD LD range is within ± 1.0 dOD.</p> <p><i>Note: [FL scale] is not displayed on the J-1100.</i></p>
[D.I.T.]	<p>Selects the data acquisition time (integration time) per point.</p> <p>J-1100: 5 msec to 8 sec</p> <p>J-1500/1700: 0.1 msec to 30 sec</p>
[Bandwidth]	<p>Selects the spectral bandwidth.</p> <p>J-1100: 1 nm (fixed)</p> <p>J-1500/1700: 0.01 nm to 16.00 nm</p>

[Slit width]	<p>Specifies the slit width within the range from 1 to 4000 μm. Selecting [Slit width] activates the text box for entering the slit width.</p> <p><i>Note: [Slit width] is not displayed on the Basic Mode.</i></p>
[Wavelength]	<p>Specifies the measurement start wavelength.</p> <p>J-1100: 180 to 600 nm J-1500: 163 to 1600 nm J-1700: 163 to 2500 nm</p>
[Measuring time]	<p>Specifies the measurement time.</p> <p>Input range: 2 to 20002 sec</p>
[Data pitch]	<p>Selects the measurement interval for data acquisition.</p> <p>J-1100: 5 msec to 60 min J-1500/1700: 0.1 msec to 60 min</p>
[Start mode]	<p>Selects the measurement start condition.</p> <p>Options: [Immediately], [Wait for ext trigger] Use [Wait for ext. trigger] when an optional accessory such as a stopped flow system is used.</p>
[NIR D. I. T.]	<p>Selects the data acquisition time (integration time) per point for NIR region.</p> <p>J-1700: 16 msec to 30 sec</p>
[NIR bandwidth]	<p>Selects the spectral bandwidth for NIR region.</p> <p>J-1700: 5nm to 50 nm</p>
[NIR slit width]	<p>Specifies the slit width within the range from 1 to 4000 μm. Selecting [Slit width] activates the text box for entering the slit width. (only for J-1700)</p> <p><i>Note: [NIR slit width] is not displayed on the Basic Mode.</i></p>
Accumulation/cycle	
[No. of cycles]	<p>Specifies the number of measurements for each sample. When 2 or more measurements are specified, the [Cycle interval] text box is displayed.</p> <p>Input range: 1 to 999</p> <p><i>Note: When the [Accumulation] check box is selected, [No. of cycles] is not displayed.</i></p>
[Accumulation]	<p>Specifies that cumulative averaging should be used to measure the spectrum. Selecting [Accumulation] displays [No. of accumulations].</p>
[No. of accumulations]	<p>Specifies the number of measurements used to average the spectrum. If measurement is stopped, all completed accumulations are averaged.</p> <p>Input range: 1 to 999.</p>

Note: "Accumulation/cycle" is not displayed on the Basic Mode.

8.1.2 [Control] tab

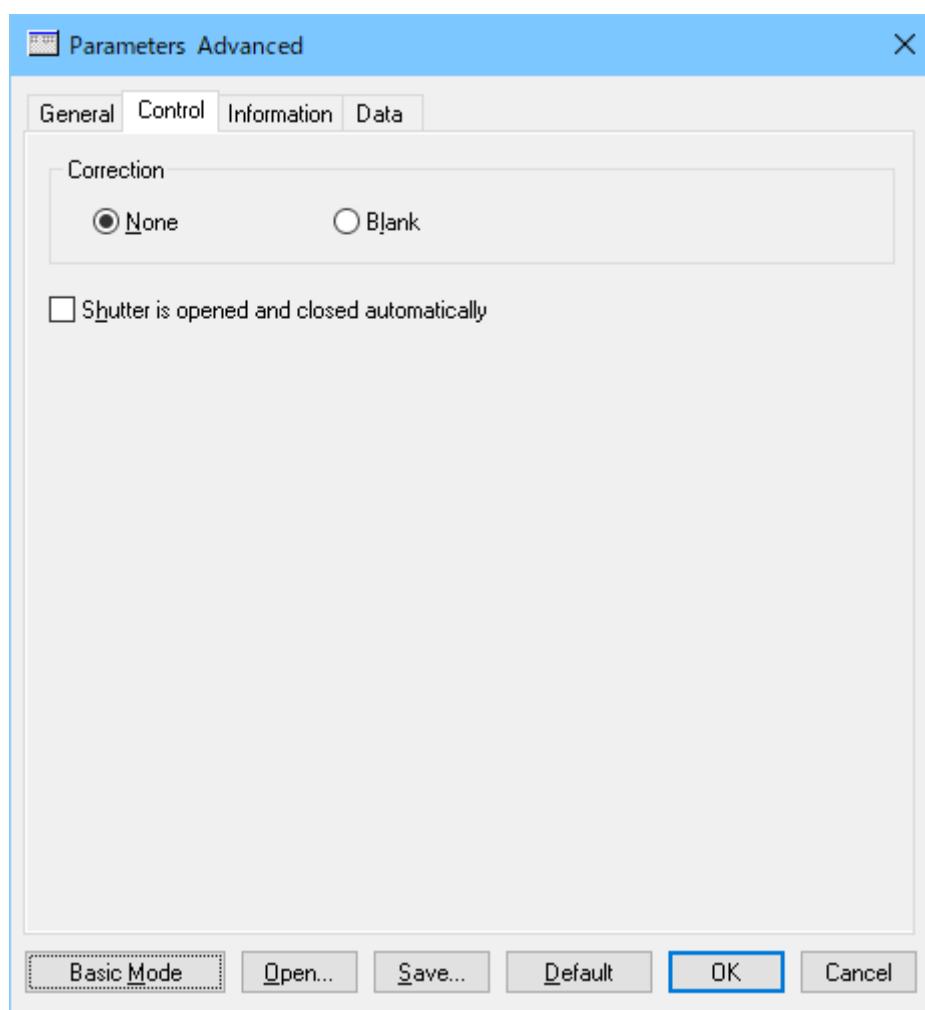


Figure 8.3 [Control] tab in [Parameters] dialog box

Correction	
[None]	Selects the use of no correction.
[Blank]	Selects blank correction using a measured blank.
[Shutter is opened and closed automatically]	Automatically opens the shutter just before the measurement starts and closes it after the measurement finishes, regardless of the current shutter status. Even when this option is selected, if cycle measurement is selected, the shutter is not closed between measurements.

8.1.3 [Information] tab

Parameters Advanced

General Control **Information** Data

Sample name:

Username:

Division:

Comment:

Concentration: mol/L

Solvent:

☐ Display the [Comments] dialog box before measurement

Basic Mode Open... Save... Default **OK** Cancel

Figure 8.4 [Information] tab in [Parameters] dialog box

[Sample name]	Enters a sample name (maximum 63 characters).
[Username]	Enters a username (maximum 63 characters).
[Division]	Enters a division name for the operator (maximum 127 characters).
[Comment]	Enters a comment (maximum 127 characters).
[Concentration]	Enters a sample concentration. There is no character limit.
[Solvent]	Enters a solvent name (maximum 127 single-byte characters).
[Display the [Comments] dialog box before measurement]	Displays the [Comments] dialog box for each measurement.

Note: In the CFR version, the [Username] and [Division] text boxes cannot be changed by the user.

8.1.4 [Data] tab

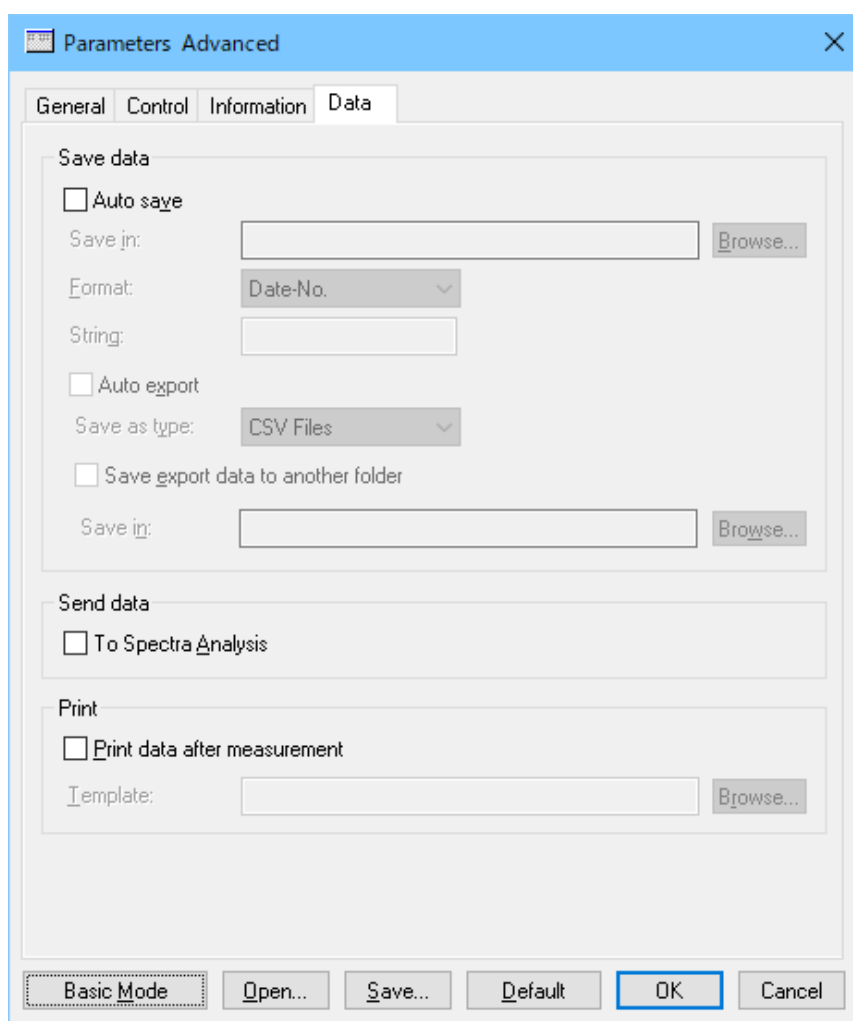


Figure 8.5 [Data] tab in [Parameters] dialog box

Save data

Selects to automatically save data. A file name is automatically allocated to the measured data and the data is saved to the specified directory.

Note: In the CFR version, when [Auto save] is selected, the following dialog box is displayed when the sample measurement starts. Start the measurement after typing a signature.

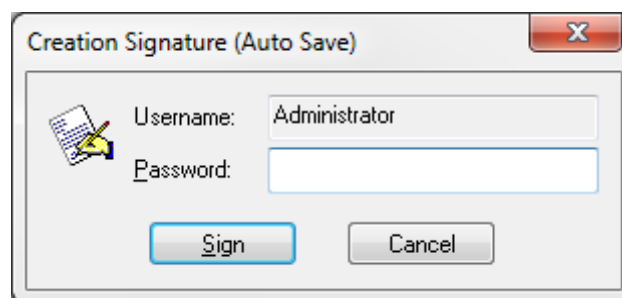


Figure 8.6 [Creation Signature (Auto Save)] dialog box

[Auto save]	Automatically saves measured data
[Save in]	Displays the drive and folder name for saving data.
<Browse...>	Browses to the [Save in] drive and folder.
[Format]	Selects the file name to be assigned during auto saving. The file format is Standard Files (*.jws).
[Date-No.]	Sets [Date] + [3-digit serial number] as the file name. The date and serial number are connected with a hyphen "-".
[Sample-No.]	Sets [Sample name] + [3-digit serial number] as the file name. [Sample name] is the name entered in the [Information] tab.
[Comment-No.]	Sets [Comment] + [3 digit serial number] as the file name. [Comment] is the comment entered in the [Information] tab.
[Sample-Comment-No.]	Sets [Sample name] + [Comment] + [3-digit serial number] as the file name. [Sample name] and [Comment] are the name and comment entered in the [Information] tab, respectively.
[String-No.]	Sets [String-No.] + [3-digit serial number] as the file name. Type the character string in the text box below the drop-down list (maximum 63 single-byte characters).
[String]	Enters a character string when [String-No.] is selected for [Format] (maximum 63 single-byte characters).
[Auto export]	Exports the measured data.
[Save as type]	Selects the file type for the exported data. Options: CSV Files, ASCII Files
[Save export data to another folder]	Selects to save the exported data to a separate folder.
[Save in]	Displays the drive and folder name for saving exported data. <div style="border: 1px solid black; padding: 5px; margin-top: 5px;"><i>Note: In the CFR version, export data can be saved in a folder other than the workgroup folder.</i></div>
<Browse...>	Browses to the [Save in] drive and folder.
Send data	Selects to automatically send measured data to [Spectra Analysis] program after measurement.
[To Spectra Analysis]	Automatically sends measured data to the [Spectra Analysis] program after measurement.
Print	Selects to automatically print data after measurement.
[Print data after measurement]	Automatically prints data after measurement.
[Template]	Displays the drive and file name of a saved template file.
<Browse...>	Selects the drive and file name of a saved template file.

9. [Data Monitor] Program Reference

The [Data Monitor] program is used to provide an enlarged display of the photometric values when the CD spectrometer is being tested or calibrated. This chapter describes specific measurement parameters for the [Data Monitor] program.

Note: In the CFR version, [Data Monitor] program is not available.

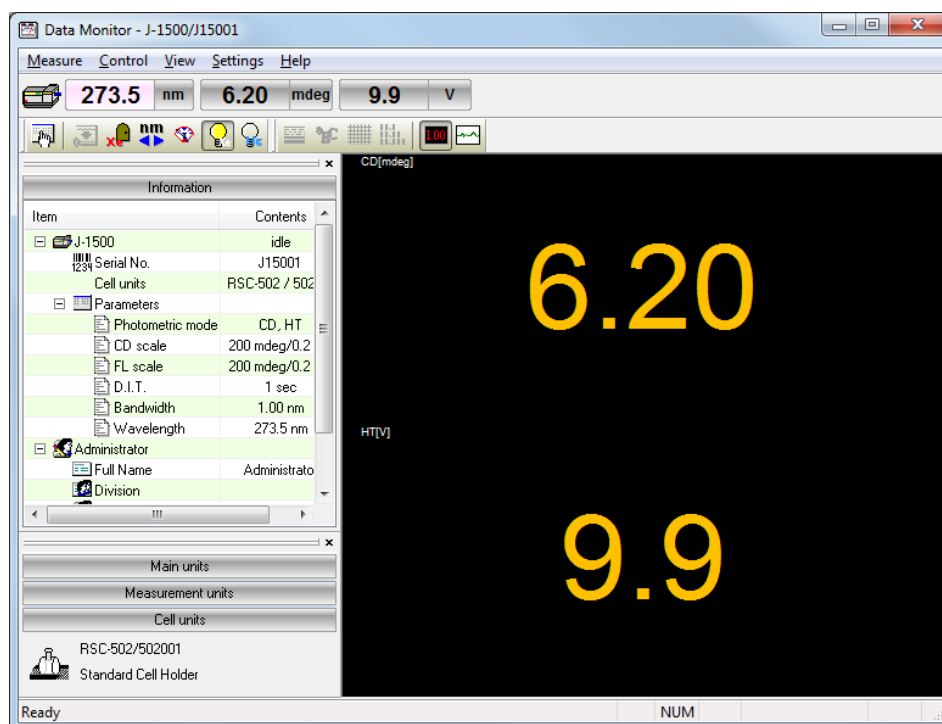


Figure 9.1 [Data Monitor] digital-mode display

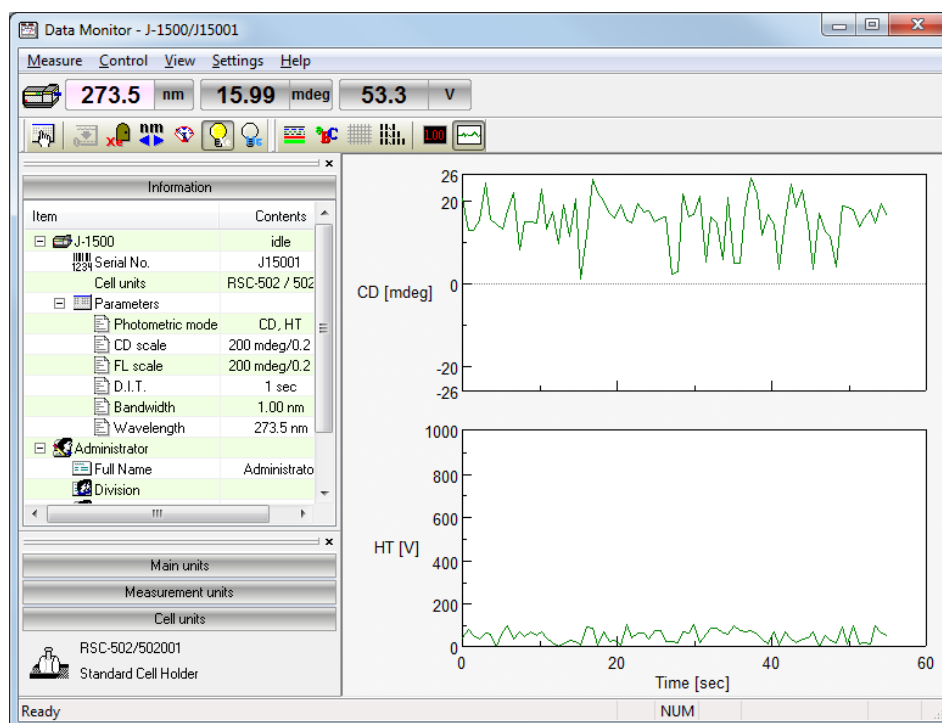



Figure 9.2 [Data Monitor] graph-mode display


9.1 [Parameters...]

Specifies the measurement parameters.

<OK>	Applies the changes to the [Parameters] settings and closes the dialog box.
<Cancel>	Closes the dialog box without applying any changes to the [Parameters] settings.


Note: This operation can also be performed by clicking the tool button  [Parameters].

9.1.1 [Parameters]

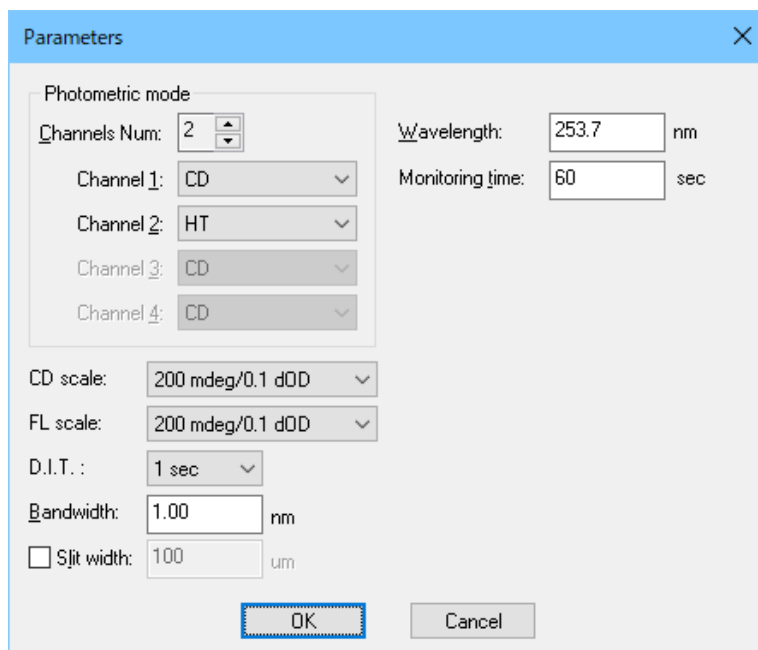

The Parameters dialog box for the J-1100/1500 model. It features a 'Photometric mode' section with 'Channels Num' set to 2. Below this are four channels: Channel 1 (CD), Channel 2 (HT), Channel 3 (CD), and Channel 4 (CD). To the right, 'Wavelength' is set to 253.7 nm and 'Monitoring time' is set to 60 sec. Further down, 'CD scale' and 'FL scale' are both set to 200 mdeg/0.1 dOD. 'D.I.T.' is set to 1 sec. 'Bandwidth' is set to 1.00 nm. There is an unchecked checkbox for 'Slit width' set to 100 um. At the bottom are 'OK' and 'Cancel' buttons.

Figure 9.3 (a) J-1100/1500 [General] tab in [Parameters] dialog box

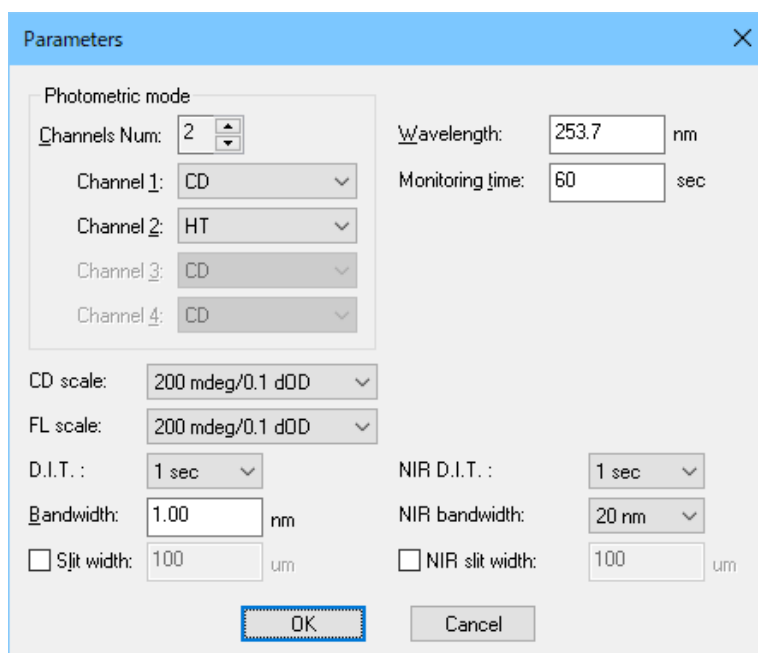

The Parameters dialog box for the J-1700 model. It includes the same 'Photometric mode' section as Figure 9.3 (a). In addition, it has 'NIR D.I.T.' set to 1 sec, 'NIR bandwidth' set to 20 nm, and an unchecked checkbox for 'NIR slit width' set to 100 um. The 'OK' and 'Cancel' buttons are at the bottom.

Figure 9.3 (b) J-1700 [General] tab in [Parameters] dialog box

Photometric mode	Selects the desired measurement parameters for the photometric mode.
[Channels Num]	Click the arrow (▼ or ▲) and select [Channels Num:] to display in the time-course data display field. Options: 1 to 4
[Channel <i>n</i>]	Selects the photometric mode for each channel. Click the drop-down list to select the photometric mode.
[CD scale]	<p>Selects the CD/LD scale range for the detector in the transmission configuration.</p> <p>[20 mdeg/0.05 dOD]: High sensitivity mode. Available when the CD range is within ± 20 mdeg and the LD range is within ± 0.05 dOD.</p> <p>[200 mdeg/1.0 dOD]: Standard sensitivity mode. Available when the CD range is within ± 200 mdeg and the LD range is within ± 1.0 dOD.</p> <p>[2000 mdeg/1.0 dOD]: Wide range mode. Available when the CD range is within ± 2000 mdeg and the LD range is within ± 1.0 dOD.</p> <p>[10000 mdeg/1.0 dOD]: Maximum range mode. Available when the CD range is within ± 10000 mdeg and the LD range is within ± 1.0 dOD.</p>
[FL scale]	<p>Selects the FDCD/FDLD scale range for the C detector in the fluorescence configuration.</p> <p>[20 mdeg/0.05 dOD]: High sensitivity mode. Available when the FDCD range is within ± 20 mdeg and the FDLD range is within ± 0.05 dOD.</p> <p>[200 mdeg/1.0 dOD]: Standard sensitivity mode. Available when the FDCD range is within ± 200 mdeg and the FDLD range is within ± 1.0 dOD.</p> <p>[2000 mdeg/1.0 dOD]: Wide range mode. Available when the FDCD range is within ± 2000 mdeg and the FDLD range is within ± 1.0 dOD.</p> <p>[10000 mdeg/1.0 dOD]: Maximum range mode. Available when the FDCD range is within ± 10000 mdeg and the FDLD range is within ± 1.0 dOD.</p> <div style="border: 1px solid black; padding: 2px; margin-top: 10px;"> <p><i>Note: [FL scale] is not displayed on the J-1100.</i></p> </div>
[D.I.T.]	<p>Selects the data acquisition time (integration time) per point.</p> <p>J-1100: 5 msec to 8 sec</p> <p>J-1500/1700: 0.1 msec to 30 sec</p>

[Bandwidth]	Selects the spectral bandwidth. J-1100: 1 nm (fixed) J-1500/1700: 0.01 nm to 16.00 nm
[Slit width]	Specifies the slit width within the range from 1 to 4000 μm . Selecting [Slit width] activates the text box for entering the slit width.
[Wavelength]	Specifies the measurement start wavelength. J-1100: 180 to 600 nm J-1500: 163 to 1600 nm J-1700: 163 to 2500 nm
[Monitoring time]	Specifies the monitoring time when graph mode is selected. Input range: 10 to 30000 sec
[NIR D. I. T.]	Selects the data acquisition time (integration time) per point for NIR region. J-1700: 16 msec to 30 sec
[NIR bandwidth]	Selects the spectral bandwidth for NIR region. J-1700: 5 nm to 50 nm
[NIR slit width]	Specifies the slit width within the range from 1 to 4000 μm . Selecting [Slit width] activates the text box for entering the slit width. (Only J-1700)

10. Appendix

10.1 Setting the Measurement Parameters

The parameters selected can seriously affect measurement results. The following summarizes how to select parameters.

10.1.1 Measurement Wavelength Range

Most CD spectrum peaks appear in a specific wavelength region (short wavelength region as in the case of ultraviolet-visible absorption spectra for most organic compounds). If the optically active absorption band position for the sample is known, take measurements beginning at a wavelength of 50 to 100 nm before the leading edge of the spectrum and ending at the trailing edge of the spectrum where it converges at zero, although this method varies slightly depending on the width and shape of the spectrum.

10.1.2 Data Acquisition Interval (Resolution)

10.1.2.1 Spectra measurement

Usually, a value of about 0.2 nm is sufficient. However, if the scanning mode is set to [Step], approximately 1 nm is more appropriate. Make the resolution coarse if the spectrum peaks are broad and fine if they are sharp.

10.1.2.2 Time-course measurement

Make the resolution fine if rapid changes with time are expected.

10.1.3 Scanning Speed

The scanning speed depends on the D. I. T. Under normal conditions, set the scanning speed at 20 to 100 nm/min and D. I. T. at 0.25 to 2 sec. Table 10.1 shows suggested D.I.T. values for different scanning speeds.

Table 10.1 D.I.T. values for different scanning speeds

Scanning speed (nm/min)	D.I.T. (recommended)	Upper limit
10000	1.0 to 2.0 msec	32.0 msec
5000	2.0 to 4.0 msec	64.0 msec
2000	4.0 to 8.0 msec	0.25 sec
1000	32.0 msec to 0.125 sec	0.5 sec
500	64.0 msec to 0.25 sec	1.0 sec
200	0.125 to 0.5 sec	2.0 sec
100	0.25 to 1.0 sec	4.0 sec
50	0.5 to 2.0 sec	8.0 sec
20	1.0 to 2.0 sec	16.0 sec
10	2.0 to 4.0 sec	16.0 sec
5	2.0 to 8.0 sec	16.0 sec
2	4.0 to 16.0 sec	16.0 sec
1	8.0 to 16.0 sec	16.0 sec

10.1.4 Number of Accumulations

The following relationship exists between the signal-to-noise ratio and the number of accumulations: $S/N \propto \sqrt{\text{Number of accumulations}}$.

Increasing the number of accumulations improves the signal-to-noise ratio for a spectrum. Determine the appropriate number of accumulations based on the spectrum (sample) condition and the overall measurement time (increasing the number of accumulations prolongs the measurement time).

If the sensitivity is lower than Standard, a single accumulation will suffice. Table 10.2 shows the suggested number of accumulations when a sample with an absorbance of 1 or less is measured with a D.I.T. of 2 sec.

Table 10.2 No. of accumulations and sensitivity (D.I.T. 2 sec)

Sensitivity	Number of accumulations
High	4 to 50
Standard, Low	1

Note: For time-course measurements, set the number of accumulations to 1. Cumulative measurement can be used in stopped flow measurement (see the instruction manual for the stopped flow unit).

10.1.5 Response (D.I.T.)

10.1.5.1 Spectra measurement

The following relationship exists between D.I.T. and the signal-to-noise ratio:

$$S/N \propto \sqrt{\text{Response}}.$$

The signal-to-noise ratio for a spectrum measured for 4 seconds is twice that for a spectrum measured for 1 second.

The spectral value at any given point in time is the average of values sampled before and after that time (D.I.T./2). Therefore, increasing D.I.T. leads to a greater effect of previous data on the current data. Do not increase D.I.T. excessively if scanning is carried out at high speed or if the resolution is low.

Table 10.3 shows suggested D.I.T. values for obtaining a spectrum with a high signal-to-noise ratio at a desired sensitivity. If the sensitivity is set at 10 mdeg and D. I. T. is set at 1 to 2 sec to improve the signal-to-noise ratio, increasing the number of accumulations is very effective because the spectrum is simultaneously averaged.

Table 10.3 D.I.T. values for obtaining high signal-to-noise ratio

CD scale (mdeg)	D.I.T. (sec)
10	0.5 to 16.0
200	0.125 to 0.5
2000	0.025 to 0.064

Note: A D.I.T. of 0.5 to 125 msec should be selected for high speed measurement using a stopped flow unit.

The above table also applies to a step scan. In this case, the measurement time is the sum of the wavelength stepping time and the data acquisition time at each step. If the D.I.T. is on the order of seconds, the number of accumulations will dominate the measurement time and the measurement time will be almost proportional to D.I.T. For a total wavelength measurement range of W (nm) and a data interval of Step (nm), the following relation holds between the total integration time and D.I.T.

$$\text{Total Integration Time} = (W/\text{Step} + 1) \times \text{D.I.T.}$$

(W/Step + 1): Number of data points

10.1.5.2 Time-course measurement

For time-course measurement, select a D.I.T. that is appropriate given the decay half-life of the CD signal due to a reaction. For normal reactions, set it to 1/10 of the decay half-life of the CD signal or smaller. For example, if the decay half-life is 20 msec, a D.I.T. of 0.5 to 2 msec is appropriate. If the decay is very slow, select a D.I.T. of 1 to 16 sec, taking into account the rate of change of the signal and the signal-to-noise ratio. When used as a monitor for HPLC, a D.I.T. of about 1 sec is suitable.

10.1.6 Bandwidth and slit width

In most cases, set the measurement bandwidth to 1.0 nm. Decrease it when measuring a sharp spectrum such as a MCD (magnetic circular dichroism) spectrum, or increase when performing measurements using a micro cell, when the signal-to-noise ratio is low. When a longer wavelength detector (400 to 1100 nm) is used, increase the bandwidth and set the slit width to about 60 μm to prevent the photomultiplier voltage (HT) from rising abnormally. However, ensure that the HT voltage does not become lower than 170 V.

Note 1: If the slit width is increased until the HT voltage drops below 170 V, the photomultiplier may be damaged.

Note 2: Most organic substances emit fluorescence. Therefore, measuring a sample having an absorbance of 2 or more at a bandwidth of 2 nm or more may distort the spectrum. Also, photolysis of a sample can occur in the ultraviolet region. If the sample is susceptible to photolysis, reduce the bandwidth to 0.5 nm or less.

JASCO Corporation
2967-5, Ishikawa-machi, Hachioji
Tokyo 192-8537, Japan

Printed in Japan